


Apr 10, 2023

High Spatial Resolution MALDI Imaging Mass Spectrometry Data Acquisition

DOI

<https://dx.doi.org/10.17504/protocols.io.n92ldpq99l5b/v1>

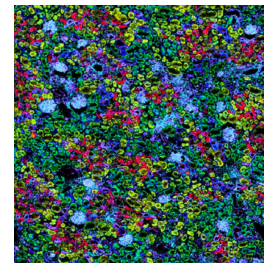

Katerina Djambazova¹, Angela Kruse¹, Jamie Allen¹, allison.b.esselman¹, Madeline E. Colley¹, David Anderson¹, ali.zahraei¹, Olof Isberg¹, Melissa Farrow¹, Jeff Spraggins¹, Martin Dufrense¹

¹Vanderbilt University

VU Biomolecular Multimodal Imaging Center / Spraggins Research Group
Tech. support email: jeff.spraggins@vanderbilte.du



Katerina V Djambazova

Vanderbilt University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ldpq99l5b/v1>



Protocol Citation: Katerina Djambazova, Angela Kruse, Jamie Allen, allison.b.esselman , Madeline E. Colley, David Anderson, ali.zahraei , Olof Isberg, Melissa Farrow, Jeff Spraggins, Martin Dufrense 2023. High Spatial Resolution MALDI Imaging Mass Spectrometry Data Acquisition. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.n92ldpq99l5b/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 08, 2023

Last Modified: October 18, 2023

Protocol Integer ID: 76683

Keywords: MALDI IMS, Lipid Imaging, maldi imaging mass spectrometry experiment, imaging mass spectrometry, mass spectrometry, negative ionization mode data acquisition in qtof mode, negative ionization mode data acquisition, spatial resolution imaging, bruker timstof flex instrument, imaging, instrument parameter, qtof mode

Abstract

This protocol provides the workflow and instrument parameters for setting up a MALDI imaging mass spectrometry experiment on a Bruker timsTOF Flex instrument. This protocol is intended for 10µm spatial resolution imaging of lipids (m/z 400 - 2000) in positive and negative ionization mode data acquisition in qTOF mode of operation.

Materials

Agilent ESI-L Tune Mix (Agilent Technologies, Santa Clara, CA)



- 1 Place slides in a MTP 2-slide holder. Scan slides using flatbed scanner, allowing for sufficient contrast to visualize the tissue boundary. Ensure the slides have fiducials.

Note

Fiducials are markings on the slide that can be used as reference points for teaching the slides (Step 4).

- 2 Load the 2-slide holder in the instrument. Open timsControl software and select appropriate MALDI IMS data acquisition method. Generate target profile by ensuring "Use Target Profile" is enabled. Check the height at a few different locations on the slide and ensure the difference is $<10\ \mu\text{m}$.

Note**Example Positive/Negative Ionization Mode Parameters for $10\ \mu\text{m}$ MALDI IMS:**

m/z 400 - 2000

Laser Power: 35 - 45%, Frequency: 10 000 Hz

Shots: 100 - 150 per pixel

Spot size: $10\ \mu\text{m}$ with Beam scanning function disabled

Source temperature: $\sim 50\ ^\circ\text{C}$

- 3 Calibrate the TOF with an external calibration using red phosphorus. Proceed if all calibration points are $<1\text{ppm}$ standard deviation error and the calibration score is $>99\%$.

**Note**

Red phosphorus is commonly used to perform calibration, however Agilent ESI-L Tune Mix can also be used for electrospray ionization.

List of Red Phosphorus Mass Calibration Peaks

	A	B
	Positive mode	Negative mode
	278.7633	278.7644
	402.6583	402.6594
	464.6058	464.6069
	526.5534	526.5545
	650.4484	650.4495
	774.3435	774.3445
	898.2385	898.2396
	960.1860	960.1870
	1022.1335	1022.1346
	1146.0286	1146.0297
	1393.8187	1393.8198
	1517.7137	1517.7148
	1641.6088	1641.6099
	1765.5038	1765.5049
	1889.3989	1889.4000

- 4 Use the FlexImaging software to load the optical image of the slides. Teach the target position with three teaching points. Check the accuracy of the teaching by moving the stage to a fourth location.
- 5 Select desired tissue measurement region. Start data acquisition in the FlexImaging software.