


Sep 08, 2020

Version 1

High-throughput SARS-CoV-2 RNA extraction from nasopharyngeal swabs using repurposed 3D printers. V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.bkynkxve>

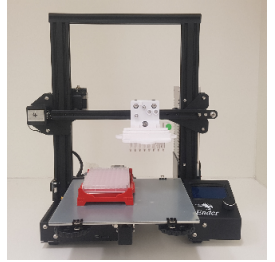

Koen Vandelannoote¹

¹Institut Pasteur du Cambodge

Coronavirus Method De...

XPRIIZE Rapid Covid Tes...

1 more workspace



Koen Vandelannoote

Institut Pasteur du Cambodge

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bkynkxve>

Protocol Citation: Koen Vandelannoote 2020. High-throughput SARS-CoV-2 RNA extraction from nasopharyngeal swabs using repurposed 3D printers.. protocols.io <https://dx.doi.org/10.17504/protocols.io.bkynkxve>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 07, 2020

Last Modified: September 08, 2020

Protocol Integer ID: 41710

Keywords: Ender VX-500, RNA extraction, 3D printer, SARS-CoV-2, nasopharyngeal swabs, COVID-19, rna extraction, extraction of sar, human upper respiratory sample, nasopharyngeal swab, rna, paramagnetic bead, throughput sar, respiratory, extraction

Abstract

Here, we describe a semi-automated workflow for paramagnetic bead-based extraction of SARS-CoV-2 RNA from human upper respiratory samples that uses low-cost repurposed *Creality3D* Ender-3 3D-printers and readily accessible reagents. The protocol described here is for the use of four instruments in parallel, each processing 24 samples, 96 samples per batch.

Attachments



20200908_Ender_VX-

50...

41KB



Icon

for

RNA ex v

RNA_ex_v10.gcode

47KB

Materials

REAGENTS:

- Lysis buffer ([M] 6 Mass Percent Guanidinium thiocyanate , [M] 150 millimolar (mM) Tris-Cl , [M] 30 millimolar (mM) EDTA , [M] 90 millimolar (mM) DTT , 3% (v/v) Triton X-100).

- SPRI bead mix prepared for RNA binding as described in https://openwetware.org/wiki/SPRI_bead_mix

Note

This SPRI bead mix is a drop-in substitute for RNAClean XP beads (Beckman Coulter) at about 1/100 of the cost.

- 100% Ethanol
- 80% Ethanol solution
- Nuclease free water (NFW)

PLASTIC CONSUMABLES:

- 96-well plates with V-shaped wells (Greiner – Ref. 651201)

Note

It is highly recommended to work with the specified 96-well plates. The use of alternative plates requires careful modifications to the Ender VX-500 "print protocol".

- 8-tube PCR strips (SSlbio - Ref. 3247-00)

Note

It is highly recommended to work with the specified strip tubes. The use of alternative strips requires careful modifications to the Ender VX-500 "print protocol".

- 96 deep-well plates (Axygen product number P-DW-11-C-S)



Safety warnings


- ⚠ Nasopharyngeal swab samples that may contain respiratory viruses should be handled in a Class II BSC. Standard PC2 laboratory personal protective equipment is required. This includes a long sleeved laboratory gown, gloves and safety glasses. Additionally, a face mask should be worn.

The lysis buffer contains guanidine thiocyanate which may produce hazardous gases (including hydrogen cyanide and chlorine) when combined with bleach (sodium hypochlorite) and/or strong acids.

Ethanol is a highly flammable liquid.

Before start


This protocol provides directions for the use of a 3D Printer RNA extraction apparatus we named the "Ender VX-500". The platform essentially consists of a *Crealty3D Ender-3* printer repurposed by installing three different custom designed interlocking attachments on the printer's extruder mounting mechanism.

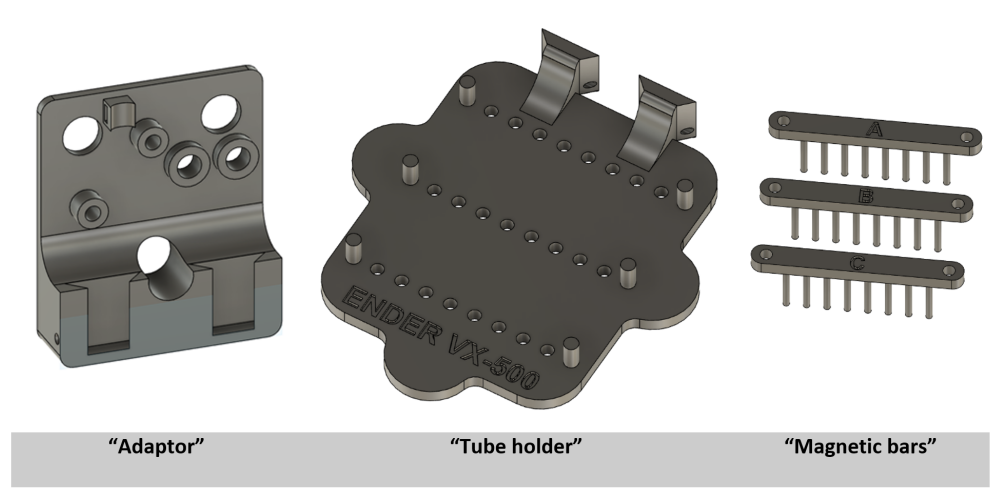

These models can be obtained from https://www.thingiverse.com/koen_vdl/designs

A 3D print "job" can be executed on the printers which will emulate the movements required to make the magnetic combs on the mounted attachments run a paramagnetic bead-based RNA extraction protocol.

Here we describe how we can efficiently process 91 nasopharyngeal swab samples in parallel on 4 Ender VX-500 machines. The following is a graphical overview of the extraction procedure.

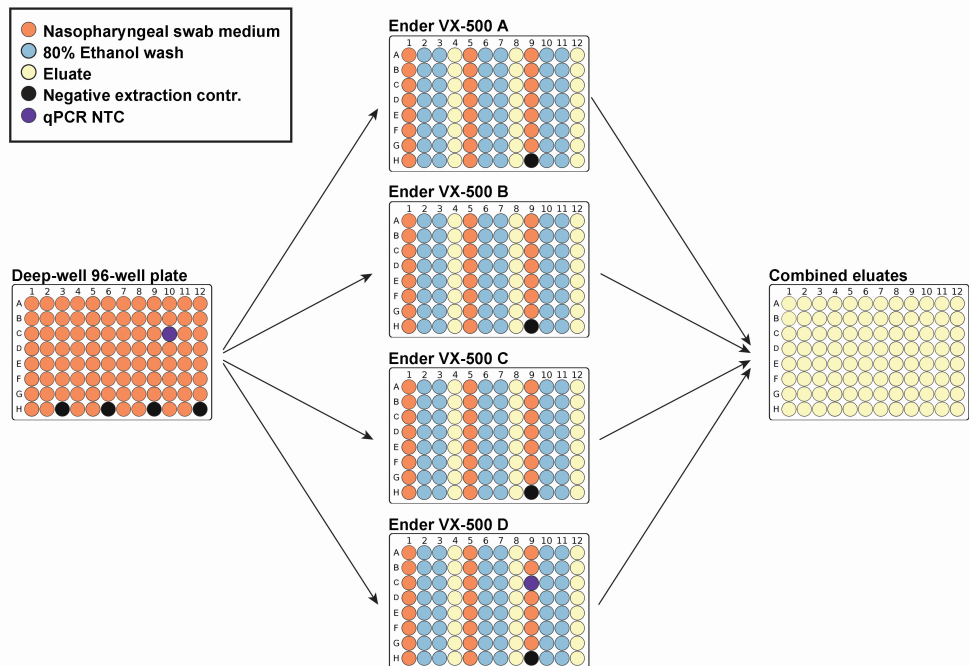

Pre-process sample information

1

Scan the barcodes of 91 nasopharyngeal swabs using a barcode scanner and the "20200908_Ender_VX-500_extraction_template.xlsx" spreadsheet.

- Populate the "SampleID" column of the "Sample log" sheet
- Randomly include a negative extraction control for every set of 24 swabs in the sheet.
- Randomly include a PCR no-template-control (NTC) control in the sheet.

Note

The spreadsheet will automatically parse 5 plate dispensing charts that can be printed for convenience:

- 4 charts represent the plate layouts of the 4 Ender VX-500 extraction runs (of 24 samples each).
- 1 chart represents the plate layout of the 96-well plate on which the eluates of the 4 extraction runs will be combined.

Note

The "20200908_Ender_VX-500_extraction_template.xlsx" spreadsheet can be found as an attachment under the abstract tab.

Pre-dispense the samples into a deep-well 96 well plate

- 2 Transfer 100 uL of swab transport medium into a deep-well 96 well plate in a Class II BSC using the plate layout charts as a guide.

Prepare four Ender VX-500 machines for extraction

- 3 Attach three 8-tube PCR strips onto the "Tube holder".

Note

- The caps of the strip fit tightly into the openings on the "Tube holder".
- Only hold the strips by the caps and make sure not to touch the outside of the tubes.

- 4 Seat the "Tube holder" into the "Adaptor".



- 5 Lock the "Tube holder" in place by inserting the provided Allan hex wrench. This will ensure the "Tube holder" is always inserted in the same position.
- 6 Place three "Magnetic bars" in the openings of the "Tube holder".

Prepare four 96-well plates for extraction

- 7 Add  18.5 μL Lysis buffer to columns 1, 5 and 9 of a 96-well plates with V-shaped wells using a multichannel pipette.

Note

Resuspend any salted out guanidine thiocyanate crystals before dispensing the buffer.

Incubate  65 °C for  00:10:00

- 8 Add  80 μL NFW to columns 4, 8 and 12 using a multichannel pipette.

- 9 Add  150 μL  80 % (v/v) Ethanol to columns 2, 3, 6, 7, 10 and 11 using a multichannel pipette.

Note

- The ethanol concentration will quickly lower in the air flow of the BSC.
- Always use a freshly made up Ethanol solution.

- 10 Transfer / pipet-mix  37 μL of nasopharyngeal swab transport medium from the deep-well plate into columns 1, 5 and 9 using using a multichannel pipette and the relevant plate dispensing chart
- 11 Incubate the plate  Room temperature for  00:01:00 .



- 12 Transfer / pipet-mix  30 μL SPRI bead mix into columns 1, 5 and 9 using a multichannel pipette.

Note

Shake the SPRI bead mix before use (the beads settle quickly).

- 13 Transfer / pipet-mix  85.5 μL [M] 100 % (v/v) Ethanol to columns 1, 5 and 9 using a multichannel pipette.

- 14 Incubate the plate  Room temperature for  00:05:00 .

Note

The prepared plates need to be extracted on the Ender VX-500 machines immediately after this incubation step

Extract 4 plates on 4 Ender VX-500 machines

- 15 Insert the four 96-well plates onto four Ender VX-500 machines prepared for extraction

Note

Ensure the microplates are correctly oriented by checking the position of the A1 marker.

- 16 Power on the Ender VX-500 machines.

- 17 Use the knob on the Ender VX-500 to start the extraction by selecting "Print from TF" and selecting "RNA_ex_v10.gcode".

Note

The "RNA_ex_v10.gcode" print job which emulates the RNA extraction can be found as an attachment under the abstract tab.

18 The Ender VX-500 will:

- Collect bead-RNA complex from the lysate (⌚ 00:05:00)
- Wash 1: vortex (⌚ 00:01:00) & collect bead-RNA complex (⌚ 00:01:30)
- Wash 2: vortex (⌚ 00:01:00) & collect bead-RNA complex (⌚ 00:01:30)
- Air dry (⌚ 00:05:00)
- Elute: vortex (⌚ 00:01:00) & collect beads (⌚ 00:01:00)
- Raise the magnetic bars out of the eluates

Note

After ⌚ 00:20:00 the machine will beep to indicate the run completed successfully.

19 Remove the 96-well plates from the Ender VX-500 machines.

20 Power off the Ender VX-500 machines.

21 Place the 96-well plates on a 96 well magnetic rack and use a multichannel pipette to transfer  50 μ L of eluate onto a fresh 96-well plate that will combine the eluates of the 4 extraction runs.

Note

 30 μ L of eluate is sacrificed to prevent any bead SPRI bead carryover.

Machine clean-up

22 Carefully remove the "Tube holder" from the EnderVX-500 "Adaptor".



- 23 Use paper tissues to release the three 8-tube PCR strips from the "Tube holder" and decontaminate the attachment in a bath containing a 1% sodium hypochlorite solution for min  00:10:00 .