

Jan 23, 2024

Histone extraction HLB protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.x54v9p49pg3e/v1>

Sigrid Verhelst¹

¹Laboratory of Pharmaceutical Biotechnology, Ghent University, Ghent, Belgium



Sigrid Verhelst

ProGenTomics, Laboratory of Pharmaceutical Biotechnology, Gh...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x54v9p49pg3e/v1>

Protocol Citation: Sigrid Verhelst 2024. Histone extraction HLB protocol. protocols.io

<https://dx.doi.org/10.17504/protocols.io.x54v9p49pg3e/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 22, 2024

Last Modified: January 23, 2024

Protocol Integer ID: 93870

Keywords: acid extraction of histone, histone protein, hypotonic lysis buffer for isolation, hypotonic lysis buffer, histone, extraction, acid extraction, hlb protocol protocol, lysis buffer

Funders Acknowledgements:

BOF

Grant ID: Mandate 01D23313

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Protocol to extract histone proteins using a hypotonic lysis buffer for isolation of nuclei prior to acid extraction of histones that is easily amenable to label-free MS owing to the lack of detergents in the lysis buffer.

Isolate nuclei

- 1 Start from a dry (snap-frozen) cell pellet
- 2 Add hypotonic lysis buffer (HLB) to cell pellet: **200 μ L for 1×10^6 cells** and resuspend softly

Note

HLB buffer: 10 mM Tris-HCl pH 8.0, 1 mM KCl, 1.5 mM MgCl₂ supplemented with 1 mM DTT, 1 mM PMSF, Halt Protease and Phosphatase Inhibitor Cocktail 100x (78440) and phosphatase inhibitor cocktails II and III (P5726 and P0044, Sigma-Aldrich, 1 mL of cocktail for 100 mL of buffer)

- 3 Rotate for 30 min at 4°C to promote lysis of cell membrane (mechanical shear)
- 4 Pellet the nuclei by centrifugation at 10.000g for 10 min at 4°C
- 5 Discard the supernatant

Extract histones

- 6 Resuspend the pellet in 0.4N hydrochloric acid (HCl) by soft pipetting until no clumps left in solution by adding **125 μ L HCl for 1×10^6 cells** (if necessary: vortex)
- 7 Incubate for **30 min** in acid on rotator at 4°C to promote lysis of nuclei and solubilization of histones
- 8 Spin down for 10 min at 4°C and 16.000 g
- 9 Transfer supernatant to new Eppendorf (histones are present in the acid since they are alkaline proteins)



Isolate histones

10 Add, drop by drop, trichloroacetic acid (TCA) until a final concentration of 33% is reached to promote precipitation of histones and invert the tube several times (results in a milky solution)

11 Incubate on ice for 30min

12 Spin for 10 min at 4°C and 16.000 g to pellet the histones

13 Remove the supernatant

Note

Be careful: the pellet is not always visible



Wash steps

14 Add ice-cold acetone (do not resuspend the pellet) to remove TCA, make sure the pellet is fully covered with acetone

15 Spin for 5 min at 4°C and 16000 g

16 Remove the supernatant

17 Add cold acetone again (do not resuspend the pellet) to remove TCA

18 Spin for 5 min at 4°C and 16000 g

19 Remove the supernatant

20 Dry at room temperature for 30 min (until no acetone left)

**Note**

Samples can be stored at -80°C or -20°C until further use (propionylation/digestion) or part of the sample can be prepared to perform sodium dodecyl sulfate polyacrylamide gel electrophoresis (SDS-PAGE)

Preparation for SDS-PAGE

21 Resuspend in MilliQ water (50 µl for 1×10^6 cells)

22 Transfer 400.000 cells to a new Eppendorf tube for gel-electrophoresis (optionally)

Note

If there are still clumps left: Spin for 10 min at 4°C and 16000 g and transfer the supernatant in a fresh Eppendorf

23 Vacuum dry the samples (SpeedVac)

Note

Store both Eppendorf tubes (for propionylation/digestion and SDS-PAGE) at -20°C or -80°C until further use