


Sep 25, 2019

Hornwort RNA extraction

DOI

<https://dx.doi.org/10.17504/protocols.io.p28dqhw>

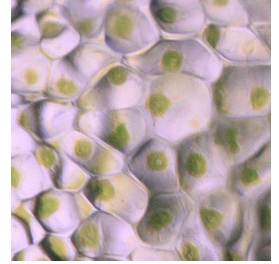

Eftychis Frangedakis¹

¹University of Cambridge



Eftychis Frangedakis

University of Cambridge, Plant Sciences

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.p28dqhw>

Protocol Citation: Eftychis Frangedakis 2019. Hornwort RNA extraction. protocols.io

<https://dx.doi.org/10.17504/protocols.io.p28dqhw>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 11, 2018

Last Modified: September 25, 2019

Protocol Integer ID: 12096

Keywords: rna extraction the same protocol, rna extraction, overnight rna precipitation step with licl, rna extraction with the addition, overnight rna precipitation step, dna extraction, rna, extraction

Abstract

The same protocol used for DNA extraction is used for RNA extraction with the addition of an overnight RNA precipitation step with LiCl.

Materials

MATERIALS

 Polyvinylpyrrolidone **Merck MilliporeSigma (Sigma-Aldrich) Catalog #PVP40**

 Chloroform **Merck MilliporeSigma (Sigma-Aldrich) Catalog #319988**

Extraction buffer:

100mM Tris-HCL pH 8 (2M)	25ml
1.4M NaCl (5M)	140ml
20mM EDTA pH 8 (0.5M)	20ml
2% CTAB	10gr
0.3% β -mercaptoethanol	150 μ l/50ml

Safety warnings

 All solutions used during this protocol were prepared using DEPC water!!



- 1 Grind 0.5-2 g of tissue using mortar and pestle in the presence of liquid nitrogen until finely ground. Transfer frozen ground tissue to a 30 ml tube.
- 2 Add 10 ml of 60 °C extraction buffer and 100 mg PVP-40/g tissue. Mix by inversion and incubate in 60 °C in a water bath for 30 min (or incubate on ice for 15 min).
- 3 Remove from heat, and let cool to room temperature for 4 to 6 min.
- 4 Add 12 ml of chloroform:IAA (24:1) and mix by inversion to form an emulsion.
- 5 After mixing thoroughly, spin at 10,000 rpm for 10 min at RT.
- 6 Transfer aqueous phase to a new 30 ml centrifuge tube. Repeat chloroform:IAA extraction to remove cloudiness (PVP) in aqueous phase. Spin at 7,000 rpm for 10 min at RT.
- 7 Add ½ volume of 5M NaCl to the final aqueous solution recovered. Mix well. Add two volumes of cold (-20 °C) 95% ethanol. Mix by inversion. Place in freezer (-20 °C) for 10 min to accentuate precipitation. Spin at 13,000 rpm for 6 min.
- 8 Optional: Resuspend in 2 ml of TE and repeat step 7.
- 9 Pour off supernatant. And wash pellet with cold (0 to 4 °C) 70% v/v ethanol. Dry pellet.
- 10 Dissolve in 400 µl TE and add 100ul 8M LiCl. Incubate o/n at 4 °C.
- 11 Spin 10 min to pellet RNA . Remove supernatant (contains DNA).
- 12 Dissolve pellet with 400ul DEPC dH₂O. Add 40ul NaoAc (pH 4.5) and 1ml 95% EtOH and freeze at -80 oC for 10 min. (optional: wash pellet with 95% EtOH).
- 13 Spin for 10 min at 13K and resuspend pellet in 40ul DEPC dH₂O.



- 14 If necessary (pellet not dissolved): incubate at 60 °C for 10 min and then on ice for 30 min. Pellet debris by spinning tube for 10 min at 13K. Transfer supernatant (RNA) to a new tube. Extraction buffer: 100mM Tris-HCL pH 8 (2M) 25ml.