

May 19, 2023

Version 2

HotSHOT genomic DNA extraction V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlkw945l5r/v2>

ZEBRAFISH UCL



FishFloorUCL¹

¹University College London

FishFloorUCL



Francois Kroll

Sorbonne Université

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlkw945l5r/v2>

Protocol Citation: FishFloorUCL 2023. HotSHOT genomic DNA extraction. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.j8nlkw945l5r/v2> Version created by **Talia Pittman**



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: March 31, 2023

Last Modified: May 18, 2026

Protocol Integer ID: 79829

Keywords: DNA, extraction, HotShot, danio rerio, zebrafish, hotshot genomic dna extraction, genomic dna from larvae, dna extraction, hotshot method, genomic dna, larvae, extraction, dna

Abstract

How to extract genomic DNA from larvae or finclips using the HotSHOT method.



Materials

1 Prepare the BASE stock solution (50X):

- 14.03 g KOH crystals (1.25M final concentration)
- 4 mL of 0.5M EDTA (10 mM final concentration)
- ddH₂O to 200 mL total volume

Can be stored at room temperature for up to four years

2 Prepare the NEUTRALISATION stock solution (50X)

- 63.04 g Tris-HCL (2M final concentration), also called Trizma HCl
- ddH₂O to 200 mL total volume

Can be stored at room temperature for up to four years

Procedure

30m

3 Prepare fresh 1X BASE and 1X NEUTRALISATION solution in nuclease-free H₂O.

4 Transfer finclips or culled larvae to individual wells of a 96-well plate.

Note

(Finclips) Transfer directly to a 96-well plate during fin-clipping

(Larvae) After culling:

- Pipette technique: Use a pipette set to 5µL to transfer larvae with as little fish water as possible
- Forceps technique: Transfer individual larvae using blunt round-ended forceps. This technique is only possible if you have already loaded the plate with 50µL 1X BASE solution (step 5)

5 Add 50 µL of 1X BASE solution into each well of the plate

6 Seal the plate and place on PCR block 95 °C 00:30:00

30m

7 Cool at Room temperature



- 8 Add  50 μL of 1x NEUTRALISATION solution into each well of the plate.
- 9 Note, the extracted DNA concentration is often *too high* for downstream applications like PCR or KASP.

Storage

- 10 Store at  4 °C if you will use the DNA in the next few weeks. Beware, the samples will slowly evaporate from a sealed plate. Alternatively,  -20 °C for long-term storage. This will also prevent the samples from evaporating.