



Aug 28, 2020 Version 2

# 🌐 HTAPP\_Depletion of CD45+ cells from ovarian cancer ascites single cell suspensions for single-cell RNA-Seq V.2

DOI

[dx.doi.org/10.17504/protocols.io.bhbij2ke](https://dx.doi.org/10.17504/protocols.io.bhbij2ke)

Benjamin Izar<sup>1,2</sup>, Parin Shah<sup>1,2</sup>, Mei-Ju Su<sup>1,2</sup>, Isaac Wakiro<sup>1</sup>, Sara Napolitano<sup>3</sup>, Jingyi Wu<sup>1,2</sup>, Sébastien Vigneau<sup>1</sup>, Asaf Rotem<sup>1,2</sup>, Orit Rozenblatt-Rosen<sup>4,2</sup>, Bruce Johnson<sup>1,2</sup>, Aviv Regev<sup>4,5,6,7</sup>

<sup>1</sup>Dana-Farber Cancer Institute; <sup>2</sup>Human Tumor Atlas Pilot Project; <sup>3</sup>University of California, Santa Barbara;

<sup>4</sup>Broad Institute; <sup>5</sup>Human Tumor Atlas Project; <sup>6</sup>Massachusetts Institute of Technology;

<sup>7</sup>Howard Hughes Medical Institute

NCIHTAN



Judit Jane-Valbuena

The Broad Institute

OPEN  ACCESS



DOI: [dx.doi.org/10.17504/protocols.io.bhbij2ke](https://dx.doi.org/10.17504/protocols.io.bhbij2ke)

**Protocol Citation:** Benjamin Izar, Parin Shah, Mei-Ju Su, Isaac Wakiro, Sara Napolitano, Jingyi Wu, Sébastien Vigneau, Asaf Rotem, Orit Rozenblatt-Rosen, Bruce Johnson, Aviv Regev 2020. HTAPP\_Depletion of CD45+ cells from ovarian cancer ascites single cell suspensions for single-cell RNA-Seq. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bhbij2ke>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** June 08, 2020

**Last Modified:** August 28, 2020

**Protocol Integer ID:** 37962



## Abstract

Many ovarian cancer ascites samples have a high proportion of CD45-expressing cells (>90%) (CD45+) and a low proportion (<10%) of malignant cells. In such cases, depletion of CD45+ immune cells can be used to increase the proportion of malignant cells, as well as other non-immune cell types, without relying on the expression of epithelial or tumor-specific markers.

For the Human Tumor Atlas Pilot Project (HTAPP) this protocol was used to deplete CD45+ immune cells from single-cell suspensions obtained from ovarian cancer ascites, and is included as a step in the "HTAPP\_Processing human ovarian cancer ascites to a single-cell suspension for single-cell RNA-seq" protocol, also available in [protocols.io](#). It is adapted from the CD45 MicroBeads MACS Separation [protocol](#) from Miltenyi Biotec (Cat No 130-045-801).

Description of this protocol and guidance for testing and selecting methods for processing different tumor and sample types can be found in [Slyper et al.](#)

## Guidelines

- Keep sample on ice and use cold reagents unless noted otherwise.
- Report information as indicated in the protocol.

## Materials

### MATERIALS

- ⊗ LS Columns **Miltenyi Biotec Catalog #130-042-401**
- ⊗ PBS pH 7.4 **Thermo Fisher Scientific Catalog #10010049**
- ⊗ ULTRAPURE 0.5M EDTA pH 8.0 **Thermo Fisher Scientific Catalog #15575020**
- ⊗ ANTI-HUMAN CD45 MICROBEADS **Miltenyi Biotec Catalog #130-045-801**
- ⊗ BSA **Cell Signaling Technology Catalog #9998S**
- ⊗ MIDIMACS SEPARATOR **Miltenyi Biotec Catalog #130-042-302**
- ⊗ MACS MULTISTAND **Miltenyi Biotec Catalog #130-042-303**
- ⊗ Falcon® 15 mL High Clarity PP Centrifuge Tube Conical Bottom with Dome Seal Screw Cap Sterile 50/Rack 500/Case **Corning Catalog #352097**

## Safety warnings

- ! ▪ Follow general lab safety and institutional guidelines when working with sharps and human derived samples.



## Before start

- Set centrifuge to 4°C.
- Prepare MACS buffer (PBS with 0.5% BSA and 2 mM EDTA), sterile filter, degas, and keep on ice. This solution can be prepared in advance and stored at 4°C for several weeks.
- Label a 15 mL conical tube as "Wash".
- Label a 15 mL conical tube as "CD45 negative fraction".

## Quality Control

- 1 Mix 5  $\mu\text{L}$  of red blood cell-free ascites single-cell suspension with 5  $\mu\text{L}$  Trypan blue and load on hemocytometer. The ascites cell suspension can be obtained following the "HTAPP\_Processing human ovarian cancer ascites to a single-cell suspension for single-cell RNA-seq" protocol.
- 2 Count and report the number of viable single cells, dead single cells, cell doublets or aggregates, and whether debris are present, then calculate additional quality control metrics below. Take picture if possible.


|   | Quality Control |
|---|-----------------|
| <b>Number of Viable Single Cells Counted</b>  |                 |
| <b>Number of Dead Single Cells Counted</b>  |                 |
| <b>Number of Cell Aggregates or Doublets Counted</b>                                    |                 |
| <b>Concentration of Viable Single Cells (cells/<math>\mu\text{L}</math>)</b>            |                 |
| <b>Concentration of Dead Single Cells (cells/<math>\mu\text{L}</math>)</b>              |                 |
| <b>Concentration of Cell Aggregates or Doublets (doublets/<math>\mu\text{L}</math>)</b> |                 |
| <b>Volume of Single Cell Suspension (<math>\mu\text{L}</math>)</b>                      |                 |
| <b>Total Number of Viable Single Cells</b>  |                 |
| <b>Proportion of Single Cells that are Viable (%)</b>                                   |                 |
| <b>Proportion of Cell Aggregates or Doublets (%)</b>                                    |                 |
| <b>Description of debris (if any)</b>   |                 |

### Note











Insert Picture:

## Magnetic Labeling

- 3 Centrifuge at 580 g for 5 minutes in 4°C pre-cooled centrifuge.

 00:05:00 Centrifugation

 4 °C Wet Ice

- 4 Remove supernatant as completely as possible without disrupting the cell pellet.  
 4 °C Wet Ice
- 5 If cell number is lower than 10 million, resuspend the cell pellet in 80  $\mu$ L MACS buffer. If cell number is higher than 10 million cells, resuspend the cell pellet in 80  $\mu$ L MACS buffer for every 10 million cells.  
 4 °C Wet Ice
- 6 Add 20  $\mu$ L CD45 microbeads per 80  $\mu$ L MACS buffer.  
 4 °C Wet Ice
- 7 Incubate on ice or in 4°C refrigerator for 15 minutes.  
 00:15:00 Labeling  
 4 °C Wet Ice
- 8 During the incubation, prime an LS column: insert the column into a MidiMACS separator and let 3 mL of MACS buffer flow through the column, collecting it in the 15 mL "Wash" tube. Once the column is primed, discard the 15 mL "Wash" tube and position the 15 mL "CD45 negative fraction" tube under the column.
- 9 Following incubation (Step 7), resuspend the labeled cell suspension with 900  $\mu$ L cold MACS buffer per 100  $\mu$ L of suspension.  
 4 °C Wet Ice
- 10 Spin the suspension at 500 g for 4 minutes in 4°C pre-cooled centrifuge.  
 00:04:00 Centrifugation  
 4 °C Centrifuge
- 11 Remove supernatant as completely as possible without disrupting the labeled cell pellet.  
 4 °C Wet Ice
- 12 Resuspend the pellet (up to 100 million cells) in 500  $\mu$ L cold MACS buffer.  
 4 °C Wet Ice

## Magnetic Separation

- 13 Transfer the labeled cell suspension to the primed LS column on the MidiMACS separator and collect flow-through in the 15 mL "CD45 negative fraction" tube. This CD45 depleted cell suspension can be used in the "HTAPP\_Processing human ovarian cancer ascites to a single-cell suspension for single-cell RNA-seq" protocol.



*Tip: Cell recovery may be increased by washing the column up to three times with 3 mL of MACS buffer.*

🌡 Room temperature