

Apr 12, 2023

Version 2

HTTPM : Illumina library preparation V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.n2bvj8oowgk5/v2>

Antoine Champie¹, Amélie De Grandmaison¹

¹Université de Sherbrooke



Antoine Champie

Université de Sherbrooke

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n2bvj8oowgk5/v2>

External link: <https://doi.org/10.1371/journal.pone.0283990>

Protocol Citation: Antoine Champie, Amélie De Grandmaison 2023. HTTPM : Illumina library preparation. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.n2bvj8oowgk5/v2>

Manuscript citation:

Champie A, Grandmaison AD, Jeanneau S, Grenier F, Jacques P, Rodrigue S (2023) Enabling low-cost and robust essentiality studies with high-throughput transposon mutagenesis (HTTPM). PLoS ONE 18(4): e0283990. doi: [10.1371/journal.pone.0283990](https://doi.org/10.1371/journal.pone.0283990)

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 24, 2022

Last Modified: April 12, 2023

Protocol Integer ID: 69139

Keywords: illumina library preparation part of the http protocol, sequencing library, illumina library preparation part, preparation of illumina, sequencing, http protocol, illumina, library preparation

Abstract

Part of the HTTP protocol dedicated to the preparation of Illumina sequencing libraries.

Materials

Preparation of Nextera Adaptaters :

Nextera (NxT) adapters are prepared by hybridisation of the following primers :

	A	B
	Nxt-XTv2-B-N701-T	CAAGCAGAAGACGGCATACGAGATTCGCCTTAGTCTCGTGGGCTCGGAGATGTGTATAAGAGACAGT
	Nxt-XTv2-B-3R-ac3-5phos	/5Phos/CTGTCTCTTATACACATCTCCGAGCCCACGAGAC/3InvdT/

■ **Preparation of the 5X annealing buffer (5X Tris NaCl buffer : 50 mM Tris, pH 7.5-8, 250 mM NaCl) :**

- 500 µl Tris-HCl 1M pH 7.5
- 500 µl NaCl 5M
- 9 ml H₂O mol.-grade

■ **Preparation of the adapters (40 µM 50 µL) :**

- Resuspend both primers in water to obtain 100 µM stocks
- Mix 20 µl of each (Nxt-XTv2-B-N7XX-T and Nxt-XTv2-B-3R-ac3-phos5')
- Add 10 µl of 5X annealing buffer
- Annealing reaction in a thermocycler (decrease temperature from 98 to 4C (-0.1C/cycle(10s/cycle)))

Primers used for the first PCR :

	A	B
	Nxt_A	AATGATACGGCGACCACCGAGATCTACAC
	Nxt_B	CAAGCAGAAGACGGCATACGAGAT

Primers template for barcoding PCR :

	A	B
	Nxt_i5_barcoding	AATGATACGGCGACCACCGAGATCTACAC [8 Nu Index] TCGTCCGCGAGCGTCAGATGTGTA



	A	B
	Nxt_i7_barcode g	CAAGCAGAAGACGGCATACGAGAT [8 Nu Index] GTCTCGTGGGCTCGGAGATGTGTATAAG

Before start

All steps and master mixes need to be kept on ice as much as possible. Thermocyclers need to be cooled at 4C before inserting sample plate.



Libraries

1h 34m

1 Transfer 2.5µl of DNA from the DNA extraction plate to a new PCR plate.

2 Prepare a fragmentation master mix with :

	A	B
	NEB Ultra II FS buffer	77 µl
	NEB Ultra II FS enzyme	22 µl
	Molecular grade water	11 µl

3 Add  1 µL of the fragmentation master mix to each well.

4 Incubate in a thermocycler with the following protocol :

45m

-  00:15:00 at  37 °C
-  00:30:00 at  65 °C

5 Add  1 µL of 4µM Nextera (NxT) adaptors to each well.

6 Prepare a ligation master mix with :

	A	B
	NEB Ultra II ligation master mix	377.4 µl
	NEB Ultra II ligation enhancer	12.1 µl

7 Add  3.5 µL of ligation master mix to each well.



8 Incubate in a thermocycler with the following protocol :

40m

- 00:30:00 at 20 °C
- 00:10:00 at 65 °C

9 Prepare a PCR master mix with :

A	B
NxT_A primer 20 µM	880 µl
Nxt_B primer 20 µM	880 µl
Molecular grade water	8360 µl
PCR Mix 2X	11000 µl

10 Add 192 µL of PCR master mix to each well.

11 Split the PCR reaction into 4 different plates (50µl per plate).

12 Incubate each plate in a thermocycler with the following cycles :

3m 15s

- 00:00:30 at 98 °C
- 00:00:15 at 98 °C
- 00:00:30 at 72 °C
- Repeat from step 2 for 20~25 cycles*
- 00:02:00 72 °C

13 Pool the 4 PCR replicates together in a.

14 Transfer 2 µL of DNA from the pool plate to a new PCR plate.



15 Add  2 μL of each barcoding primer to the DNA :

- Nxt_i5_barcoding
- Nxt_i7_barcoding

16 Prepare a PCR master mix with :

	A	B
	Molecular grade water	2090 μl
	PCR mix 2X	2750 μl

17 Add  44 μL of the PCR master mix to each well of the plate.

18 Incubate in a thermocycler with the following protocol :

3m 45s

-  00:00:30 at  98 °C
-  00:00:15 at  98 °C
-  00:01:00 at  72 °C (no anneal step)
- Repeat from step 2 for 5 cycles
-  00:02:00 at  72 °C

19 Pool together  2 μL of each sample.

20 Purify with SPRI beads using a 0.8 ratio. Resuspend with  50 μL of molecular grade water.



21 Proceed with QC and sequencing.