

Mar 18, 2025

Version 2

HuBMAP Lung TMC UCSD & URM C Protocol Overall for Multimodal Single Cell/Nucleus Assays V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l27qpqg1y/v2>

Gloria S Pryhuber¹, Elizabeth Duong², Blue Lake², Cory Poole¹, Heidie Huyck¹, Ravi Misra¹, Xin Sun², Jim Hagood³, Dinh H Diep², Kun Zhang², Zoey Zhao²

¹University of Rochester Medical Center; ²University of California San Diego;

³University of North Carolina at Chapel Hill

Human BioMolecular Atl...

URMC Pryhuber Lab



Gloria S Pryhuber

University of Rochester Medical Center

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l27qpqg1y/v2>

Protocol Citation: Gloria S Pryhuber, Elizabeth Duong, Blue Lake, Cory Poole, Heidie Huyck, Ravi Misra, Xin Sun, Jim Hagood, Dinh H Diep, Kun Zhang, Zoey Zhao 2025. HuBMAP Lung TMC UCSD & URM C Protocol Overall for Multimodal Single Cell/Nucleus Assays . **protocols.io** <https://dx.doi.org/10.17504/protocols.io.4r3l27qpqg1y/v2> Version created by **Gloria S Pryhuber**

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 18, 2025

Last Modified: March 18, 2025

Protocol Integer ID: 124593

Keywords: hubmap lung tmc ucscd, lung cell type mapping, lung tissue mapping center, urmc protocol overall for multimodal single cell, use for the single nuclear rna, pipeline for the human biomolecular atlas program, single nuclear rna, human biomolecular atlas program, nucleus assay, final data to the hubmap hive, tmc urmc, tissue sample, lung tissue, datasets for hubmap, hubmap hive, frozen lung tissue, tmc ucscd, urmc protocol overall, hubmap, multimodal single cell, rna, human kidney, nucleus, isolation of nuclei, processing tissue

Funders Acknowledgements:

HuBMAP

Grant ID: U54 HL145608; U54HL165443

Abstract

This protocol is an overview of the protocols in use for the single nuclear RNA sequencing and ATACseq datasets for HubMAP-Lung Tissue Mapping Center, TMC UCSD / TMC URM C. It includes protocols.io links to each of the individual protocols that make up this project workflow, from processing tissue, sectioning, H&E imaging the tissue samples, isolation of nuclei from frozen lung tissue, running of multimodal single cell/nucleus assays.uploading the final data to the HuBMAP HIVE.

See also: Human Kidney, Urinary Tract, and Lung Cell Type Mapping Pipeline for the Human Biomolecular Atlas Program (HuBMAP) V.2 dx.doi.org/10.17504/protocols.io.bvvin64e

Materials

Required materials are provided within each linked protocol.



- 1 Screen and accept donor organs, gross and preserve the tissues into samples for analysis using the protocols:

602.2 Donor Acceptance Criteria for URMHC HTC HuBMAP and LungMAP Inclusion
dx.doi.org/10.17504/protocols.io.bjuxkxn
603.3 & 604.5_URMC_HTC_Whole Lung and Lobe Processing
dx.doi.org/10.17504/protocols.io.biz7kf9n
611.2 URMHC HTC Formalin-Inflated, Paraffin-Embedded Human Lung Tissue
dx.doi.org/10.17504/protocols.io.bjtknnn
615.1 URMHC HTC Non-Inflated Fresh-Frozen Embedded Lung and Associated Tissue
dx.doi.org/10.17504/protocols.io.bjtnknme
615.2 URMHC HTC Non-Inflated Fresh-Frozen Embedded Lung and Associated Tissue V2
dx.doi.org/10.17504/protocols.io.rm7vz84k2vx1/v2
- 2 Create the cryosections to be used in the single nuclear isolation preparation. Generally placing 10 × 40 micron sections in a single cryovial.
624.1_URMC_HTC_Cryosectioning Non-Inflated Fresh-Frozen Embedded Lung for Nuclear Isolation or Multi-plexed In Situ Hybridization
dx.doi.org/10.17504/protocols.io.3byl4j91jlo5/v1
- 3 Register via the HuBMAP UUID generator at <https://ingest.hubmapconsortium.org/>, the Donor, the Organ Lobe, the Tissue Block, and the set of tissue sections (40 micron curls) from that block used in single nuclear isolation. common coordinate region information. Record and keep locally the key for the generated UUIDs.
- 4 Register the Tissue Blocks, just registered in the UUID generator, in the Common Coordinate Framework via the Registration User Interface at <https://hubmapconsortium.github.io/ccf-ui/rui/>
- 5 Approximately every 10th section, make a 10 micron section just before and just after, the set of 40 micron curls. Stain these thinner sections by H&E.
621.1.HTC_H&E Stain (Paraffin or Cryosections)
dx.doi.org/10.17504/protocols.io.36wggjng3vk5/v1
- 6 Isolate single nuclei from the tissue curls using protocol:
Frozen lung tissue dissociation for single-nucleus RNA & Chromatin assays
dx.doi.org/10.17504/protocols.io.bh26j8he
- 7 Running of multimodal single cell/nucleus assays.
10X snRNA-seq: dx.doi.org/10.17504/protocols.io.86khzcw
10X snATAC-seq: dx.doi.org/10.17504/protocols.io.bvssn6ee
SNARE-Seq2: dx.doi.org/10.17504/protocols.io.bzdsp26e



- 8 To Be Added:
10X Chromium Single Cell *Multiome* ATAC + Gene Expression
MERFISH: dx.doi.org/10.17504/protocols.io.rm7vzjy1rlx1/v1
DARTFISH