

Dec 12, 2019

Hydra Medium 4.0

DOI

<https://dx.doi.org/10.17504/protocols.io.z4jf8un>

Rob Steele¹

¹University of California, Irvine



Rob Steele

University of California, Irvine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.z4jf8un>

Protocol Citation: Rob Steele 2019. Hydra Medium 4.0. protocols.io <https://dx.doi.org/10.17504/protocols.io.z4jf8un>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 16, 2019

Last Modified: December 12, 2019



Protocol Integer ID: 22379

Keywords: Hydra, medium, culture, hydra medium, hydra strain, recipe for the medium, detailed description of the medium, steele lab, recipe

Abstract

This is the recipe for the medium that the Steele Lab currently uses for culturing all of its Hydra strains. A detailed description of the medium is in the following document:



Hydra_Medium.pdf

Guidelines

Because we are culturing quite a few different strains of Hydra, some in large quantities, we make Hydra Medium up in 100 liter batches. This also avoids the need to make up stock solutions of the reagents. We simply add the solid reagents to 100 liters of water. There is no need to adjust the pH of the medium; the sodium bicarbonate provides buffering to the appropriate pH.

Safety warnings



The components of Hydra Medium are not dangerous.

Before start

Be sure you have a source of high purity water, such as MilliQ or Nanopure.



Fill Tank

- 1 Fill Nalgene tank with 100 liters of high purity water (Nanopure, MilliQ, or equivalent). The volume markings on the side of the tank are accurate enough for determining how much water you have added.

Add ingredients

- 2 Add anhydrous magnesium sulfate  3.97 g
Mix well with plastic paddle
- 3 Add potassium chloride  0.22 g
Mix well with plastic paddle
- 4 Add sodium bicarbonate  4.2 g
Mix well with plastic paddle
- 5 Add calcium chloride dihydrate  14.7 g
Mix well with plastic paddle