

Feb 05, 2026

Version 1

Identification of inhibitory midbrain neurons in FFPE human midbrain sections V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.dm6gp153pgzp/v1>

Anastasia Filimontseva^{1,2}, YuHong Fu^{1,2}, Glenda Halliday^{1,2}

¹Brain and Mind Centre & Faculty of Medicine and Health School of Medical Sciences, The University of Sydney, Sydney, NSW 2050, Australia;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 20815



Anastasia Filimontseva

Brain and Mind Centre & Faculty of Medicine and Health Schoo...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.dm6gp153pgzp/v1>

Collection Citation: Anastasia Filimontseva, YuHong Fu, Glenda Halliday 2026. Identification of inhibitory midbrain neurons in FFPE human midbrain sections. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.dm6gp153pgzp/v1>

License: This is an open access collection distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this collection and it's working

Created: February 05, 2026

Last Modified: February 06, 2026

Collection Integer ID: 242661

Keywords: ASAPCRN, identification of inhibitory midbrain neuron, inhibitory midbrain neuron, ffpe human midbrain sections protocols for standard processing, ffpe human brainstem section, human midbrain sections protocol, immunohistochemical staining for tyrosine hydroxylase, human brainstem section, tyrosine hydroxylase, immunohistochemical staining, ffpe

Funders Acknowledgements:

National Health and Medical Research Council Investigator Grant

Grant ID: 1176607

Aligning Science Across Parkinson's

Grant ID: ASAP-020505

Abstract

Protocols for standard processing of FFPE human brainstem sections and immunohistochemical staining for tyrosine hydroxylase (TH) and GAD67.



Files

 SEARCH

Protocol

NAME

Human Brain Tissue Processing

VERSION 1

CREATED BY



Anastasia Filimontseva

Brain and Mind Centre & Faculty of Medicine and Health School of Medical Sciences, The University of Sydney, Sydney, NSW 2050, Australia

OPEN →

Protocol

NAME

Immunohistochemical staining for GAD67 and TH in paraffin sections.

VERSION 1

CREATED BY



Anastasia Filimontseva

Brain and Mind Centre & Faculty of Medicine and Health School of Medical Sciences, The University of Sydney, Sydney, NSW 2050, Australia

OPEN →