

Jun 11, 2019

IHC on fixed frozen cryo sections

DOI

dx.doi.org/10.17504/protocols.io.32egqbe

Grace Burgin¹, Inbal Benhar¹

¹Broad Institute



Grace Burgin

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.32egqbe

Protocol Citation: Grace Burgin, Inbal Benhar 2019. IHC on fixed frozen cryo sections. [protocols.io](https://dx.doi.org/10.17504/protocols.io.32egqbe)
<https://dx.doi.org/10.17504/protocols.io.32egqbe>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 11, 2019

Last Modified: June 11, 2019

Protocol Integer ID: 24358

Guidelines

All incubation steps in humidifier tray (slide box with PBS soaked paper towels)

All wash steps in coplin jars (no agitation)



- 1 30-60 min Protein Block
5% Normal Horse (or Donkey) Serum in PBS + 0.3% Triton-X
- 2 Overnight (~18-22hrs) Primaries at 4C in PBS + 0.3% Triton-X
- 3 Wash 2x5 min w/ PBS
- 4 2hr Secondaries at room temp in PBS + 0.3% Triton-X
- 5 Wash 2x5 min w/ PBS
- 6 Coverslip