


Apr 14, 2022

Imaging of Calcium Dynamics in Vasoactive Intestinal Peptide-expressing Neurons of Enteric Nervous System

DOI

dx.doi.org/10.17504/protocols.io.14egn76xpv5d/v1

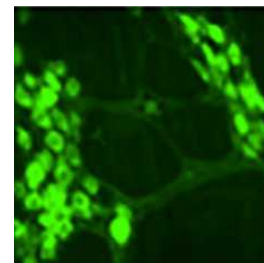

Joseph Margiotta¹, Marthe Howard¹

¹Department of Neurosciences, University of Toledo College of Medicine and Life Sciences, Toledo, Ohio

SPARC

Tech. support email: info@neuinfo.org



Marlena S Pela

University of California, San Diego

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.14egn76xpv5d/v1>

Protocol Citation: Joseph Margiotta, Marthe Howard 2022. Imaging of Calcium Dynamics in Vasoactive Intestinal Peptide-expressing Neurons of Enteric Nervous System. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.14egn76xpv5d/v1>

**Manuscript citation:**

<https://doi.org/10.1016/j.jcmgh.2021.08.016>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 11, 2022

Last Modified: April 14, 2022

Protocol Integer ID: 59375

Keywords: vasoactive intestinal peptide, neurons of enteric nervous system, enteric nervous system, imaging of calcium dynamic, gcamp reporter mouse, calcium dynamic, laboratory animal care, approved animal care, association of laboratory animal care, toledo health science campus animal care, animal care, expressing neuron, animal

Abstract

The protocol was published on behalf of the investigators by the SPARC project team.

The University of Toledo Health Science Campus animal care and use committee approved animal care, breeding procedures, and experimental protocols. Animals were housed in an Association of Laboratory animal Care-approved facility with a 12-hour light cycle with food (standard chow) and water ad libitum.

One male VIP-GCaMP reporter mouse, age 4 months, was used in this study. The mouse was purchased from the Jackson Laboratories, and all genotyping was performed using primers recommended by the Jackson Laboratories according to their protocols.

Materials

Mouse strains:

A	B	C	D	E	F
Formal strain name	Commonly used name	Mouse genome informatics ID	Jax Lab stock number	RRID	Strain background
Vip-tm1(cre)Zjh-J	Vip-IRES-cre	4431361	031628	IMSR_JAX:031628	C57BL/6J
B6J.Cg-Gt(ROSA)26Sortm95.1(CAGGCaMP6f)Hze/MwarJ	Ai95(RCL-GCaMP6f)-D(C57BL/6J) or Ai95D(C57BL/6J)	5558090	028865	IMSR_JAX:028865	C57BL/6J

Mice that contained a floxed GCaMP6f construct within the Gt(ROSA)26 locus were bred with mice that expressed Cre recombinase under the control of the Vip promoter.

Artificial cerebrospinal fluid:

NaCl - catalogue #: [S9888](#), Sigma, St. Louis, MO
 KCL - catalogue #: [P3911](#), Sigma, St. Louis, MO
 NaHCO₃ - catalogue #: [S6014](#), Sigma, St. Louis, MO
 NaH₂PO₄ - catalogue #: [S8282](#), Sigma, St. Louis, MO
 MgSO₄•7H₂O - catalogue #: [230391](#), Sigma, St. Louis, MO
 CaCl₂ - catalogue #: [C1016](#), Sigma, St. Louis, MO
 D-glucose - catalogue #: [G5767](#), Sigma, St. Louis, MO
 sodium butyrate - catalogue #: [B5887](#), Sigma, St. Louis, MO
 sodium acetate - catalogue #: [S2889](#), Sigma, St. Louis, MO

Equipment:

Picospritzer II, [Parker Instrumentation Corp](#), Barnstable, UK
[DM6000 FS Leica](#) fluorescence microscope - Leica, Buffalo Grove, IN
[Prime 95B](#) CMOS camera - Teledyne Photometrics, Tuscon, AZ
[Metamorph](#) software - Molecular Devices, Downingtown, PA

Troubleshooting



1 On the day of the experiment prepare solutions of:

1. *Artificial cerebrospinal fluid (CSF)* containing:

- 117.9 mmol/L NaCl
- 4.7 mmol/L KCl
- 25 mmol/L NaHCO₃
- 1.3 mmol/L NaH₂PO₄
- 1.2 mmol/L MgSO₄•7H₂O
- 2.5 mmol/L CaCl₂
- 11.1 mmol/L D-glucose
- 2 mmol/L sodium butyrate
- 20 mmol/L sodium acetate

2. *Artificial CSF with test reagents*, e.g., agonists.

2 Euthanize the mouse.

3 Remove the colon.

4 Open the excised proximal colon segment along the mesenteric border.

5 Pinn the colon segment mucosal side down onto a Sylgard surface lining a glass coverslip attached to the bottom of a plastic imaging chamber containing artificial CSF perfused and bubbled with Carbogen (95% O₂/5% CO₂) to oxygenate and achieve

 7.4 .

Note

In experiments that involve recording neural activity leading up to spontaneous Colonic Motor Complexes (CMCs) heat the artificial CSF to  35 °C .

6 Detect GCaMP-positive neurons by their basal Ca²⁺ fluorescence.

**Note**

To do this we used an upright DM6000 FS Leica fluorescence microscope (Leica, Buffalo Grove, IL) and Prime 95B Scientific CMOS camera (Photometrics, Tucson, AZ).

Images were collected with Metamorph software (Molecular Devices, Downingtown, PA) at a 40-Hz sampling rate and 25-ms exposure time.

- 7 Record spontaneous and evoked changes in GCaMP6f-mediated Ca^{2+} fluorescence intensity.

Note

We acquired 12-bit images using a 1.44-megapixel CMOS camera capable of capturing at 40 frames/sec (Prime 95B; Teledyne Photometrics, Tuscon, AZ) controlled by Meta-Morph software (version 7.10.1.161; Molecular Devices, Silicon Valley, CA).

- 8 At second 20 focally deliver an agonist by pressure microperfusion (10psi, 10 sec; via Picospritzer II; Parker Instrumentation Corp, Barnstable, UK) from blunt glass micropipettes (diameter, 5–10 mm) delivered within 50 μm of an adjacent myenteric ganglion.

Note

Here we used dimethylphenyl-piperazinum (DMPP) 1M 10 micromolar (μM) .