



Mar 15, 2024

## Immunocytochemistry

 [Science Advances](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvor7pdv4o/v1>

anita.adami<sup>1</sup>

<sup>1</sup>Laboratory of Molecular Neurogenetics, Department of Experimental Medical Science, Wallenberg Neuroscience Center and Lund Stem Cell Center, BMC A11, Lund University, 221 84 Lund, Sweden.



Anita Adami

Lund University

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.5qpvor7pdv4o/v1>

**External link:** <https://doi.org/10.1126/sciadv.adh9543>

**Protocol Citation:** anita.adami 2024. Immunocytochemistry. [protocols.io](#)  
<https://dx.doi.org/10.17504/protocols.io.5qpvor7pdv4o/v1>

**Manuscript citation:**

Raquel Garza, Diahann A. M. Atacho, Anita Adami, Patricia Gerdes, Meghna Vinod, PingHsun Hsieh, Ofelia Karlsson, Vivien Horvath, Pia A. Johansson, Ninoslav Pandiloski, Jon Matas-Fuentes, Annelies Quaegebeur, Antonina Kouli, Yogita Sharma, Marie E. Jönsson, Emanuela Monni, Elisabet Englund, Evan E. Eichler, Molly Gale Hammell, Roger A. Barker, Zaal Kokaia, Christopher H. Douse, Johan Jakobsson (2023) LINE-1 retrotransposons drive human neuronal transcriptome complexity and functional diversification. Science Advances doi: [10.1126/sciadv.adh9543](https://doi.org/10.1126/sciadv.adh9543)

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** February 02, 2023

**Last Modified:** March 15, 2024

**Protocol Integer ID:** 76288

**Keywords:** immunocytochemistry on 2d, immunocytochemistry this protocol, immunocytochemistry, fixed cell, cell, 2d

**Funders Acknowledgements:**

**Aligning Science Across Parkinson's through the Michael J. Fox Foundation for Parkinson's Research**

Grant ID: ASAP-000520

**Swedish Research Council**

Grant ID: 2018-02694

**Swedish Brain Foundation**

Grant ID: FO2019-0098

**Cancerfonden**

Grant ID: 190326

**Barncancerfonden**

Grant ID: PR2017-0053

**NIHR Cambridge Biomedical Research Centre**

Grant ID: NIHR203312

**Swedish Society for Medical Research**

Grant ID: S19-0100

**National Institutes of Health**

Grant ID: HG002385

**Swedish Research Council**

Grant ID: 2021-03494

**Swedish Research Council**

Grant ID: 2020-01660

## Abstract

This protocol describes how to perform immunocytochemistry on 2D fixed cells



## Fixing

- 1 The cells were washed three times with DPBS (GIBCO) and fixed for 10 minutes with 4% paraformaldehyde (Merck Millipore), followed by three more rinses with DPBS.

## Blocking and permeabilisation

1h

- 2 The cells were blocked for 01:00:00 in blocking solution (KPBS 0.25% Triton X-100 (Fisher Scientific) and 5% normal donkey serum) at Room temperature .

1h

## Staining

2h

- 3 The primary antibody was added to the blocking solution and incubated Overnight at Room temperature (follow manufacturer's instructions for dilutions).
- 4 The cells were then washed the next day three times with KPBS 0.25%.
- 5 The secondary antibody was added with DAPI (1:1000; Sigma-Aldrich) to the blocking solution and incubated at Room temperature for 01:00:00 .
- 6 The cells were finally rinsed 2-3 times with KPBS and visualised on a Leica microscope (model DMI6000B).

1h

1h