

Nov 14, 2025

Immunocytochemistry Protocol for Progenitor Cells in 96-Well Plates

DOI

dx.doi.org/10.17504/protocols.io.4r3l216jgg1y/v1

Guochang Lyu¹, Ernest Arenas¹

¹Karolinska Institute

Ernest Arenas: Deceased 15.09.2024

SOX6 mDA differentiation



Roberto Garcia Swinburn

Karolinska Institute Stockholm

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l216jgg1y/v1>

Protocol Citation: Guochang Lyu, Ernest Arenas 2025. Immunocytochemistry Protocol for Progenitor Cells in 96-Well Plates. [protocols.io https://dx.doi.org/10.17504/protocols.io.4r3l216jgg1y/v1](https://dx.doi.org/10.17504/protocols.io.4r3l216jgg1y/v1)

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 09, 2025

Last Modified: November 14, 2025

Protocol Integer ID: 226843

Keywords: immunocytochemistry protocol for progenitor cell, canonical markers with fluorophore, immunocytochemistry protocol, progenitor cell, canonical marker, fluorophore, plates this protocol

Abstract

This protocol was used to tag canonical markers with fluorophores on *in vitro* differentiated KOLF2.1J cells.

Troubleshooting

Cell Seeding

- 1 Seed ~300,000 progenitor cells per well in a 96-well plate.
- 2 Culture as needed before fixation.

Fixation

37m

- 3 Remove medium from each well. 2m
- 4 Add 4% PFA to each well. Incubate for 20 minutes at 4°C. 20m

- 5 Wash wells 3× with PBS to remove residual fixative (5 minutes each wash). 15m

Permeabilization & Blocking

1h

- 6 Prepare blocking/permeabilization buffer:
 - 5% normal donkey serum
 - 0.3% Triton X-100 in PBS (PBST)
- 7 Add buffer to each well. Incubate for 1 hour at 🌡️ Room temperature 1h



Primary Antibody Incubation

16h

- 8 Dilute **primary antibodies** in PBST + 1% donkey serum. *Remember to mix antibodies with hosts that will not interact undesiredly with secondary antibodies. Example: goat primary antibody with goat secondary antibodies.*
- 9 Add to each well. Incubate **overnight at 4°C.** 🕒 Overnight 🌡️ 4 °C 16h



Secondary Antibody Incubation

45m

10 Wash wells 3× with PBST (5 minutes each wash at room temperature).

15m

11 Add **Alexa Fluor-conjugated secondary antibodies** (e.g., Alexa488, Alexa555, or Alexa647) diluted in PBST. Incubate for **30 minutes at room temperature**, protected from light.  Room temperature  00:30:00 . *Remember not to mix secondary antibodies that will interact with each other, example goat anti-donkey and donkey anti-rabbit.*

30m

Nuclear Staining

15m

12 Add **DAPI** to each well. Incubate for **15 minutes** at room temperature, **protected from light**.

15m

13 Wash **3× with PBS (5 minutes each)**.

15m

Imaging

14 Store samples in PBS (short term, <48h) or fluorescence mounting media (mid-long term).

15 Image using a fluorescence microscope (appropriate channels for Alexa dyes and DAPI).

