



Jul 03, 2024

Immunofluorescence for adherent cells



Science Advances



In 2 collections

DOI

<https://dx.doi.org/10.17504/protocols.io.ewov19z5olr2/v1>

Alexandros C Kokotos^{1,2}, Tim Ryan^{1,2}

¹Department of Biochemistry, Weill Cornell Medicine, New York, NY 10065, USA;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, Maryland 20815, USA



Alexandros C Kokotos

Weill Cornell Medicine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov19z5olr2/v1>

Protocol Citation: Alexandros C Kokotos, Tim Ryan 2024. Immunofluorescence for adherent cells. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.ewov19z5olr2/v1>

**Manuscript citation:**

Phosphoglycerate kinase is a central leverage point in Parkinson's Disease driven neuronal metabolic deficits

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 03, 2024

Last Modified: July 03, 2024

Protocol Integer ID: 102821

Keywords: ASAPCRN, immunofluorescence, immunostaining, immunocytochemistry, antibodies, immunofluorescence for adherent cell, adherent cell, immunofluorescence, immunostaining, cell

Funders Acknowledgements:

ASAP

Grant ID: 000580

Abstract

This protocol describes a method to perform immunostaining for adherent cells.

Guidelines

Cells should be cultured and handled following all local biosafety regulations.

Materials

Equipment

Inverted tweezers (like Roboz RS-5020)

Consumables

PBS (Sigma Aldrich P3813) supplemented with 0.01 % sodium azide (Sigma Aldrich S-8032)

4 % PFA in PBS solution (Thermo Scientific J19943-K2)

NH₄Cl (Sigma Aldrich A4514)

Triton X-100 (Sigma Aldrich T8787)

BSA (Sigma Aldrich A4503)

Parafilm (Electron Microscopy Sciences 70990)

Glass slides (VWR 16004-368)

Prolong Diamond mountant (Thermo Scientific P36970)

Clear nail polish



Before start

This protocol describes a method to perform immunofluorescence on adherent cells on coverslips. It can also be used for immunocytochemistry by changing the secondary antibodies and also adjusted for cells growing on plastic. This protocol has been verified for use with both cell line cells and primary cells, like neurons.



Immunofluorescence

4h

- 1 Fix cells with [IM] 4 % (v/v) PFA in PBS for ⌚ 00:10:00 . 10m
Note: Discard PFA in a special container and not in the drain.
- 2 Quench PFA by using a [IM] 50 millimolar (mM) NH_4Cl solution in PBS for ⌚ 00:10:00 . 10m
- 3 Permeabilize cells with [IM] 0.25 % (v/v) Triton X-100 in PBS for ⌚ 00:10:00 . 10m
- 4 Block with [IM] 4 % (v/v) BSA in PBS for ⌚ 01:00:00 . 1h
Note: At this step the protocol can be paused. Blocking can also be done overnight. Other blocking solutions can be used, such as goat or donkey serum matching with the species of the secondary antibodies used.
- 5 Incubate primary antibodies in [IM] 4 % (v/v) BSA in PBS for ⌚ 01:00:00 . 1h
For adherent cells on glass coverslips, the antibody solution is drop seeded on parafilm and the coverslip is placed upside down on the antibody solution using the inverted tweezers. For the 22×22 mm coverslips, 100 ul of antibody solution is sufficient for the whole coverslip to be incubated with antibody. For cells growing on plastic the antibody solution can be added in the well/dish.
- 6 Wash cells with PBS for ⌚ 00:05:00 three times. 15m
- 7 Incubate secondary antibodies in [IM] 4 % (v/v) BSA in PBS for ⌚ 01:00:00 . 1h
For immunofluorescence, we usually use Alexa dye secondary antibodies (Thermo Scientific) at a 1:500 dilution.
- 8 Wash cells with PBS for ⌚ 00:05:00 three times. 15m
- 9 Wash cells with ddH₂O to remove excess salt and protein crystals.
For cells on coverslips, the coverslips are dipped into a beaker with ddH₂O four times and excess ddH₂O is removed by touching the edge of the coverslips on absorbent paper.
- 10 Add the coverslip on a glass slide that has been previously drop seeded with Prolong Diamond mountant. Place coverslips inverted and avoid trapping any air in the process.



Allow to air dry overnight in a dark environment.

For the 22×22 mm coverslips, 50 ul of mountant is sufficient. Pipette mountant with a cut tip to avoid bubbles.

- 11 Seal the coverslips the next day using clear nail polish.