

Jul 09, 2024

Immunofluorescence of ATP13A2 and CD63

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l6xy3klqe/v1>

Stephanie Vrijzen^{1,2}, Peter Vangheluwe^{1,2}

¹Laboratory of Cellular Transport Systems, Department of Cellular and Molecular Medicine, KU Leuven, B-3000 Leuven, Belgium;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network Chevy Chase, MD 20815, USA



Stephanie Vrijzen

Laboratory of Cellular Transport Systems

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l6xy3klqe/v1>

Protocol Citation: Stephanie Vrijzen, Peter Vangheluwe 2024. Immunofluorescence of ATP13A2 and CD63. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bp2l6xy3klqe/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: December 05, 2023

Last Modified: July 09, 2024

Protocol Integer ID: 91848

Keywords: ASAPCRN, immunofluorescent visualization of atp13a2, immunofluorescence of atp13a2, atp13a2, immunofluorescent visualization, immunofluorescence, cd63 this protocol, cd63, staining of sample

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-000458

Abstract

This protocol describes the seeding, fixation and staining of samples for immunofluorescent visualization of ATP13A2 and CD63.

Materials

- 4% paraformaldehyde (Thermo Fisher Scientific): J61899
- BSA (Roth): 8076.4
- Saponin (Sigma): S-2149
- anti-ATP13A2 (Sigma): A3361
- anti-CD63 (ExBio): 11-343-C100
- alexa-488 goat anti-rabbit IgG (Invitrogen): R37116
- alexa-594 goat anti-mouse IgG (Invitrogen): A11005
- DAPI (Sigma): D9542
- FluorSave (Merck): 345789-20ml



- 1 Seed cells: 75 000 cells/well (SH-SY5Y) in a 12-well plate with coverslips. Allow the cells to attach and grow for  48:00:00 . 2d
- 2 Wash the cells with ice-cold PBS.
- 3 Fix the cells with 4% paraformaldehyde  00:30:00 at 37 °C. 30m
- 4 Wash the cells twice with ice-cold PBS.
- 5 Permeabilization and blocking with a mixture of 5% BSA (Roth) and 0.5% saponin (Sigma) (diluted in PBS) (referred to as blocking buffer) for  01:00:00 on a shaker. 1h
- 6 Incubate with primary antibody (anti-ATP13A2, A3361, Sigma; anti-CD63, 11-343-C100, ExBio) (diluted 1/200 in blocking buffer) for  02:00:00 (40 µl antibody mixture is sufficient per coverslip in a wet chamber). 2h
- 7 Wash with DPBS for  00:05:00 on a shaker, repeat this step 3x. 5m
- 8 Incubate the samples in secondary antibody (alexa-488 goat anti-rabbit IgG R37116, Invitrogen; alexa-594 goat anti-mouse IgG, A11005, Invitrogen) (diluted 1/1000 in blocking buffer) for  00:30:00 on a shaker. 30m
- 9 Wash with DPBS for  00:05:00 on a shaker, repeat this step 3x. 5m
- 10 Incubate the samples in 200 ng/ml DAPI (D9542, Sigma) for  00:15:00 . 15m
- 11 Wash with DPBS for  00:05:00 on a shaker, repeat this step 3x. 5m
- 12 Mount the samples with FluorSave (345789-20ml, Merck).



- 13 Acquire images using a LSM880 microscope (Zeiss) with a 63x objective and Airyscan detector.