

Aug 13, 2025

IMMUNOFLUORESCENCE PROTOCOL

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl46joogo5/v1>

Yuanzhuo Zhou^{1,2}, Kohei Asai^{1,2}, Hirohisa Kyogoku^{1,3}, Tomoya S. Kitajima^{1,2}

¹Laboratory for Chromosome Segregation, RIKEN Center for Biosystems Dynamics Research (BDR), Kobe, Japan;

²Graduate School of Biostudies, Kyoto University, Kyoto, Japan;

³Graduate School of Agricultural Science, Kobe University, Kobe, Japan

Designing protein-base...



Yuanzhuo Zhou

Laboratory for Chromosome Segregation, RIKEN Center for Bios...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3byl46joogo5/v1>

Protocol Citation: Yuanzhuo Zhou, Kohei Asai, Hirohisa Kyogoku, Tomoya S. Kitajima 2025. IMMUNOFLUORESCENCE PROTOCOL. protocols.io <https://dx.doi.org/10.17504/protocols.io.3byl46joogo5/v1>



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 13, 2025

Last Modified: August 13, 2025

Protocol Integer ID: 224589

Keywords: staining of mouse oocyte, immunofluorescence protocol, immunofluorescence protocol this protocol, mouse oocyte, paraformaldehyde solution

Abstract

This protocol describes the fixation and immunofluorescent staining of mouse oocyte in paraformaldehyde solution.

Materials

1.6% formaldehyde, 100 mM PIPES (pH 7.0), 1 mM MgCl₂, 0.1% Triton X-100, PBS, 3% BSA, primary antibodies, secondary antibodies, Hoechst 33342 (Invitrogen)



Protocol

- 1 Prepare fixation buffer consisting of 1.6% formaldehyde in 100 mM PIPES (pH 7.0), 1 mM MgCl₂, and 0.1% Triton X-100.
- 2 Fix oocytes in fixation buffer at room temperature for 30 minutes.
- 3 Wash oocytes 4 times in PBS containing 0.1% Triton X-100 (PBT).
- 4 Incubate oocytes in PBT at 4 °C overnight.
- 5 Prepare blocking solution PBT containing 3% BSA.
- 6 Block oocytes in blocking solution at room temperature for 1 hour.
- 7 Prepare primary antibody solution in 3% BSA-PBT.
- 8 Incubate oocytes with primary antibodies at 4 °C overnight.
- 9 Wash oocytes 4 times in 3% BSA-PBT.
- 10 Prepare secondary antibody solution in 3% BSA-PBT containing 5 µg/ml Hoechst 33342 (Invitrogen).
- 11 Incubate oocytes in secondary antibody solution at room temperature for 2 hours or at 4 °C overnight.