

Jan 09, 2026

Immunofluorescence staining of cultured mouse hippocampal neurons

DOI

<https://dx.doi.org/10.17504/protocols.io.x54v95x5pl3e/v1>

Akio Mori¹, Robert Edwards¹

¹University of California San Francisco



Akio Mori

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x54v95x5pl3e/v1>

Protocol Citation: Akio Mori, Robert Edwards 2026. Immunofluorescence staining of cultured mouse hippocampal neurons. protocols.io <https://dx.doi.org/10.17504/protocols.io.x54v95x5pl3e/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 20, 2025



Last Modified: January 09, 2026

Protocol Integer ID: 233094

Keywords: Neuron culture, Immunofluorescence, immunofluorescence staining of cultured mouse hippocampal neuron, cultured mouse hippocampal neuron, cultured neuron, immunofluorescence, neuron, hippocampal neuron

Abstract

This protocol describes the immunofluorescence staining of cultured neurons.

Materials

Reagents

- 4% Paraformaldehyde (PFA) in 0.1 M phosphate buffer, pH 7.4
- PBS (phosphate-buffered saline)
- Normal goat serum (NGS) (G9023 Sigma-Aldrich)
- Triton X-100 (BP151, Fisher Scientific)
- Primary antibodies (user-defined)
- Secondary antibodies (Alexa Fluor series or others)
- ProLong Gold Antifade Mountant (P36930, Thermo Fisher Scientific)

Consumables

- Glass coverslips
- Culture dishes

Equipment

- Nikon Ti microscope, CSU-W1 spinning disk confocal unit, Andor Zyla 4.2 sCMOS camera
- 60× oil immersion objective (PlanApo, NA 1.40)
- Micro-Manager software (v2.0-gamma)
- ImageJ / Fiji (NIH)



Fixation

- 1 Remove culture medium and rinse coverslips briefly with PBS.
- 2 Fix neurons with 4% PFA in 0.1 M phosphate buffer for 10–15 min at room temperature.
- 3 Wash coverslips 3 times with PBS.

Permeabilization and Blocking

- 4 Incubate coverslips for 15–30 min in PBS containing: 4% NGS, 0.1% Triton X-100

Primary Antibody Incubation

- 5 Prepare primary antibodies in PBS containing: 1% NGS, 0.025% Triton X-100
- 6 Incubate coverslips overnight at 4°C in a humidified chamber.
- 7 Wash coverslips 3 times with PBS.



Secondary Antibody Incubation

- 8 Prepare secondary antibodies in PBS containing: 1% NGS, 0.025% Triton X-100
Incubate coverslips for 1 h at room temperature, protected from light.
- 9 Wash coverslips 3 times with PBS, protecting from light.



Mounting

- 10 Mount coverslips on slides using ProLong Gold Antifade Mountant.



Imaging

- 11 Visualize fluorescence using a Nikon Ti + CSU-W1 spinning disk confocal system equipped with a 60× oil- immersion (NA 1.4) objective.
- 12 Process images using ImageJ (NIH).

