




Aug 31, 2025

Immunohistochemical detection of porcine epidemic diarrhea virus (PEDV) in formalin-fixed, paraffin-embedded (FFPE) pig tissues

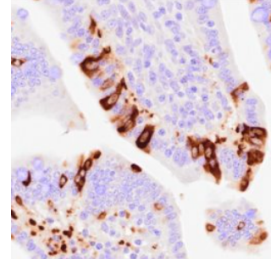

Forked from a private protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.8epv5k2d6v1b/v1>

Jayne Wiarda¹, Eraldo Zanella¹

¹National Animal Disease Center, ARS, USDA

Jayne Wiarda



Jayne E Wiarda

National Animal Disease Center, ARS, USDA

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8epv5k2d6v1b/v1>

Protocol Citation: Jayne Wiarda, Eraldo Zanella 2025. Immunohistochemical detection of porcine epidemic diarrhea virus (PEDV) in formalin-fixed, paraffin-embedded (FFPE) pig tissues. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.8epv5k2d6v1b/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 31, 2025

Last Modified: August 31, 2025

Protocol Integer ID: 226105

Keywords: NP103, immunohistochemical detection of porcine epidemic diarrhea virus, porcine epidemic diarrhea virus, pig tissues an immunohistochemistry, pig tissue, immunohistochemical detection, immunohistochemistry, pedv, tissue, staining protocol

Funders Acknowledgements:

USDA-ARS

Grant ID: CRIS #5030-32000-230-000-D

Disclaimer

All opinions expressed in this paper are the authors' and do not necessarily reflect the policies and views of USDA, ARS, DOE, or ORAU/ORISE. Mention of trade names or products is for information purposes only and does not imply endorsement by the USDA. USDA is an equal opportunity employer and provider.

Abstract

An immunohistochemistry (IHC) staining protocol for *in situ* identification of porcine epidemic diarrhea virus (PEDV) in formalin-fixed, paraffin-embedded (FFPE) tissue.

Attachments



IHC_Pig_PEDV_ManualA.

⌵

4.3MB

Guidelines

	A	B	C	D	E	F	G	H
	Antibody	Undiluted	100ug/ml	10ug/ml	1ug/ml	0.1ug/ml	0.01ug/ml	0.001ug/ml
	MPEDV-5A-2F7	3.6 mg/ml	1:36	1:360	1:3,600	1:36,000	1:360,000	1:3,600,000
	MPEDV-5A-3D6	5.4 mg/ml	1:54	1:540	1:5,400	1:54,000	1:540,000	1:5,400,000
	MPEDV-5A-5B9	2.62 mg/ml	1:26.2	1:262	1:2,620	1:26,200	1:262,000	1:2,620,000
	MPEDV-5A-6H10	1.95 mg/ml	1:19.5	1:195	1:1,950	1:19,500	1:195,000	1:1,950,000
	MPEDV-5A-7A12	1.75 mg/ml	1:17.5	1:175	1:1,750	1:17,500	1:175,000	1:1,750,000

Primary antibody dilutions

Assay Controls

- Controls will be required to ensure there is minimal/no cross-reactivity with off-target epitopes in the tissue and to ensure stains are specific to the intended antibody targets
- IHC controls:
 - Secondary only control
 - This slide receives only diluent in place of primary antibody. All other reagents are still applied to the slide.

Assay Variations

- Parameters may need to be further optimized for different experiments, tissues, targets, or species.

Materials

Equipment

- Pipettes/pipette tips
- Drying oven (able to reach & hold 60°C)
- Fume hood
- Slide staining tray (e.g. Simport M920-2)
- EZ-Batch Slide Holder (Advanced Cell Diagnostics 321716)
- EZ-Batch Wash Tray (Advanced Cell Diagnostics 321717)
- Tissue-Tek Vertical 24 slide rack (American Master Tech Scientific LWS2124)
- Tissue-Tek Staining Dishes (American Master Tech Scientific LWS20WH)
- Tissue-Tek Clearing Agent Dishes, xylene resistant (American Master Tech Scientific LWS20GR)
- Brightfield microscope and/or slide imaging platform

Reagents/Supplies

- Distilled water (obtained in-house)
- Phosphate-buffered saline (PBS), pH 7.2-7.4 (made in-house)
- 0.05% PBS-Tween (PBS-T), pH 7.35 (made in-house)
- Xylenes (Macron Fine Chemicals 8668-16)
- 100% ethanol (Pharmco 111000200)
 - Dilute with distilled water to make 95%, 85%, and 70% concentrations
- Pro-Par Clearant (Anatech 510)
- Fixative
 - 10% NBF (Cancer Diagnostics, Inc. 111) or 4% PFA (Electron Microscopy Sciences 15713)
- ImmEdge Hydrophobic Barrier Pen (Vector H-4000)
- Ready to use Proteinase-K (Dako S302080-2/S302030-2)
- BLOXALL Endogenous Blocking Solution (Vector SP-6000)
- ImmPRESS-VR Horse Anti-Mouse IgG Polymer Detection Kit, Peroxidase (Vector MP-6402-15)
 - Normal Horse Serum, 2.5%
 - ImmPRESS-VR Horse Anti-Mouse IgG Polymer Reagent
- Mouse anti-PEDV antibody
 - Clone #2F7 (Immunology Consultants Laboratory MPEDV-5A-2F7, stock concentration 3.6mg/mL)
 - Clone #3D6 (Immunology Consultants Laboratory MPEDV-5A-3D6, stock concentration 5.4mg/mL)
 - Clone #5B9 (Immunology Consultants Laboratory MPEDV-5A-5B9, stock concentration 2.62mg/mL)
 - Clone #6H10 (Immunology Consultants Laboratory MPEDV-5A-6H10, stock concentration 1.95mg/mL)
 - Clone #7A12 (Immunology Consultants Laboratory MPEDV-5A-7A12, stock concentration 1.75mg/mL)
- 1% bovine serum albumin (BSA) in PBS (made in-house)
- ImmPACT DAB EqV Substrate Kit, Peroxidase (Vector SK-4103)
 - ImmPACT DAB EqV Reagent 1 (Chromogen)
 - ImmPACT DAB EqV Reagent 2 (Diluent)
- Gill's Hematoxylin I (American Master Tech Scientific HXGHE1LT)
- Refrax Mounting Medium (Anatech 711)
- #1 thickness cover glass (e.g. Fisherbrand 12-545-F)



Safety warnings

For all reagents, refer to MSDS to determine appropriate precautions, personal protective equipment (PPE), and disposal methods before use

Before start

Starting specimens

Starting samples = FFPE tissues cut to 4 micron thickness and adhered to positively-charged microscopy slides (e.g. SuperFrost Plus Slides; Fisher Scientific 12-550-15). It is crucial that tissues are adequately fixed to prevent tissue degradation. Tissues no thicker than 0.5 centimeters should be freshly harvested and placed into 10% neutral-buffered formalin (NBF) or 4% paraformaldehyde (PFA) at a ratio of at least 20 volumes fixative per one volume tissue. Fix tissues between 16-30 hours at room temperature (RT), followed by immediate transfer to 70% ethanol and processing into FFPE tissue blocks. Fixation times should be optimized for individual tissues and experiments.



Baking

20m

1 Before starting the assay:

- Preheat a dry oven to 60°C
- Load slides for assay into vertical slide rack

2 • **Bake slides 20 min 60°C** ⌚ 00:20:00

20m

2.1 While slides bake:

- o Prepare 0.05% PBS-T

2.2 Immediately before deparaffinizing:

- o Add ~200 mL xylenes to each of three clearing agent dishes in a fume hood
- o Add ~200 mL 100% ethanol to each of two staining dishes in a fume hood
- o Add ~200 mL 95% ethanol to a staining dish in a fume hood
- o Add ~200 mL 85% ethanol to a staining dish in a fume hood
- o Add ~200 mL 70% ethanol to a staining dish in a fume hood
- o Add ~200 mL distilled water to a staining dish in a fume hood
- o Add ~200 mL PBS-T to a staining dish

Deparaffinizing & Rehydrating

20m

3 • Submerge slide rack in fresh **xylenes 5 min RT** ⌚ 00:20:00

20m

- Submerge slide rack in fresh **xylenes 5 min RT**
- Submerge slide rack in fresh **xylenes 5 min RT**
- Submerge slides rack in fresh **100% ethanol 1 min RT**
- Submerge slides rack in fresh **100% ethanol 1 min RT**
- Submerge slides rack in fresh **95% ethanol 1 min RT**
- Submerge slides rack in fresh **85% ethanol 1 min RT**
- Submerge slides rack in fresh **70% ethanol 1 min RT**
- Submerge slides rack in fresh **distilled water 3 min RT**
- Submerge slides rack in fresh **PBS-T for transport**

3.1 While slides deparaffinize/rehydrate:

- o Turn off dry oven
- o Prepare humidified slide staining tray by adding water to bottom & placing lid on top
- o Add ~200 mL PBS-T to each of two wash trays

Hydrophobic Barrier

20m



- 4
- **Apply hydrophobic barrier** around each tissue  00:20:00 20m
 - o One by one, unload slides from vertical rack submerged in PBS-T. Dry off only the area around the tissue where a barrier will be drawn with a hydrophobic barrier pen. Keep tissue area wet the whole time. Draw barrier and place slides into the EZ-Batch slide holder placed inside the slide staining tray. Using a pipette, apply a small amount of PBS-T within the barrier (just enough to keep the tissue wet while drawing barriers on remaining slides). Slides will remain locked in the EZ-Batch slide holder throughout the protocol until being transferred back to a vertical rack for counterstaining.
 - Leave slide holder in slide staining tray

Proteinase Digestion

10m

- 5
- Decant slide holder and again place flat in slide staining tray  00:10:00 10m
 - Incubate with **Dako Ready to use Proteinase-K 3 min RT**
 - o Apply drops to completely cover tissues
 - o Let incubate in slide staining tray with lid closed
 - Decant slide holder and transfer to wash trays for PBS-T washes
 - Submerge slide holder in fresh **PBS-T 2 min RT**
 - Submerge slide holder in fresh **PBS-T 2 min RT**
- 5.1 While slides incubate with Proteinase:
- o Discard deparaffinizing & rehydrating reagents
 - o Add ~200 mL PBS-T to each of two wash trays

Tissue Quenching

15m

- 6
- Decant slide holder and again place flat in slide staining tray  00:15:00 15m
 - Incubate with **BLOXALL Endogenous Blocking Solution 10 min RT**
 - o Apply to completely cover tissues; let incubate in slide staining tray with lid closed
 - Decant slide holder and transfer to wash trays for PBS-T washes
 - Submerge slide holder in fresh **PBS-T 2 min RT**
 - Submerge slide holder in fresh **PBS-T 2 min RT**
- 6.1 While slides incubate with enzyme block:
- o Discard epitope retrieval reagents
 - o Add ~200 mL PBS-T to each of two wash trays

Protein Blocking

30m

- 7
- Decant slide holder and again place flat in slide staining tray  00:30:00 30m
 - Incubate with **Normal Horse Serum 2.5% 30 min RT**



- o Serum is a component of the secondary antibody reagent kit
- o Apply to completely cover tissues; let incubate in slide staining tray with lid closed

7.1 While slides incubate with serum:

- o Discard tissue quenching reagents
- o Prepare diluted primary antibody by adding antibody stock to 1% BSA in PBS at a dilution of 1:10,000. Total volume to use is dependent on tissue sizes. Make sure to mix reagents before pipetting. Mix diluted antibody well before use.

Primary Antibody

5m

- 8
- Decant slide holder and again place flat in slide staining tray  Overnight
 - o Do not wash with PBS-T between decanting serum and applying secondary antibody
 - Incubate with **diluted primary antibody 4C overnight**
 - o Apply to completely cover tissues; let incubate in slide staining tray with lid closed
 - • Decant slide holder and transfer to wash trays for PBS-T washes
 - Submerge slide holder in fresh **PBS-T 2 min RT**
 - Submerge slide holder in fresh **PBS-T 2 min RT**

5m

8.1 While slides incubate with primary antibody:

- o Discard protein blocking reagents
- The Next Day:
- o Add ~200 mL PBS-T to each of two wash trays

Secondary Antibody

35m

- 9
- Decant slide holder and again place flat in slide staining tray  00:35:00
- Incubate with **Horse anti Mouse IgG (HRP-conjugated polymer) 30 min RT**
 - o Secondary antibody polymer is a component of the secondary antibody reagent kit
 - o Apply to completely cover tissues; let incubate in slide staining tray with lid closed
 - Decant slides and transfer to vertical slide rack for PBS-T washes
 - Submerge slide rack in fresh **PBS-T 2 min RT**
 - Submerge slide rack in fresh **PBS-T 2 min RT**

35m

9.1 While slides incubate with secondary antibody:

- o Discard primary antibody reagents
- o Add ~200 mL PBS-T to each of two wash trays



9.2 Immediately before chromogen detection:

- o Prepare diluted DAB chromogen by adding equal volumes of DAB Reagent 1 (Chromogen) and DAB Reagent 2 (Diluent) together. Total volume to use is dependent on tissue sizes. Make sure to mix reagents thoroughly. Store in the dark due to light sensitivity.

Chromogenic Detection (DAB)

12m

- 10
- Decant slide holder and again place flat in slide staining tray  00:12:00
 - Incubate with diluted **DAB chromogen 7 min RT**
 - o Apply to completely cover tissues; let incubate in slide staining tray with lid closed
 - Decant slide holder and transfer to vertical slide rack for PBS-T washes
 - Submerge slide rack in fresh **PBS-T 2 min RT**
 - Submerge slide rack in fresh **PBS-T 2 min RT**

12m

10.1 While slides incubate with chromogen:

- o Discard remaining secondary antibody reagents
- o Add ~200 mL PBS-T to each of two wash trays
- o Add ~200 mL hematoxylin to one staining dish
- o Add ~200 mL distilled water to each of three staining dishes

Counterstaining

2m

- 11
- Transfer slides to vertical slide rack  00:02:00
 - o Do quickly to avoid drying out slides or alternatively place vertical slide rack in a staining dish containing PBS-T and then transfer slides
 - Submerge slide rack in **hematoxylin 1 min RT**
 - Submerge slide rack in fresh **distilled water, dunking 3-5 times**
 - Submerge slide rack in fresh **distilled water, dunking 3-5 times**
 - Submerge slide rack in fresh **distilled water, dunking 3-5 times**
 - o Water should no longer appear purple in the third water dish used

2m

Mounting

45m

- 12
- Submerge slide rack in fresh **95% ethanol 1 min RT**  00:45:00
 - Submerge slide rack in fresh **100% ethanol 1 min RT**
 - Submerge slide rack in fresh **100% ethanol 1 min RT**
 - Submerge slide rack in fresh **100% ethanol 1 min RT**
 - Submerge slide rack in fresh **Pro-Par 5 min RT**
 - Submerge slide rack in fresh **Pro-Par 5 min RT**
 - Submerge slide rack in fresh **Pro-Par 5 min RT**

45m



- **Mount slides** by adding 2-4 drops of mounting media to each slide, followed by application of a cover glass. Remove bubbles from tissue by applying pressure to cover glass.
- Place slides flat in a dry, dark space to air dry at RT overnight
- Assess staining with a brightfield microscope

12.1 While slides are air drying:

- o Discard chromogen detection, counterstaining, and slide mounting reagents

Protocol references

- Staining protocol was developed by Dr. Jayne Wiarda
- Staining protocol was optimized and executed by Dr. Jayne Wiarda, Eraldo Zanella, and Colin Stoy
- We thank Adrienne Shircliff for slide sectioning and imaging