



Aug 15, 2024

Immunohistochemical staining of wholemount major pelvic ganglia (MPG) for analysis of myelinated bladder afferents

Forked from a private protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.yxmvme3nbg3p/v1>

Janet R Keast¹, Peregrine B Osborne¹, Nicole Wiedmann¹

¹University of Melbourne

SPARC

Tech. support email: info@neuinfo.org



John-Paul Fuller-Jackson

University of Melbourne

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.yxmvme3nbg3p/v1>

Protocol Citation: Janet R Keast, Peregrine B Osborne, Nicole Wiedmann 2024. Immunohistochemical staining of wholemount major pelvic ganglia (MPG) for analysis of myelinated bladder afferents. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.yxmvme3nbg3p/v1>



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 23, 2024

Last Modified: August 15, 2024

Protocol Integer ID: 100336

Keywords: myelinated bladder afferent, immunohistochemical staining of wholemount major pelvic ganglia, myelinated axon, major pelvic ganglia, wholemount major pelvic ganglia, immunohistochemical staining, neurofascin, peripheral neuron, bladder afferents this protocol, immunohistochemical procedure, bladder, neuron, bladder afferent

Funders Acknowledgements:

NIH SPARC

Grant ID: 3OT2OD023872

Abstract

This protocol describes immunohistochemical procedures applied to wholemount major pelvic ganglia (MPG) for the visualization of myelinated axons filled with cholera toxin subunit B (CTB) and/or neurons transfected with adeno-associated virus (AAV) expressing TdTomato. In this protocol, samples were obtained from rats in which CTB was microinjected into the bladder, and AAV-PHP.S was intravenously administered to preferentially and sparsely label peripheral neurons. In this context, neurofascin (paranode marker) identifies myelinated axons.

Materials

- ☒ Horse serum **Merck MilliporeSigma (Sigma-Aldrich) Catalog #12449C**
- ☒ Triton X-100 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #T8787-50ML**
- ☒ Anti-RFP antibody (guinea pig) **Synaptic Systems Catalog #390004**
- ☒ Anti-cholera toxin subunit B antibody (goat) **List Labs Catalog #703**
- ☒ Anti-neurofascin antibody (rabbit) **Alomone Labs Catalog #AIP-025**
- ☒ Cy3 Donkey anti-guinea pig IgG **Jackson ImmunoResearch Laboratories, Inc. Catalog #706-165-148**
- ☒ AF488 Donkey anti-goat IgG **Jackson ImmunoResearch Laboratories, Inc. Catalog #705-545-147**
- ☒ AF647 Donkey anti-rabbit antibody **Invitrogen Catalog #A32795**

Solutions:

- PBS: phosphate-buffered saline, 0.1 M, pH 7.2
- PBS containing 0.1% sodium azide
- PB: phosphate-buffer, 0.1M, pH7.2
- Blocking solution: PBS containing 10% normal horse serum and 0.5% triton X-100
- PBS containing 0.1% sodium azide, 2% normal horse serum and 0.5% triton X-100

Primary Antibodies:

	A	B	C	D	E
	Abbreviation	Synonym	RRID	Host Species	Dilution
	RFP	Red fluorescent protein	AB_2737052	Guinea pig	1:500
	CTB	Cholera toxin subunit B	AB_10013220	Goat	1:10,000
	Neurofascin	NF155	AB_2756657	Rabbit	1:1000

Secondary Antibodies:

	A	B	C	D
	Tag-antibody	Host Species	RRID	Dilution



	A	B	C	D
	Cy3 anti-guinea pig	Donkey	AB_2340460	1:2000
	AF488 anti-goat	Donkey	AB_2336933	1:1000
	AF647 anti-rabbit	Donkey	AB_2762835	1:1000

Immunohistochemistry

- 1 Wash whole MPGs in phosphate buffer (PB; 0.1M; pH 7.2) (3 × 30 min)
- 2 Incubate sections in blocking solution (PB; 10% horse serum; and 0.5% Triton-X) at room temperature for 2 h
- 3 Incubate sections in appropriate dilutions of primary antibodies (or combinations of primary antibodies) for 72h. Antibodies are diluted in PBS containing 0.1% sodium azide, 2% horse serum, and 0.5% triton-X.
- 4 Wash tissue in PBS (3 × 30 min)
- 5 Incubate sections in appropriate dilutions of secondary antibodies (or combinations of secondary antibodies) 24 h in the dark. Antibodies are diluted in PBS containing 2% horse serum, and 0.5% triton-X.
- 6 Wash tissue in PBS (3 × 30 min)
- 7 Mount tissue onto glass slides and coverslip in preferred anti-fade mountant.