

Mar 14, 2023

(3) Immunohistochemistry of liver tissue sections



Forked from Immunohistochemistry of liver tissue sections

DOI

dx.doi.org/10.17504/protocols.io.36wgqjpkxvk5/v1

Presha Rajbhandari¹, Taruna Neelakantan¹, Brent R. Stockwell¹

¹Columbia University

Human BioMolecular Atlas Program (HuBMAP) Method Development Community Tech. support email: Jeff.spraggins@vanderbilt.edu



Presha Rajbhandari

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account





DOI: https://dx.doi.org/10.17504/protocols.io.36wgqjpkxvk5/v1

Protocol Citation: Presha Rajbhandari, Taruna Neelakantan, Brent R. Stockwell 2023. Immunohistochemistry of liver tissue sections. **protocols.io** https://dx.doi.org/10.17504/protocols.io.36wgqjpkxvk5/v1

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: March 14, 2023

Last Modified: March 14, 2023

Protocol Integer ID: 78790

Keywords: IHC, immunohistochemistry, frozen human liver tissue sample, standard immunohistochemistry for antibody validation, liver tissue sections this protocol, standard immunohistochemistry, liver tissue section, antibody validation, molecular pathology core facility at columbia university, molecular pathology core facility, antibody

Funders Acknowledgements:

NIH

Grant ID: 4UH3CA256962

Abstract

This protocol outlines the steps used to perform standard immunohistochemistry for antibody validation in frozen human liver tissue samples, performed at Molecular Pathology Core facility at Columbia University.

Troubleshooting



- 1 Cryosection the frozen liver tissue at 5µm thickness and place it on charged slide
- Air dry the sections for 3 minutes 3 03:00:00
- Fix the tissue sections in cold acetone for 15 minutes 00:15:00. Alternatively, fix with 1% paraformaldehyde at 4C for 15min followed by ice-cold methanol at -20C for 5min.
- Air dry the sections at room temperature for 2 minutes 00:02:00
- Incubate the slides in 0.3% hydrogen peroxide in PBS for 5 minutes 00:05:00 to block peroxidase activity
- Wash slides with PBS 3 times, for 5 minutes each 00:05:00 X3
- 7 1. Block the tissue sections in 10% normal goat serum or 5%Horse serum with 0.1%BSA for 20 minutes 00:20:00
- 8 Incubate the tissue sections with primary antibody diluted in DAKO antibody diluent at room temperature for 1.5-2 hours 01:30:00 02:00:00
- 9 Wash the slides with PBS 3 times, for 5 minutes each 60 00:05:00 each 3 times
- Incubate the tissue sections with biotinylated secondary antibody diluted in PBS at room temperature for 30-45 minutes 00:30:00 00:45:00
- 11 Wash the slides with PBS 3 times, for 5 minutes each 00:05:00 each 3 times
- 12 Incubate the tissue sections with ABC (Avidin-Biotin complex) peroxidase solution at room temperature for 30 minutes 00:30:00

15m



- 13 Wash the slides with PBS 3 times, 5 minutes each 00:05:00 each - 3 times
- 14 Incubate the tissue sections with DAB (3,3'-diaminobenzidine) peroxidase substrate solution until desired color intensity is reached and immerse slides in distilled water
- 15 Counterstain with hematoxylin and rinse with distilled water
- 16 Dehydrate the sections using 95% ethanol followed by 100% ethanol
- 17 Clear with xylene and mount coverslip using mounting medium