

Nov 27, 2023

Immunohistochemistry Protocol for Paraffin-Embedded Sections - General

DOI

<https://dx.doi.org/10.17504/protocols.io.kqdg3p7mql25/v1>

Julianna Tomlinson¹, Nathalie Lengacher²

¹Ottawa Hospital Research Institute; ²OHRI



Julianna Tomlinson

Ottawa Hospital Research Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg3p7mql25/v1>

Protocol Citation: Julianna Tomlinson, Nathalie Lengacher 2023. Immunohistochemistry Protocol for Paraffin-Embedded Sections - General. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.kqdg3p7mql25/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 17, 2022

Last Modified: June 01, 2024

Protocol Integer ID: 62744

Keywords: Immunohistochemistry Protocol, Antigen retrieval, ASAPCRN, immunohistochemistry protocol for paraffin, immunohistochemistry protocol, general immunohistochemistry procedure, immunohistochemistry, paraffin, general this protocol detail, embedded section, protocol detail

Abstract

This protocol details general immunohistochemistry procedure.

Attachments



[gegpbqbbhp.pdf](#)

183KB

Materials

For DAB:

Solution-based: Sigma Cat# D5637; prepared in Tris buffer as per manufacturer's instructions

SIGMAFAST DAB Tablets: Sigma Cat# 4293

SIGMAFAST DAB with Metal Enhancer - Sigma Cat# 0426



Day 1 - Deparaffinize

- 1 Dip slides in 100% Citrisolv 3 × 5 mins.
- 1.1 Dip slides in 100% Citrisolv 3 x  00:05:00 (1/3). 5m
- 1.2 Dip slides in 100% Citrisolv 3 x  00:05:00 (2/3). 5m
- 1.3 Dip slides in 100% Citrisolv 3 x  00:05:00 (3/3). 5m

Hydration

- 2 Dip in 100% EtOH 2 × 2 mins.
- 2.1 Dip in 100% EtOH 2 x  00:02:00 (1/2). 2m
- 2.2 Dip in 100% EtOH 2 x  00:02:00 (2/2). 2m
- 3 Dip in 95% EtOH 2 × 2 mins.
- 3.1 Dip in 95% EtOH 2 x  00:02:00 (1/2). 2m
- 3.2 Dip in 95% EtOH 2 x  00:02:00 (2/2). 2m
- 4 Dip in 80% EtOH 2 × 2 mins.
- 4.1 Dip in 80% EtOH 2 x  00:02:00 (1/2). 2m



4.2 Dip in 80% EtOH 2 x 00:02:00 (2/2).

2m

5 Dip in ddH₂O 2 × 2 mins.

5.1 Dip in ddH₂O 2 x 00:02:00 (1/2).

2m

5.2 Dip in ddH₂O 2 x 00:02:00 (2/2).

2m

Quench endogenous peroxides

6 Incubate slides in 0.3% H₂O₂ in methanol for 00:15:00 (500 µL of 30% H₂O₂ in 50 mL MetOH, in white containers).

15m



Antigen retrieval

7

5m

Note

Use citrate buffer 6 or Tris-EDTA buffer 9 based on antigen desired.

Pre-heat antigen retrieval solution in microwave in white containers (~ 50 mL), for 00:05:00 at full power. Use large plastic container and fill halfway with water.

8 Transfer slides to white containers with AR solution. Microwave for 00:10:00 at power level 1.

10m

9 Rest slides at Room temperature for 00:20:00 , change water in bottom of container.

20m



Wash

10 Remove AR solution and wash slides with ddH₂O for 00:05:00 .

5m



11 Draw bubble with hydrophobic pen during this wash.



12 Rinse slides with PBS for 00:05:00 .

5m



Blocking

13 Block with 5% normal goat serum in PBS + 0.1% Triton-x-100 + 0.05% Tween-20, 00:30:00 @ Room temperature (300 µL /slide).

30m

Primary antibody

14 Dilute antibody in 5% normal goat serum in PBS + 0.1% Triton-x-100 + 0.05% Tween-20, Overnight at 4 °C (300 µL /slide).

5m

Day 2 - Wash

15 Rinse with PBS; 3 × 5 minutes



15.1 Rinse with PBS; 3 x 00:05:00 (1/3).

5m

15.2 Rinse with PBS; 3 x 00:05:00 (2/3).

5m

15.3 Rinse with PBS; 3 x 00:05:00 (3/3).

5m

Secondary antibody



16 Add secondary antibody diluted 1:225 in PBS + 0.1% triton-X + 0.05% tween-20 with 5% normal goat serum (~  300 μ L /slide).



16.1 Commonly: biotinylated anti-rabbit IgG (H+L) made in goat (Vector Labs,BA-1000) or biotinylated anti mouse IgG (H+L) made in goat (Vector Labs, BA-9200).

17 Incubate at  Room temperature for  01:00:00 -  02:00:00 .

3h



Wash

18 Rinse with PBS; 3 $\times$ 5 minutes.



18.1 Rinse slides with PBS; 3 $\times$  00:05:00 (1/3).

5m

18.2 Rinse slides with PBS; 3 $\times$  00:05:00 (2/3).

5m

18.3 Rinse slides with PBS; 3 $\times$  00:05:00 (3/3).

5m

Amplification

19 Prepare amplification solution at least  00:30:00 before use.

30m

20 Amplification with VECTASTAIN® Elite® ABC HRP Kit.

20.1  9 μ L solution A per mL PBS, vortex; +  9 μ L solution B per mL PBS, vortex; (
 300 μ L /slide).

20.2 Incubate at  Room temperature for  01:00:00 -  02:00:00 .

3h





Wash

21 Rinse slides with PBS; 3 × 5 minutes

21.1 Rinse slides with PBS; 3 x  00:05:00 (1/3).

5m

21.2 Rinse slides with PBS; 3 x  00:05:00 (2/3).

5m

21.3 Rinse slides with PBS; 3 x  00:05:00 (3/3).

5m

Substrate development

22

Note

Use DAB or eDAB depending on antigen being detected.

For regular DAB:

22.1 Add  1 mL of DAB to  150 mL of [M] 50 millimolar (mM) Tris +  50 µL 30% 
H₂O₂.

22.2 Dip slides into DAB solution, develop until colour appears ~  00:03:00 .

3m

23 For eDAB (SIGMAFAST™ DAB with Metal Enhancer):

23.1 Add 1 tablet DAB/Cobalt and 1 tablet Urea Hydrogen Peroxide to  5 mL ddH₂O, 
vortex until dissolved.

**Note**

Prepare immediately before use.

23.2 Use ~  300 μ L /slide, develop until colour appears ~  00:03:00 .

3m



24 Rinse with PBS to stop reaction.



25 Rinse with ddH₂O for  00:05:00 .

5m

**Counterstain**

26 Counterstain with 1:2 Hematoxylin:ddH₂O, dip slides in quickly.

27 Remove excess hematoxylin in ddH₂O; 2X.

Dehydration

28 5 dips, 1 minute each:

28.1 Dip in 50% EtOH  00:01:00

1m

28.2 Dip in 80% EtOH x2  00:01:00 .

1m

28.3 Dip in 95% EtOH  00:01:00 .

1m

28.4 Dip in 100% EtOH x2  00:01:00 .

1m



28.5 Dip in 1:1 Citrisolv: 100% EtOH  00:01:00 .

1m

28.6 Dip in 100% Citrisolv x2  00:01:00 .

1m

Coverslip

29 Coverslip with permount solution, allow slides to dry before viewing on microscope.