



Dec 16, 2024

## Immunohistochemistry with Rodent Brain Sections

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l292kpv1y/v1>

Mukesh Kumar<sup>1</sup>, Yumei Wu<sup>2,3,4</sup>, Pietro De Camilli<sup>2,3,4</sup>, Timothy A. Ryan<sup>1,2</sup>

<sup>1</sup>Department of Biochemistry, Weill Cornell Medicine, New York, NY 10065;

<sup>2</sup>Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, Maryland 20815, USA;

<sup>3</sup>Department of Neuroscience, Howard Hughes Medical Institute, Program in Cellular Neuroscience, Neurodegeneration and Repair, Yale University School of Medicine, New Haven, Connecticut 06520, USA;

<sup>4</sup>Department of Cell Biology, Yale University School of Medicine, New Haven, Connecticut 06520, USA



**Mukesh Kumar**

Weill Cornell Medical College

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.4r3l292kpv1y/v1>

**Protocol Citation:** Mukesh Kumar, Yumei Wu, Pietro De Camilli, Timothy A. Ryan 2024. Immunohistochemistry with Rodent Brain Sections. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.4r3l292kpv1y/v1>



**License:** This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** November 27, 2024

**Last Modified:** December 16, 2024

**Protocol Integer ID:** 114543

**Keywords:** ASAPCRN, immunohistochemistry with rodent brain section, rodent brain section, detailed stepwise procedure of immunohistochemical analysis, immunohistochemical analysis, rodent brain tissue, immunohistochemistry, regions of the hippocampus, hippocampus, detailed visualization of target antigen, brain, target antigen, antigen

**Funders Acknowledgements:**

NIH

Grant ID: NS036942

NIH

Grant ID: NS11739

ASAP

Grant ID: ASAP-024404

## Abstract

This protocol describes a detailed stepwise procedure of immunohistochemical analysis of rodent brain tissue, enabling detailed visualization of target antigens within regions of the hippocampus.

## Guidelines

This protocol needs prior approval by the users' Institutional Animal Care and Use Committee (IACUC) or equivalent ethics committee.



## Materials



Ketamine hydrochloride/xylazine hydrochloride solution **Merck MilliporeSigma (Sigma-Aldrich) Catalog #K113**

Fixative: 4% Formaldehyde + 0.125% Glutaraldehyde in 0.1 M phosphate buffer



Formaldehyde solution **Merck MilliporeSigma (Sigma-Aldrich) Catalog #F8775-25ML** +



Glutaraldehyde 50% EM Grade Aqueous **Electron Microscopy Sciences Catalog #16320** in 0.1 M phosphate buffer (PB), warmed to  37 °C

Normal goat serum:



Normal Goat Serum **Jackson ImmunoResearch Laboratories, Inc. Catalog #005-000-121**



Bovine Serum Albumin **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A4503**



Triton X-100 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #T8787**

Primary antibodies:  DDHD2 Polyclonal antibody **Proteintech Catalog #25203-1-AP** ,



Synapsin1/2 antibody **Synaptic Systems Catalog #106 004**

Alexa-conjugated secondary antibodies (Thermofisher Scientific)



ProLong™ Gold Antifade Mountant **Invitrogen - Thermo Fisher Catalog #P36934**

Nail polish (clear)

## Equipments:

- Dissection tools
- Perfusion setup
- Vibratome



## Procedure

4h 30m

- 1 Prepare the anesthetic cocktail (Ketamine/xylazine) and administer it to the mouse via intraperitoneal injection to achieve deep anesthesia.

### Note

Confirm the absence of reflexes before proceeding.

- 2 Perform transcardial perfusion with  37 °C warmed fixative (4% formaldehyde + 0.125% glutaraldehyde in  0.1 Mass Percent PB) until the liver and lungs blanch, indicating successful perfusion. 
- 3 Carefully dissect the brain and immerse it in the same fixative  Overnight at  4 °C . 
- 4 Wash the fixed brain in  0.1 Mass Percent PB to remove excess fixative. 
- 5 Use a vibratome to cut coronal sections at 25 µm thickness. Collect sections in  0.1 Mass Percent PB.
- 6 Prepare the blocking solution containing 2% normal goat serum, 2% BSA, 0.1 M PB, and 0.4% Triton X-100 and incubate brain sections in the blocking solution for  01:00:00 at  Room temperature with gentle agitation.   
1h
- 7 Dilute the primary antibodies in the blocking buffer and incubate the sections with the primary antibody solution  Overnight at  4 °C .   
1h
- 8 Wash the sections 3 times in 0.1 M PB, 10 minutes per wash, at room temperature. 



- 8.1 Wash the sections in [IM] 0.1 Mass Percent PB for ⌚ 00:10:00 at 10m  
🌡 Room temperature (1/3).
- 8.2 Wash the sections in [IM] 0.1 Mass Percent PB for ⌚ 00:10:00 at 10m  
🌡 Room temperature (2/3).
- 8.3 Wash the sections in [IM] 0.1 Mass Percent PB for ⌚ 00:10:00 at 10m  
🌡 Room temperature (3/3).
- 9 Incubate the sections with Alexa-conjugated secondary antibodies prepared in the blocking buffer for ⌚ 02:00:00 at 🌡 Room temperature with gentle agitation in the dark to protect fluorescent dyes from photobleaching. 2h
- 10 Rinse the sections briefly in [IM] 0.1 Mass Percent PB and mount on microscope slides.
- 11 Add ProLong Gold Antifade Reagent to each section and cover with a coverslip.
- 12 Seal the edges of the coverslip with nail polish to prevent drying.
- 13 Images were acquired with the Yokogawa spinning disk field scanning confocal system with microlensing (63x, oil immersion objective lens, CSU-W1 SoRa, Nikon) controlled by NIS elements (Nikon) software (RRID:SCR\_014329).

#### Note

- Ensure all steps are performed with care to avoid tissue damage.
- Optimize primary and secondary antibody dilutions for specific targets.
- Protect fluorescently labeled sections from light exposure during and after staining.
- Perform controls to verify antibody specificity and avoid background staining.