


May 27, 2025

🌐 Immunolabeling and Quantitative Analysis of Synapses in Human iPSC-Derived Dopaminergic Neurons

DOI

<https://dx.doi.org/10.17504/protocols.io.eq2lyq7opvx9/v1>

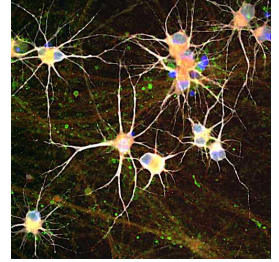

Roxanne Larivière¹, Edward A. Fon¹

¹Montreal Neurological Institute - McGill University



Roxanne Larivière

Montreal Neurological Institute - McGill University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.eq2lyq7opvx9/v1>

Protocol Citation: Roxanne Larivière, Edward A. Fon 2025. Immunolabeling and Quantitative Analysis of Synapses in Human iPSC-Derived Dopaminergic Neurons. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.eq2lyq7opvx9/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: May 02, 2025

Last Modified: May 27, 2025

Protocol Integer ID: 210771

Keywords: iPSC , Dopaminergic neurons, Synapse, Immunofluorescence, Synaptophysin, PSD-95, cellular assay, labeling synapse, derived dopaminergic neurons this protocol, quantitative analysis of synapse, derived dopaminergic neuron, synaptophysin, dopaminergic neuron, synapse, induced pluripotent stem cell, pluripotent stem cell, immunofluorescence, using immunofluorescence, neuron


Funders Acknowledgements:

Michael J. Fox Foundation for Parkinson's research

Grant ID: MJFF-025378

Abstract

This protocol outlines the procedure for labeling synapses with pre-synaptic and post-synaptic markers, synaptophysin and PSD-95 respectively, and microtubule-associated protein 2 (MAP2) in induced pluripotent stem cell (iPSC)-derived dopaminergic neurons using immunofluorescence.

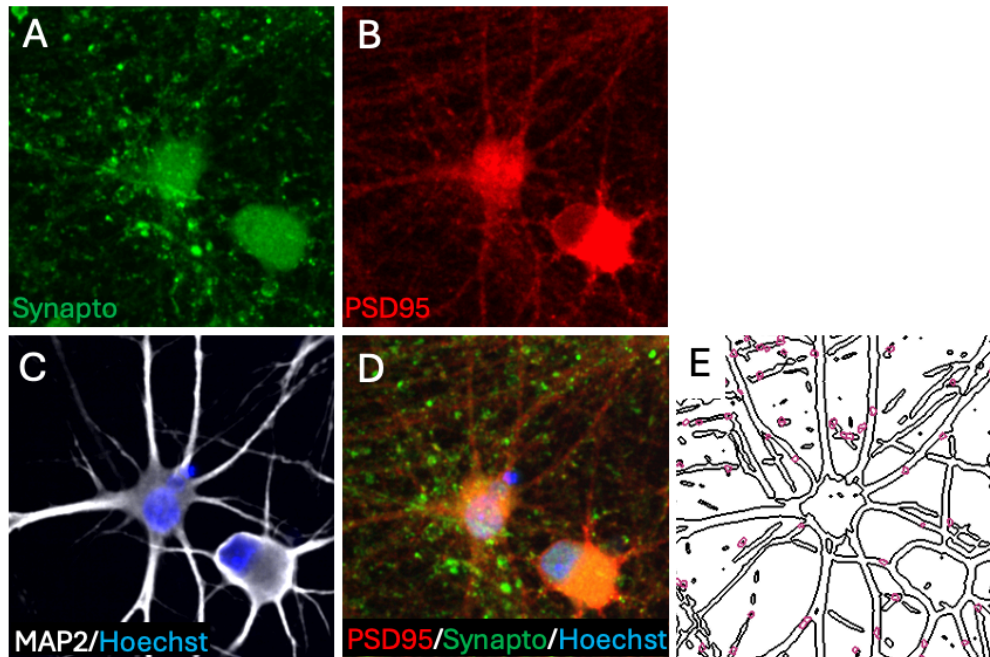

Immunofluorescence labeling of neurons differentiated for 6 weeks was performed using antibodies against Synaptophysin (A), PSD-95 (B), MAP2, and Hoechst (C). (D) shows the merged image of Synaptophysin, PSD-95, and Hoechst staining. (E) displays representative colocalization of Synaptophysin and PSD-95 (highlighted with magenta circles) overlaid on the MAP2 mask following analysis with CellProfiler.



Materials

Reagents	Company	Cat. Number
Rabbit Synaptophysin antibody	Synaptic Systems	101002
Mouse PSD-95 antibody	Millipore	MABN68
Chicken MAP2 antibody	Encor Biotech	CPCA-MAP2
Goat anti-Chicken IgY (H+L) Secondary Antibody, Alexa Fluor 555	ThermoFisher Scientific	A21437
Goat anti-Mouse IgG (H+L) Cross-Adsorbed Secondary Antibody, Alexa Fluor 647	ThermoFisher Scientific	A21235
Goat anti-Rabbit IgG (H+L) Cross-Adsorbed Secondary Antibody, Alexa Fluor 488	ThermoFisher Scientific	A11008
Hoechst 33342	SIGMA	H3570
16% Formaldehyde solution (w/v) Methanol-free	ThermoScientific	28908
Normal goat serum	SIGMA	G9023
Triton X-100	SIGMA	X-100

Before start

iPSC-derived dopaminergic neurons are cultured on 96-well plates for up to 4 to 6 weeks of differentiation. We use black Costar 96-well assay plates (Corning cat. no. 3904), which are compatible with imaging using the Opera Phenix high-content screening system.

Be extremely gentle when dispensing into the wells, as dopaminergic neurons are prone to detaching with forceful pipetting.



Fixation of Dopaminergic Neurons in 96-well plates

35m

- 1
 - Carefully aspirate the culture media, leaving $50\ \mu\text{L}$ in each well.
 - Gently add $50\ \mu\text{L}$ of 8% paraformaldehyde (PFA) to each well to reach a final concentration of 4 % (v/v) PFA.
 - Incubate at room temperature for 00:20:00 .
 - Remove the fixative and wash each well three times with $75\ \mu\text{L}$ of 1X PBS. Allow each wash to incubate for 00:05:00 .

35m

Permeabilization

25m

- 2
 - Prepare the permeabilization solution by diluting Triton X-100 to a final concentration of 0.5 % (v/v) in 1X PBS.
 - Add $60\ \mu\text{L}$ of the permeabilization solution to each well and incubate for 00:10:00 at room temperature.
 - Remove the solution and wash wells three times with $75\ \mu\text{L}$ of 1X PBS, allowing each wash to incubate for 00:05:00 .

25m

Blocking

1h

- 3
 - Prepare the blocking buffer by mixing 5 % (v/v) normal goat serum and 0.02 % (v/v) Triton X-100 in 1X PBS.
 - Add $60\ \mu\text{L}$ of the blocking buffer to each well and incubate for 01:00:00 at Room temperature .

1h

Primary Antibody Incubation

16h

- 4
 1. Prepare the primary antibody solution in blocking buffer with the following dilutions:
 - Rabbit anti-Synaptophysin: 1:500
 - Mouse anti-PSD-95: 1:250
 - Chicken anti-MAP2: 1:1500

16h



2. Add  60 μL of the antibody mix to each well.
3. Incubate  Overnight at  4 °C with gentle shaking.
4. Remove antibody mix and wash four times with  75 μL 1X PBS, allowing each wash to incubate for  00:05:00 .

Secondary Antibody Incubation

2h 15m

- 5
 1. Prepare the secondary antibody solution in  1 % (v/v) normal goat serum diluted in 1X PBS with the following components:
 - Alexa Fluor 488 anti-Rabbit: 1:1000
 - Alexa Fluor 647 anti-Mouse: 1:1000
 - Alexa Fluor 555 anti-Chicken: 1:1000
 - Hoechst 33342: 1:5000
 2. Add  60 μL of the antibody mix to each well.
 3. Incubate for  02:00:00 at  Room temperature with gentle shaking, protected from light.
 4. Remove secondary antibody mix and wash four times with  75 μL 1X PBS, allowing each wash to incubate for  00:05:00 .

2h 15m

Imaging using the Opera Phenix high-content imaging system

- 6
 1. Insert the prepared plate into the Opera Phenix chamber.
 2. Utilize a 40 $\times$ water immersion objective. Capture images using the Hoechst, 488 nm, 555 nm and 647 nm laser channels. Acquire 15-18 fields of view per well

Synapse Detection and Data analysis using Cell Profiler software

- 7 Images are analyzed using a Cell Profiler automated synaptic quantification pipeline, previously published by Berryer et al., eLife, 2023. <https://doi.org/10.7554/eLife.80168>
<https://github.com/mberryer/ALPAQAS>



Protocol references

1. Martin H Berryer, Gizem Rizki, Anna Nathanson, Jenny A Klein, Darina Trendafilova, Sara G Susco, Daisy Lam, Angelica Messana, Kristina M Holton, Kyle W Karhohs, Beth A Cimini, Kathleen Pfaff, Anne E Carpenter, Lee L Rubin, Lindy E Barrett (2023) High-content synaptic phenotyping in human cellular models reveals a role for BET proteins in synapse assembly eLife 12:e80168 <https://doi.org/10.7554/eLife.80168>