

Feb 25, 2025

Immunolabeling of endogenous proteins for fluorescence microscopy in cultured mammalian cells

DOI

<https://dx.doi.org/10.17504/protocols.io.n92ldr1xxg5b/v1>

Ahmad Shami¹, Hera Saqub¹, Samantha C Lewis¹

¹UC Berkeley



Samantha C Lewis

UC Berkeley

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ldr1xxg5b/v1>

Protocol Citation: Ahmad Shami, Hera Saqub, Samantha C Lewis 2025. Immunolabeling of endogenous proteins for fluorescence microscopy in cultured mammalian cells. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.n92ldr1xxg5b/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: February 17, 2025

Last Modified: February 25, 2025

Protocol Integer ID: 120859

Keywords: Immunofluorescence, Mitochondrial protein detection, mitochondria, fluorescence microscopy, imaging endogenous protein, endogenous proteins by immunofluorescence, fluorescence microscopy in cultured mammalian cell, immunolabeling of endogenous protein, endogenous protein, fluorescence microscopy, immunofluorescence, microscopy, mitochondria, cultured mammalian cell, bound organelle, protein, cell, immunolabeling, other membrane, fluorescence, imaging

Funders Acknowledgements:

National Institutes of Health

Grant ID: R35GM147218

Abstract

This protocol details the detection of endogenous proteins by immunofluorescence. This is our standard protocol for imaging endogenous proteins associated with mitochondria and other membrane-bound organelles.

Guidelines

Introduction

This protocol provides detailed instructions for detection of endogenous proteins in fixed cells by immunofluorescence (Immunofluorescence Assay/IFA). The protocol results in a fluorescent sample, to be visualized using a high-resolution fluorescence confocal microscope. Target proteins appear as distinct signals corresponding to the fluorophore-conjugated secondary antibodies, allowing the spatial analysis of subcellular structures, protein interactions, and nucleic acids.

Materials

- Proliferating cancer cell line or fibroblasts at 40% confluency
- 1X Phosphate-buffered saline (PBS), pH 7.4
- Fixative: 4% Paraformaldehyde in PBS, pH 7.4
- Quenching Solution: 1M Glycine in PBS, pH 7.4
- Permeabilization Solution: 0.1% Triton-X 100 in PBS, pH 7.4
- Blocking Solution: 0.1% Tween in 1X TBS + 1% BSA (Bovine Serum Albumin)
- Glass bottom poly-D-lysine-coated dishes (P35GC-1.5-14-C, MatTek)
- Primary antibody (specific to the target protein)
- Secondary antibody conjugated to fluorophore of choice



Method

4d 1h 1m

- 1 Prepare the sample by plating cells in  2 mL of media at a 20% density in a glass bottom poly-D-lysine-coated dish (P35GC-1.5-14-C, MatTek).
- 2 Grow cells for  24:00:00 until they reach a confluency of approximately 40%. 1d
- 3 Aspirate media and fix the cells with  2 mL of pre-warmed 4% PFA in PBS. 
- 4 Incubate for  00:20:00 at  Room temperature on a shaker gently rotating. 20m
 
- 5 Quench the reaction by adding  200 μ L of 1M glycine in PBS to the fixed sample and incubate for  00:01:00 at  Room temperature . 1m
 
- 6 Aspirate the quenched fixative and wash the sample once with PBS.  
- 7 Add  2 mL of 0.1% Triton X-100 in PBS and incubate for  00:10:00 at  Room temperature on a shaker to permeabilize the sample. 10m
  
- 8 Aspirate permeabilized solution and wash sample once with PBS.  
- 9 Add  2 mL of blocking buffer (1% BSA + 0.1% Tween in 1X TBS) with a  00:10:00 incubation at  Room temperature on a shaker. 10m
  
- 10 Aspirate blocking buffer and add a freshly prepared  2 mL primary antibody dilution (typically a 1:1000 dilution) made in blocking buffer. 
- 11 Incubate the sample  Overnight at  4 °C in the dark.   



- 12 Aspirate the primary antibody solution and wash the sample once with  2 mL of blocking buffer.  
- 13 Aspirate wash and add  2 mL of freshly prepared secondary antibody dilution (typically at a 1:2000 dilution) with a  00:10:00 incubation at  Room temperature on a shaker in the dark.    10m
- 14 Aspirate the secondary antibody solution and wash the sample with  2 mL of blocking buffer with a  00:10:00 incubation at  Room temperature on a shaker.    10m
- 15 Remove wash and cover sample with  2 mL PBS. 
- 16 Proceed with fluorescence microscopy. 
- 17 Samples may be stored at  4 °C protected from light for up to  72:00:00 ; bring dishes to  Room temperature before imaging.  3d

Protocol references

- Begeman A, Smolka JA, Shami A, Waingankar TP, Lewis SC. A spatial atlas of mitochondrial gene expression reveals dynamic translation hubs and remodeling in stress. bioRxiv [Preprint]. 2024 Aug 17:2024.08.05.604215. doi: 10.1101/2024.08.05.604215. PMID: 39149346; PMCID: PMC11326164.
- Smolka JA, Lewis SC. In Situ Analysis of Mitochondrial DNA Synthesis Using Metabolic Labeling Coupled to Fluorescence Microscopy. Methods Mol Biol. 2023;2615:99-106. doi: 10.1007/978-1-0716-2922-2_8. PMID: 36807787.

Acknowledgements

We thank the Lewis Lab for helpful discussion.