Sep 05, 2018 Version 1

In vitro α-amylase inhibitory assay V.1

DOI

dx.doi.org/10.17504/protocols.io.s9meh46

Maira Rubi, Maira Rubi Segura Campos¹

¹Universidad Autónoma de Yucatán

Maira Rubi Segura Campos





DOI: dx.doi.org/10.17504/protocols.io.s9meh46

Protocol Citation: Maira Rubi, Maira Rubi Segura Campos 2018. In vitro α-amylase inhibitory assay. **protocols.io** <u>https://dx.doi.org/10.17504/protocols.io.s9meh46</u>

Manuscript citation:

Dineshkumar B, Mitra A, Manjunatha M. Studies on the anti-diabetic and hypolipidemic potentials of mangiferin (xanthone glucoside) in streptozotocin-induced type 1 and type 2 diabetic model rats. Int J Adv Pharm Sci. 2010; 1: 75-85.

License: This is an open access protocol distributed under the terms of the <u>Creative Commons Attribution License</u>, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working We use this protocol and it's working

Created: September 05, 2018

Last Modified: September 05, 2018

Protocol Integer ID: 15373

Abstract

The assay was carried out following the protocol reported by Dineshkumar et al. [8]. Starch (2 mg) was suspended in a tube containing 0.2 mL of 0.5 M Tris-HCI (Sigma-Aldrich, USA) buffer (pH 6.9) with 0.01 M calcium chloride as substrate. The tube was boiled for 5 min and then preincubated at 37 °C for 5 min. Plant aqueous extract (1 mg) was dissolved with 1 mL of 0.1% of dimethyl sulfoxide in order to obtain a concentration of 1,000 μ g/mL; then 0.2 mL of aqueous extract was added to the tube containing the substrate solution, 0.1 mL of porcine pancreatic amylase in Tris-HCI buffer (2 U/mL) was also added, and incubated for 10 min at 37 °C. Finally, the reaction was stopped with 0.5 mL of acetic acid (50% v/v) and centrifuged 5 min at 1,811 × g and 4 °C. The assay was performed in triplicate. The a-amylase inhibitory activity was calculated using the formula (Ac⁺) – (Ac⁻) – (As – Ab)/(Ac⁺) – (Ac⁻) × 100, where Ac⁺, Ac⁻, As, Ab are defined as the absorbance (595 nm) of 100% enzyme activity (only solvent with enzyme), respectively.