

Sep 23, 2023

In vitro kinase assay

 Nature Structural & Molecular Biology

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l225xjl1y/v1>

Elias Adriaenssens¹

¹Sascha Martens lab, University of Vienna, Max Perutz Labs - Vienna



Elias Adriaenssens

Sascha Martens lab, University of Vienna, Max Perutz Labs - ...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l225xjl1y/v1>

External link: <https://doi.org/10.1038/s41594-024-01338-y>

Protocol Citation: Elias Adriaenssens 2023. In vitro kinase assay. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.4r3l225xjl1y/v1>

Manuscript citation:

Elias Adriaenssens, Thanh Ngoc Nguyen, Justyna Sawa-Makarska, Grace Khuu, Martina Schuschnig, Stephen Shoebridge, Marvin Skulsuppaisarn, Emily Maria Watts, Kitty Dora Csalyi, Benjamin Scott Padman, Michael Lazarou, Sascha Martens (2024) Control of mitophagy initiation and progression by the TBK1 adaptors NAP1 and SINTBAD. Nature Structural & Molecular Biology doi: [10.1038/s41594-024-01338-y](https://doi.org/10.1038/s41594-024-01338-y)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 27, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 84091

Keywords: in vitro kinase assay, ASAPCRN, kinase, assay this protocol, assay

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-000350

Abstract

This protocol describes *in vitro* kinase assay.

Attachments



[755-1923.pdf](#)

51KB



Materials

Materials

- Recombinant proteins TBK1, ULK1 complex, and NAP1
- MgCl_2
- ATP
- dH_2O
- Nitrocellulose membranes (RPN132D, GE Healthcare)
- Mini Trans-Blot Cell (Bio-Rad).
- SDS-PAGE gels (NP0321BOX, NP0322BOX, or NP0323BOX, Thermo Fisher)
- PageRuler Prestained protein marker (Thermo Fisher)

Kinase buffer

	A	B
	Tris-HCl pH 7.4	20 mM
	NaCl	150 mM
	DTT	1 mM

Fixation solution

	A
	40% ethanol
	10% acetic acid
	50% dH_2O



In vitro kinase assay

25m

- 1 Mix recombinant proteins TBK1 or ULK1-complex (composed of ULK1, FIP200, ATG13, and ATG101) and NAP1 in kinase buffer. 
- 2 Use the kinases at [M] 50 nanomolar (nM) and mix with [M] 250 nanomolar (nM) NAP1.
- 3 Start the kinase reactions by the adding 2x ATP/MgCl₂ kinase buffer to a final concentration of [M] 10 millimolar (mM) MgCl₂ and [M] 100 millimolar (mM) ATP.
- 4 Prepare protein mixtures as master mixes and divide over the number of time points.
- 5 To control for potential protein instability, induce the latest time point first and then go gradually to the shortest time point.
- 6 In this way, keep all protein mixtures at  Room temperature for the same time, and terminate the reactions together.
- 7 Achieve the termination of reactions by the addition of 6x Protein Loading dye and heat inactivation at  95 °C for  00:05:00 . 
- 8 Separate the samples on 4-12% SDS-PAGE gels (NP0321BOX, NP0322BOX, or NP0323BOX, Thermo Fisher) with PageRuler Prestained protein marker (Thermo Fisher).
- 9 After the run, either stain the SDS-PAGE gel with Coomassie or transfer to nitrocellulose membranes for western blot analysis.
- 10 In the case of Coomassie staining, incubate the gel for  00:10:00 in Coomassie solution, fix for  00:10:00 with fixation solution, and then destain it  Overnight  
- 11 Cut the band corresponding to NAP1 from the gel with a fresh scalpel and submit for mass spectrometry analysis.



- 12 In the case of western blotting, transfer the proteins onto nitrocellulose membranes (RPN132D, GE Healthcare) for  01:00:00 at  4 °C using the Mini Trans-Blot Cell (Bio-Rad).
- 13 Process the membranes further for western blot analysis, as described in the western blot protocol.

1h