

Sep 07, 2023

Version 1

# In vitro LRRK2 kinase activity assay using mass-spectrometry as readout V.1

DOI

[dx.doi.org/10.17504/protocols.io.6qpvr385ovmk/v1](https://dx.doi.org/10.17504/protocols.io.6qpvr385ovmk/v1)

Verena Dederer<sup>1,2</sup>, chatterjeedeep<sup>1,2</sup>, Sebastian Mathea<sup>1,2</sup>, Stefan Knapp<sup>1,2</sup>

<sup>1</sup>Institute of Pharmaceutical Chemistry, Goethe University Frankfurt, Max-von-Laue-Straße 9, Frankfurt 60438, Germany;

<sup>2</sup>Structural Genomics Consortium, Buchman Institute for Molecular Life Science (BMLS), Max-von-Laue-Straße 15, Frankfurt 60438, Germany



Verena Dederer

Goethe University Frankfurt

## Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.6qpvr385ovmk/v1>

**Protocol Citation:** Verena Dederer, chatterjeedeep, Sebastian Mathea, Stefan Knapp 2023. In vitro LRRK2 kinase activity assay using mass-spectrometry as readout. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.6qpvr385ovmk/v1>



**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** September 06, 2023

**Last Modified:** September 07, 2023

**Protocol Integer ID:** 87458

**Keywords:** Irrk2 kinase activity, concentration series of kinase inhibitor, kinase inhibitor, substrate phosphorylation for any purified kinase substrate pair, kinase, purified kinase substrate pair, Irrk2, substrate phosphorylation, phosphorylation,  $\mu$ l reaction, inhibition curves by addition, inhibition curve

**Funders Acknowledgements:**

**Aligning Science Across Parkinson's**

Grant ID: ASAP-000519

## Abstract

This protocol can be used in this form or with small adjustments regarding concentration and reaction time to determine in vitro substrate phosphorylation for any purified kinase substrate pair.

It can be used to determine inhibition curves by addition of concentration series of kinase inhibitors.

Reaction is carried out in 50  $\mu$ L reaction mix containing final concentration of 50 nM kinase and 5 mM substrate.

## Materials

purified protein kinase

purified kinase substrate

buffer: 20 mM Hepes pH 7.4, 150 mM NaCl, 5% glycerol, 0.5 mM TCEP, 20  $\mu$ M GDP, 2.5 mM  $\text{MgCl}_2$

100 mM ATP

## Troubleshooting

## Step-by-Step Protocol

- 1 Prepare 50  $\mu\text{L}$  reaction mix I: 50 nM purified LRRK2 and 5  $\mu\text{M}$  Rab8 substrate in buffer containing 20 mM Hepes pH 7.4, 150 mM NaCl, 5% glycerol, 0.5 mM TCEP, 20  $\mu\text{M}$  GDP, 2.5 mM  $\text{MgCl}_2$ .  
Aliquot 25  $\mu\text{L}$  per tube.

### Note

If you want to run multiple reactions prepare a master mix and aliquot 25  $\mu\text{L}$  per tube.

- 2 Prepare 25  $\mu\text{L}$  reaction mix II: 2 mM ATP + 2.5 mM  $\text{MgCl}_2$  in buffer containing 20 mM Hepes pH 7.4, 150 mM NaCl, 5% glycerol, 0.5 mM TCEP, 20  $\mu\text{M}$  GDP, 2.5 mM  $\text{MgCl}_2$ .

### Note

Multiply if you want to run more than one reaction.

- 3 Prepare 25  $\mu\text{L}$  reaction mix III: same as reaction mix II but no ATP as negative control.
- 4 Start reaction by adding 25  $\mu\text{L}$  reaction II to reaction mix I.  
Note: for negative control add 25  $\mu\text{L}$  reaction mix III to second aliquot of reaction mix I.
- 5 Mix by vortexing and brief centrifugation for 30 sec with 500xg.
- 6 Incubate reaction for 3 h at room temperature.  
Note: depending on kinase this may be shorter or longer.
- 7 Stop reaction by adding 50  $\mu\text{L}$  mass spec buffer ( $\text{dH}_2\text{O}$ +0.1% formic acid).
- 8 Store at  $-80^\circ\text{C}$  or proceed directly with mass spectrometry analysis.



#### Note

For determination of inhibition curves, prepare a concentration series of inhibitor in DMSO. Prepare a master mix of kinase and substrate pair and aliquot 25  $\mu$ L per tube. Add the desired amount of inhibitor to each of the tube and incubate all reactions at room temperature.

#### Note

For analysis: substrate phosphorylation/turnover can be determined as ratio between phosphorylated and unphosphorylated peak intensity for each sample analyzed.