

Jul 10, 2024

In vivo BioID protein purification

DOI

<https://dx.doi.org/10.17504/protocols.io.8epv5r3z6g1b/v1>

Shiyi Wang¹

¹Duke University



Shiyi Wang

Duke University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8epv5r3z6g1b/v1>

Protocol Citation: Shiyi Wang 2024. In vivo BioID protein purification. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.8epv5r3z6g1b/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 10, 2024

Last Modified: August 13, 2024

Protocol Integer ID: 103177

Keywords: ASAPCRN, vivo bioid protein purification, purification, protein

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP) initiative

Grant ID: ASAP-020607

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

The **protocols.io** team notes that research involving animals and humans must be conducted according to internationally-accepted standards and should always have prior approval from an Institutional Ethics Committee or Board.

Abstract

In vivo Bioid protein purification



- 1 ****Animal Preparation**** - Breed genotype-matched animals (wild-type C57BL6 or LRRK2 G2019Ski/ki) to produce single-genotype litters. - For each genotype (WT or G2019S), prepare 6 pups for injection.
- 2 ****AAV Injection**** - Inject 1 μ L of AAVs carrying Astro-Ezrin-BioID (PHP.eB.GfaABC1D-Ezrin WT-BioID2-HA) or Astro-CYTO-BioID (PHP.eB.GfaABC1D-BioID2-HA) bilaterally into the cortex of P0-P2 mouse pups using a Hamilton syringe. - Monitor pups until they recover on a heating pad.
- 3 ****Biotin Injection**** - At P18, P19, and P20, subcutaneously inject biotin at 24 mg/kg to increase biotinylation efficiency.
- 4 ****Tissue Collection**** - At P21, remove the cerebral cortices and store at -80°C . - Pool 2 genotype-matched cortices at the time of protein isolation, yielding 3 independent replicates per BioID construct.
- 5 ****Protein Purification**** - Lyse each cortex in a buffer containing: - 50 mM Tris/HCl, pH 7.5 - 150 mM NaCl - 1 mM EDTA - Protease inhibitor mixture (Roche) - Phosphatase inhibitor mixture (PhosSTOP, Roche) - Add an equal volume of buffer containing: - 50 mM Tris/HCl, pH 7.5 - 150 mM NaCl - 1 mM EDTA - 0.4% SDS - 2% TritonX-100 - 2% deoxycholate - Protease inhibitor mixture - Phosphatase inhibitor mixture
- 6 ****Sonication and Centrifugation**** - Sonicate samples. - Centrifuge at 15,000 g for 10 minutes. - Ultracentrifuge the supernatant at 100,000 g for 30 minutes at 4°C .
- 7 ****Sample Preparation for Protein Binding**** - Add SDS detergent to the samples and heat at 45°C for 45 minutes. - Cool on ice.
- 8 ****Protein Binding**** - Incubate each sample with High-Capacity Streptavidin Agarose beads (ThermoFisher) at 4°C overnight.
- 9 ****Bead Washing**** - Wash beads serially:
 - 10 Twice with a solution containing 2% SDS.
 - 11 Twice with a buffer containing 1% TritonX-100, 1% deoxycholate, and 25 mM LiCl.
 - 12 Twice with 1 M NaCl.



- 13 Five times with 50 mM ammonium bicarbonate.
- 14 ****Protein Elution**** - Elute biotinylated proteins attached to the agarose beads in a buffer containing: - 125 mM Tris/HCl, pH 6.8 - 4% SDS - 0.2% β -mercaptoethanol - 20% glycerol - 3 mM biotin - Heat at 60°C for 15 minutes.
- 15 ****Downstream Analysis**** - Subject the 12 total samples (3 per genotype per construct) to LC-MS/MS and downstream analysis.