

Feb 17, 2026

In Vivo Fiber Photometry Recording With Concurrent Optogenetic Stimulation

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.n2bvj1jzxvk5/v1>

Cristian González-Cabrera¹, Matthias Prigge¹

¹Leibniz Institute for Neurobiology, Magdeburg, Germany



Cristian González-Cabrera

Leibniz Institute für Neurobiologie

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n2bvj1jzxvk5/v1>

Protocol Citation: Cristian González-Cabrera, Matthias Prigge 2026. In Vivo Fiber Photometry Recording With Concurrent Optogenetic Stimulation. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.n2bvj1jzxvk5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: February 16, 2026

Last Modified: February 17, 2026

Protocol Integer ID: 243363

Keywords: vivo fiber photometry recording with concurrent optogenetic stimulation, vivo fiber photometry recording, concurrent optogenetic stimulation of chrimson, optogenetic stimulation, concurrent optogenetic stimulation, wavelength excitation, vta neuron, doric fluorescence minicube system, nm signal channel, expressing neuron

Abstract

This protocol describes in vivo fiber photometry recordings from freely moving DAT-Cre mice expressing jRCaMP1f in VTA neurons. Dual-wavelength excitation (470 nm signal channel and 405 nm isosbestic channel) was sinusoidally modulated and demodulated using a Doric fluorescence minicube system. In some experiments, concurrent optogenetic stimulation of Chrimson-expressing neurons was delivered through the same ferrule.

Guidelines

- Avoid spectral overlap in filter configuration.
- Use non-opsin controls to confirm absence of optical artifacts.

Materials

****Viral Strategy and Implant****

- DAT-Cre mice.
- VTA injection: AAV5-Syn-DIO-jGCaMP8f and AAV8-CAG-FLPX-rc[ChrimsonR-tdTomato].
- PnO injection: retroAAV-Ef1a-FlpO.
- Optical fiber: 400 μm core diameter, NA 0.66, implanted above VTA.

****Recording Setup****

- Doric fluorescence minicube system.
- Silicon photodiode detector.
- Lock-in demodulation via Doric acquisition system.
- Behavioral arena: 20 $\times$ 30 cm.
- Freely moving configuration.

****Excitation Parameters****

GCaMP Signal Channel

- 470 nm excitation.
- Bandpass: 460–490 nm.
- Emission: 500–540 nm.
- Power at ferrule tip: approximately 70 μW .
- Sinusoidal modulation frequency: 208.616 Hz.

Isosbestic Control Channel

- 405 nm excitation.
- Power at ferrule tip: approximately 20 μW .
- Sinusoidal modulation frequency: 333.786 Hz.

****Signal Detection****

- Fluorescence collected through implanted 400 μm fiber.
- Signal detected via silicon photodiode.
- Channels demodulated by Doric system.
- Demodulated signals exported for downstream processing.

****Concurrent Optogenetic Stimulation (When Performed)****

- Chrimson stimulation delivered through the same ferrule.
- Wavelength: 568 nm.
- Power at ferrule tip: approximately 150 μW .
- Delivered via external TTL trigger.

****Optical Separation Safeguards****

- Separate excitation and emission bands for GCaMP.



- Dedicated return paths for green emission.
- 568 nm channel not modulated at 405 or 470 carrier frequencies.
- Non-opsin controls confirmed negligible 568 nm influence on GCaMP detection channel.



Procedure

- 1 Confirm excitation powers at the ferrule tip.
- 2 Connect mouse to patch cord.
- 3 Verify stable fluorescence baseline.
- 4 Start sinusoidal LED modulation.
- 5 Begin demodulated signal acquisition.
- 6 Deliver optogenetic stimulation if applicable.
- 7 Record synchronized stimulation TTLs.

Outputs

- 8 Demodulated 470 nm signal trace.
- 9 Demodulated 405 nm isosbestic trace.
- 10 Stimulation TTL timestamps.
- 11 Continuous behavioral session recording.

Critical Steps



- 12 Maintain stable patch cord coupling.
- 13 Confirm modulation frequencies remain distinct and stable.
- 14 Verify optical power before each session.
- 15 Avoid spectral overlap in filter configuration.
- 16 Use non-opsin controls to confirm absence of optical artifacts.