

Jun 07, 2023

Isolation of bacteria and fungi from cheese rind microbiomes

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzb7b4vx1/v1>



Emily C.P. Weiss¹

¹Arcadia Science

Arcadia Science



Arcadia Science

Arcadia Science

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzb7b4vx1/v1>

Protocol Citation: Emily C.P. Weiss 2023. Isolation of bacteria and fungi from cheese rind microbiomes. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.rm7vzb7b4vx1/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 20, 2023

Last Modified: January 30, 2024

Protocol Integer ID: 75598

Keywords: Bacteria, bacterial, fungus, fungi, microbes, isolate, isolates, cheese, microbial community, community, contaminant, contamination, separate, single, colony, colonies, cheese rind microbiomes this protocol, cheese rind microbiome, fungal isolates from cheese rind, fungal isolate, microbial community sample, cheese rind, bacteria, fungi

Abstract

This protocol describes how to harvest and stock bacterial and fungal isolates from cheese rind microbial community samples.

Image Attribution

Arcadia Science



Materials

- 1.5 mL sterile Eppendorf tubes
- Sterile razor blade
- Sterile 96 well plates (flat bottom)
- Multi-channel pipette
- Solution reservoir (sterile)
- 1 mL and 200 μ L sterile pipette tips
- Sterile toothpicks or wooden dowels
- Sterile plastic pestles
- Sterile 2 mL cryotubes
- Sterile 1 $\times$ phosphate-buffered saline + 20% glycerol
- Sterile 1 $\times$ phosphate-buffered saline + 0.05% Tween
- Sterile 80% glycerol
- LB broth
- LB agar plates with 100 μ g/mL cycloheximide
- PCAMS plates with 50 μ g/mL chloramphenicol

Recipe for 1 L of PCAMS without antibiotic:

5 g tryptone

2.5 g yeast extract

1 g dextrose

1 g whole milk powder

10 g sodium chloride

15 g agar

965.5 mL deionized water



Harvesting the cheese rind biofilm

- 1 If you are starting from a glycerol stock of a previously stored rind community, proceed directly to the dilutions in step 8. If starting from an intact cheese, start at step 2.

Note

This protocol is focused on obtaining pure bacterial and fungal isolates. For a more detailed protocol on sampling and stocking cheese rind communities (for metatranscriptomics, metagenomics, long-term storage, etc.), see <https://www.protocols.io/view/harvesting-and-stocking-cheese-rind-community-samp-ckf2utqe>.

- 2 Holding the cheese in one (gloved) hand and a sterile razor blade in the other, gently scrape the surface of the cheese to remove the rind biofilm. Avoid applying too much pressure, as you will start to dig into the paste. Scrape the rind biofilm into a weigh boat or other container to clean off the blade as needed.

Note

If the cheese surface contains nooks/crannies, use the corner of the razor blade to gently scrape the rind from these.

- 3 After removing the desired amount of rind, use a sterile wooden dowel or a pipette tip to homogenize the harvested rind.

Note

If not harvesting the entire rind, scraping from multiple locations may help to capture more of the microbial diversity present.

Resuspension and dilution

- 4 Add a lentil-sized amount of harvested cheese rind biofilm to 500 μ L of sterile 1 $\times$ phosphate-buffered saline (PBS) + 0.05% Tween in a 1.5 mL microcentrifuge tube.



- 5 Using a small sterile pestle, gently grind the cheese rind sample until it is completely resuspended in the buffer.

Equipment

Disposable Pellet Pestles

NAME

Fisher Scientific

BRAND

12-141-364

SKU

<https://www.fishersci.com/shop/products/fisherbrand-rnase-free-disposable-pellet-pestles-2/p-4246531>

LINK



- 6 Add another 500 μL of PBS + 0.05% Tween to the suspension.

- 7 Vortex thoroughly to mix the suspension.

- 8 Make 10-fold serial dilutions of the cheese rind suspension eight times in PBS + 0.05% Tween to create 10^{-1} , 10^{-2} , 10^{-3} , 10^{-4} , 10^{-5} , 10^{-6} , 10^{-7} , 10^{-8} dilutions.

Note

You can do these serial dilutions in a sterile 96-well plate. If you have multiple samples, you can use a multichannel pipette after the initial dilution to serially dilute all samples at once. If plating for both bacteria and fungi, you may want a minimum of 250 μL of each dilution.

Bacterial isolation

- 9 Use four LB + 100 $\mu\text{g}/\text{mL}$ cycloheximide agar plates to plate 100 μL each of the 10^{-5} , 10^{-6} , 10^{-7} , 10^{-8} dilutions from step 8.

Note

When working with new communities, it may take some trial and error to identify the correct dilution to plate for getting single colonies. These recommendations are based on our past experience.

Cycloheximide is an antifungal. It is not always effective against *Galactomyces*. If trying to remove *Galactomyces* fungi, you can add 21.6 µg/mL natamycin.

- 10 Leave the plates at room temperature until you can see individual bacterial colonies on at least one of the dilution plates.

Note

In our experience with cheese communities, this may take around 3–5 days. Over time, pigments will start to develop in some colonies, which will make identification of different morphologies easier. Leaving the plates exposed to light (e.g. on the bench top) can help encourage pigment production.



Bacterial colonies from a cheese rind community

- 11 Using a sterile toothpick or other sterile instrument, gently touch a single individual colony and quadrant-streak onto a fresh LB + 100 µg/mL cycloheximide agar plate.

- 12 Leave at room temperature until single colonies appear.

Note

You should store plates inside a plastic bag with a moistened paper towel to prevent them from drying out.

- 13 Using a sterile toothpick or other sterile instrument, gently touch a single pure individual colony and inoculate into 2 mL LB broth.

- 14 Shake the culture at 200 rpm at room temperature until cloudy.

- 15 Add 250 μL of culture to 750 μL of sterile 80% glycerol in a labeled 2 mL cryotube, invert to mix well, and immediately store tube at $-80\text{ }^{\circ}\text{C}$.

- 16 Repeat this process until all desired colonies from the original community plate have been stocked.

Note

We usually try to get at least one representative from each unique morphology, but this will depend on your downstream purpose.

Fungal isolation

- 17 Use four PCAMS + 50 $\mu\text{g}/\text{mL}$ chloramphenicol agar plates to plate 100 μL each of the 10^{-1} , 10^{-2} , 10^{-3} , 10^{-4} dilutions from step 8.

Note

When working with new communities, it may take some trial and error to identify the correct dilution to plate for getting single colonies. These recommendations are based on our past experience.

Chloramphenicol is an antibacterial drug.

- 18 Leave the plates at room temperature until you can see individual fungal colonies on at least one of the dilution plates.

Note

In our experience with cheese communities, this may take around 3–7 days for fungi. Morphologies may become more distinguishable over time as more color develops.



Fungal colonies from a cheese rind community

- 19 Using a sterile toothpick or other sterile instrument, gently touch a single individual colony and quadrant streak onto a fresh PCAMS + 50 µg/mL chloramphenicol agar plate.
- 20 Repeat as many times as needed onto a new plate to get a pure single colony.
- 21 From a pure single colony, use a sterile toothpick to pick up fungal spores and then streak the spores in horizontal lines back and forth across a new full plate. Then turn the plate 90° and streak in horizontal lines back and forth across the plate. The goal is to spread the fungus evenly over the surface of the plate.

Note

You should store plates inside a plastic bag with a moistened paper towel to prevent them from drying out.

- 22 Once the fungus is grown and sporulated (~5–7 days), add 1 mL of sterile PBS with 20% glycerol to the plate. Use a cell scraper to gently scrape the surface of the plate to



collect the fungal spores into the liquid.

- 23 Use a 1 mL pipette to collect the liquid with the resuspended spores. Vortex to mix thoroughly.
- 24 Transfer the liquid to a labeled 2 mL cryotube and store at -80°C .
- 25 Repeat this process until all desired colonies from the original community plate have been stocked.

Note

We usually try to get at least one representative from each unique morphology, but this will depend on your downstream purpose.