



May 03, 2022

Version 3

JAX - EZ Lysis Nuclei Isolation for 10x Genomics Assays V.3

DOI

<https://dx.doi.org/10.17504/protocols.io.ewov1n9npgr2/v3>

Shannon Bessonett¹, Gregory A Perry¹, Paul.gabriel¹, Diane Luo¹, Jessica Grassmann¹, William Flynn¹, Elise Courtois¹, Paul Robson², Sandra L Daigle²

¹The Jackson Laboratory; ²The Jackson Laboratory for Genomic Medicine

JAX Single Cell Biology

Tech. support email: scblservice@jax.org



William F F Flynn

The Jackson Laboratory

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov1n9npgr2/v3>

Protocol Citation: Shannon Bessonett, Gregory A Perry, Paul.gabriel, Diane Luo, Jessica Grassmann, William Flynn, Elise Courtois, Paul Robson, Sandra L Daigle 2022. JAX - EZ Lysis Nuclei Isolation for 10x Genomics Assays. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.ewov1n9npgr2/v3> Version created by [William F F Flynn](#)

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 03, 2022

Last Modified: May 03, 2022

Protocol Integer ID: 61887

Keywords: nuclei isolation, 10x genomics, multiome, ez lysis nuclei isolation for 10x genomics assay, 10x genomics multiome assay, protocol for single cell multiome atac, single cell multiome atac, ez lysis nuclei isolation, 10x genomics assay, sigma aldrich nuclei isolation kit, single nuclei from frozen human tissue, gene expression sequencing, single nuclei, human glioblastoma, human placenta tissue, ez prep protocol, gene expression, placenta tissue, nuclei, jax

Abstract

The purpose of this protocol is to produce single nuclei from frozen human tissues for downstream assaying with the 10x Genomics Multiome assay.

This protocol has been demonstrated using Human placenta tissue as well as human Glioblastoma.

This protocol is modified from Sigma Aldrich Nuclei Isolation Kit: EZ Prep protocol and 10x Chromium Demonstrated protocol for Single Cell Multiome ATAC + Gene Expression Sequencing (CG000365).

Guidelines

1. Assess nuclear membrane integrity (refer to 10x Genomics CG000365), using Brightfield or Trypan blue
2. Assess for debris, if large amount of debris present, filter with smaller pore size to decrease debris.
3. Use either Trypan blue and Countess (for clean average size nuclei preps) or AO/PI and Nexcellom K2 counter (if nuclei prep has debris)
4. If RNA content is of concern Imaging nuclei can be done with fluorescence, and Syto RNA stain

Materials

Nuclei Buffer (20x): 2000153/2000207 (10x Genomics)
 Digitonin: BN2006 (Thermo Fisher)
 Trizma Hydrochloride Solution, pH 7.4: T2194 (Sigma Aldrich)
 Magnesium Chloride Solution: M1028 (Sigma Aldrich)
 Sodium Chloride solution, 5M: 59222C (Sigma Aldrich)
 1M Nonidet P40 substitute: 74385 (Sigma Aldrich)
 Protector RNase inhibitor (DO NOT SUBSTITUTE): 3335399001 (Sigma Aldrich)
 Nuclei Isolation Kit: Nuclei EZ Prep: NUC101-1KT (Sigma Aldrich)
 DTT: 43816-10ML (Sigma Aldrich)
 Tween-20: 1662404 (Bio-Rad)
 PBS: 14190-144 (Gibco)
 BSA: 130-091-376 (Miltentyi)
 Countess Slides: C10228 (Invitrogen)
 Flowmi 40um filter (Bel-Art™ H136800040)
 Bel-art disposable pestle: BAF199230000-100EA (Sigma Aldrich)
 Pluristrainer mini filters: 43-10040-40 (Pluriselect)

Buffers

	A	B	C	D
		Stock	Final	1mL
	EZ lysis	-	-	972ul
	Digitonin	5%	0.01%	2ul
	DTT	1000 mM	1 mM	1 µl
	RNase inhibitor 40 U/µl	40 U/µl	1 U/µl	25ul

Lysis Buffer, Prepare fresh, maintain at 4°C

	A	B	C	D
		Stock	Final	4mL
	Tris-HCl (pH 7.4)	1 M	10 mM	40 µl
	NaCl	5 M	10 mM	8 µl
	MgCl₂	1 M	3 mM	12 µl



	A	B	C	D
	Tween-20	10%	1.00%	400 µl
	BSA	10%	0%	40 µl
	DTT	1000 mM	1 mM	4 µl
	RNase inhibitor 40 U/µl	40 U/µl	1 U/µl	100 µl
	Nuclease-free Water	-	-	3.40 ml

Wash Buffer, Prepare fresh, maintain at 4°C

	A	B	C	D
		Stock	Final	1mL
	Nuclei Buffer* (20X)	20X	1X	50 µl
	DTT	1000 mM	1mM	1 µl
	RNase inhibitor	40 U/µl	1 U/µl	25 µl
	Nuclease-free Water	-	-	924 µl

Nuclei Buffer (1X), Prepare fresh, maintain at 4°C

Before start

Familiarize yourself with 10x Genomics protocol CG000365.



Sample Prep

20m

- 1 Slightly thaw sample out on ice and place onto pre chilled Petri dish on ice.
- 2 Cut tissue into small pieces and add to a microcentrifuge tube.

Cell lysis

17m

- 3 In micro centrifuge tube containing the tissue add  100 μ L lysis buffer of lysis buffer.
- 4 Using a plastic pestle to grind tissue by pushing pestle down and twisting to break up the tissue.
Start timer once lysis buffer is added.
- 5 Add  25 μ L lysis buffer to wash off pestle and place tube  On ice for a total of  00:05:00 .
- 6 After 5 minutes of lysis  On ice ,  500 x g, 4°C, 00:05:00 .
- 7 Remove supernatant and resuspend pellet using wide bore tips in  100 μ L lysis buffer .
- 8 Incubate  On ice for a total of  00:05:00 .

Washing

15m

- 9 Add  500 μ L wash buffer and  500 x g, 4°C, 00:05:00 .
- 10  Repeat for 3 total washes.
- 11 After last wash, filter suspension through 40um filter.



- 12 Resuspend sample in 1x Nuclei buffer (10x provided) containing RNase inhibitor, and DTT.

Load

5m

- 13 Load for transposition and continue with the 10x multiome protocol.