

Sep 27, 2023

KAPP-Sen TMC: Dissociation of Pancreatic Acinar and Ducts (non-recovered)

DOI

<https://dx.doi.org/10.17504/protocols.io.eq2lyjnorlx9/v1>

Juliana Alcoforado Diniz¹, Jessica Garofalo¹, Dylan Baker¹, Paul Robson^{1,2,3}

¹The Jackson Laboratory for Genomic Medicine, Farmington, CT, USA;

²Department of Genetics and Genome Sciences, University of Connecticut School of Medicine, Farmington, CT, USA;

³Institute for Systems Genomics, University of Connecticut, Farmington, CT, USA

Cellular Senescence Net...

KAPP-Sen TMC



Ashley M Raynock

UConn Health, UConn Center on Aging

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.eq2lyjnorlx9/v1>

Protocol Citation: Juliana Alcoforado Diniz, Jessica Garofalo, Dylan Baker, Paul Robson 2023. KAPP-Sen TMC: Dissociation of Pancreatic Acinar and Ducts (non-recovered). **protocols.io** <https://dx.doi.org/10.17504/protocols.io.eq2lyjnorlx9/v1>



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 01, 2023

Last Modified: September 27, 2023

Protocol Integer ID: 85795

Keywords: SenNet KAPP-Sen TMC, dissociation of pancreatic acinar, pancreatic acinar, sen tmc, scrna, samples from brain, dead donor, cold from prodolab, kapp

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **[protocols.io](#)** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **[protocols.io](#)**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

The dispersed samples were shipped cold from **PRODOLABS**. Prior to scRNA-seq, dispersed samples from brain dead donor's pancreatic acinar and ducts were dissociated as follows.



Cell Dissociation with TrypLE

1

Note

Before beginning cell dissociation coat all the materials (pipettes, tubes, etc.) with stopping media to prevent sticking. Stopping Media: DMEM + 10% FBS + 1:100 Glutamate.

- 1.1 Distribute specimens into 50 ml conical tubes, try to make similar cell concentrations across all tubes. *** Receive acinar in ~300 ml, should distribute into 12 tubes with 25 ml each***
- 1.2 Centrifuge acinar samples for 2 min at 230 g and 4°C. Leave duct samples to settle without disturbance.
- 1.3 Wash samples 3X with 25 ml PBS.
- 1.4 Add 10 ml TrypLE E and incubate at 37°C, mixing with a pipette approximately every 5 min.
 1. Acinar: in TrypLE E at 37°C for approximately 30 mins.
 2. Ducts: in TrypLE E at 37°C for approximately 15 mins.
- 1.5 To stop reaction, add 30 ml per tube of stopping media.
- 1.6 Centrifuge acinar and ducts at 1300 rpm, 4°C, 2 mins.
- 1.7 To stop: add 30 ml per tube of stopping media.
- 1.8 Centrifuge acinar and ducts at 1300 rpm, 4C, 2 mins.
- 1.9 Aspirate TrypLE E and wash 1X with stop media. Aspirate then resuspend in stop media to (Acinar) 5 ml or (Ducts) 2 ml.
- 1.10 Transfer cell suspension through 70 µm filter then combine to 2 ml (Acinar) and 1ml (Ducts). Filter again through a 40 µm Flowmi.



1.11 Count cells using AO/PI (acridine orange/propidium iodide) Cell Viability Kit for Luna-FL automated cell counter.

1.12 Proceed to cell fixation.

Fixation of Cells & Nuclei for Chromium Fixed RNA Profiling

2 Cells were fixated prior to scRNAseq according to <https://dx.doi.org/10.17504/protocols.io.x54v9py5zg3e/v1>