

Oct 22, 2019 Version 2

## Lambda Fague Lysis V.2

DOI

[dx.doi.org/10.17504/protocols.io.8j7hurn](https://dx.doi.org/10.17504/protocols.io.8j7hurn)

Laura Sánchez<sup>1</sup>

<sup>1</sup>Universidad Complutense de Madrid, AEGIS - Madrid iGEM 2019

AEGIS - Madrid iGEM 2019



Laura Sánchez

---

OPEN  ACCESS



DOI: [dx.doi.org/10.17504/protocols.io.8j7hurn](https://dx.doi.org/10.17504/protocols.io.8j7hurn)

**Protocol Citation:** Laura Sánchez 2019. Lambda Fague Lysis . protocols.io <https://dx.doi.org/10.17504/protocols.io.8j7hurn>

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** October 22, 2019

**Last Modified:** October 22, 2019

**Protocol Integer ID:** 29023

## Abstract

Lambda phage is a bacteriophage that infect Gam- bacterias, such as E.coli, through the outer membrane protein LamB. We use the Lambda phage to characterize if the AEGIS's display system has been properly expressed.

## Guidelines

Make sure to have all work surface and material clean before the start.



## Safety warnings

- ! The phage Lambda is under the white list of microorganism. However, it extremely infective and could contaminate an entire laboratory if not handle with care.

