

Sep 22, 2019

LB Medium (1 Liter)

DOI

dx.doi.org/10.17504/protocols.io.7jshkne

Alba Balletbó¹

¹Wageningen University

iGEM Wageningen 2019



Alba Balletbó

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.7jshkne

Protocol Citation: Alba Balletbó 2019. LB Medium (1 Liter). **protocols.io** <https://dx.doi.org/10.17504/protocols.io.7jshkne>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 22, 2019

Last Modified: September 22, 2019

Protocol Integer ID: 27986

Guidelines

Mixture must be autoclaved before use.



Materials

MATERIALS



Distilled Water



NaCl Merck MilliporeSigma (Sigma-Aldrich) Catalog #53014



Tryptone Fisher Scientific Catalog #BP1421-500



Yeast Extract Thermo Fisher Catalog #211930



- 1 Add LB medium components in a 1 L autoclavable. LB medium components (1 L):

	NaCl	10 gram
	Trypt one	10 gram
	Yeast Extra ct	5 gram

- 2 Add 1 L of distilled water into the bottle.
- 3 Vigorously shake the bottle.
- 4 Place the bottle into the autoclave and sterilize.