

May 04, 2024

Version 2

LC-MS of Native Nanodiscs V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.e6nvw1k97lmk/v2>

Caroline Brown^{1,2,3}, Snehasish Ghosh^{1,2,3}, Kallol Gupta^{1,2,3}

¹Nanobiology Institute, Yale University, West Haven, CT, USA;

²Department of Cell Biology, Yale University School of Medicine, New Haven, CT, USA;

³Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 20815



Caroline Brown

Yale University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.e6nvw1k97lmk/v2>

Protocol Citation: Caroline Brown, Snehasish Ghosh, Kallol Gupta 2024. LC-MS of Native Nanodiscs. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.e6nvw1k97lmk/v2> Version created by **Caroline Brown**

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: May 03, 2024

Last Modified: May 04, 2024

Protocol Integer ID: 99224

Keywords: ms analysis of native nanodisc, ms of native nanodisc, native nanodisc, biological membrane

Abstract

This protocol provides a step-by-step guide for conducting LC-MS analysis of native nanodiscs extracted from biological membranes.



Liquid Chromatography

- 1 Resuspend peptides in water + 0.1 formic acid varying amount of water so that 500ng of peptide can be loaded in a reasonable volume.
- 2 Chromatography is subsequently conducted using home-packed C18 columns (15cm x 75uM ID) for separation of peptides by hydrophobicity.
- 3 Use a 75 minute separation gradient with a buffer system of water/acetonitrile with 0.1% formic acid beginning at 6% acetonitrile and ending at 100% acetonitrile.

Mass Spectrometry

- 4 Data Dependent Acquisition (DDA) mass spectrometry is conducted using an Orbitrap Eclipse instrument.
- 5 MS1 scans are acquired using the Orbitrap detector at a resolution of 120000 and stored in profile mode using a mass range of 350-1400 m/z and an intensity threshold of 1.0e3.
- 6 Use a cycle time of 1 second for ion isolation and MS2 fragmentation isolating ions with a 1.4 m/z window and fragmented with CID energy at 30% activation.
- 7 Product ions are detected at the ion trap set to rapid detection mode with precursor ions being dynamically excluded for 20 seconds after one instance of detection.