

Nov 15, 2019

Version 1

## *Lemna minor* (duckweed) sterilization protocol V.1

DOI

[dx.doi.org/10.17504/protocols.io.9dth26n](https://dx.doi.org/10.17504/protocols.io.9dth26n)

Jason Laurich<sup>1</sup>

<sup>1</sup>University of Toronto



Jason Laurich

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.9dth26n>

**Protocol Citation:** Jason Laurich 2019. *Lemna minor* (duckweed) sterilization protocol. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.9dth26n>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**Protocol for sterilization of duckweed.** This protocol works, but may need to be repeated multiple times.

**Created:** November 15, 2019



**Last Modified:** November 15, 2019

**Protocol Integer ID:** 29843

**Keywords:** Sterilization, Duckweed, Plants, sterilization protocol, duckweed, lemna minor

## Troubleshooting

## Duckweed Sterilization : preparation

- 1 In a tea-strainer, strain duckweed plants under running water (doesn't have to be sterile). This step gets rid of some surface algae.
- 2 Move surface-rinsed duckweed into a water bath; remove up to 50 plants from water bath and place in 50 mL Falcon tube approximately half full with autoclaved Phosphate Buffered Saline (PBS).
- 3 Vortex or shake intensely for 5 minutes

## Duckweed Sterilization : sterilization

- 4 All of the subsequent sterilization steps should take place in a safety cabinet.
- 5 Strain plants into a tea-strainer reserved for bleaching. Plants stuck to the side of the Falcon tube can be dislodge with distilled water.
- 6 Place in (1% sodium hypochlorite) bleach for 60 seconds.
- 7 Remove from bleach; rinse tea-strainer thoroughly by moving around in jar of autoclaved distilled water.
- 8 Soak in autoclaved distilled water for 5 minutes or more. Repeat this step 3 times, in a new jar of fresh autoclaved distilled water each time.
- 9 Open tea-strainer; pour autoclaved medium (e.g. Hoagland's, Krajncic medium) over tea-strainer to dislodge plants. This should be done over a clean, autoclaved jar (I use 500 mL Mason jars). Seal with lid and move to environmental chamber.

## Checking duckweed sterility

- 10 Wait at least 3 days to give plants a chance to recover from the bleaching.



- 11 In the fumehood, transfer plants to autoclaved Mason jar approximately half full of your ENRICHED media of choice. I typically add 1 g yeast extract and 10 g sucrose per L to make it enriched.
- 12 In as little as 2-3 days, you will know if your duckweed remains contaminated with bacteria — in this case, the media will become opaque as microorganisms replicate. If the media remains clear, congratulations!!! You have successfully sterilized your duckweed plants.
- 13 If the media is cloudy, wash plants and transfer to Mason jar with autoclaved, unenriched media (in safety cabinet). After plants have recovered (ie. look green and healthy) you can re-sterilize them until all bacteria are removed.