

Jan 09, 2026

Lentivirus production

DOI

<https://dx.doi.org/10.17504/protocols.io.yxmvmmb656g3p/v1>

Akio Mori¹, Robert Edwards¹

¹University of California San Francisco



Akio Mori

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.yxmvmmb656g3p/v1>

Protocol Citation: Akio Mori, Robert Edwards 2026. Lentivirus production. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.yxmvmmb656g3p/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 20, 2025



Last Modified: January 09, 2026

Protocol Integer ID: 233012

Keywords: production of lentivirus, lentivirus production, lentivirus production this protocol, lentivirus, primary neuronal culture, neuronal culture, cell

Abstract

This protocol describes the production of lentivirus in HEK293T cells for the transduction of primary neuronal cultures.

Materials

Cells

- HEK293T cells (early passage) (CRL-3216, ATCC)

Plasmids

- Gene of interest in lentiviral backbone
- Accessory plasmids: pMDLG/pRRE, pIVS-VSVG, pRSV-Rev

Reagents

- FuGENE transfection reagent (E2311, Promega)
- Neurobasal medium (21103-049, Gibco)
- Poly-L-lysine for coating 6-well plates (P2636, Sigma-Aldrich)

Consumables

- 6-well tissue culture plates
- 0.45 μm filter units



Cell seeding

- 1 Coat 6-well plates with poly-L-lysine for 1 h, and wash twice with PBS.
- 2 Seed early-passage (< P10) HEK293T cells onto coated 6-well plates to reach ~70–80% confluency the next day.

Transfection

- 3 Prepare a DNA mixture per well:
 - a. In 200 μ L Opti-MEM, combine:
 - 1.5 μ g plasmid encoding the gene of interest (pFUGW, pJHUG, or pJHUIC)
 - 0.5 μ g each of accessory plasmids (pMDLG/pRRE, pIVS-VSVG, pRSV-Rev)
 - b. Add 12 μ L FuGENE HD transfection reagent to the DNA solution.
 - c. Mix gently by vortexing.
 - d. Incubate at room temperature for 10–15 min to allow complex formation.
- 4 Remove HEK cells' medium and replace it with 1.8 mL of fresh medium per well.
- 5 Add the transfection mixture to each well and incubate overnight at 37°C, 5% CO₂.

Virus production

- 6 The next morning, remove the transfection medium and replace it with 2 mL neurobasal medium or HEK cells' medium per well.
- 7 Incubate cells for 36–48 h to allow virus production.
- 8 Collect the culture medium containing lentivirus. Filter the collected medium through a 0.45 μ m filter to remove cellular debris. Use the viral supernatant immediately for neuronal transduction, or aliquot and store at -80°C.

Infection of primary neurons

- 9 Add 20–40 μ L viral supernatant per cloning ring of primary neuronal culture to achieve efficient transduction.

