

Jul 17, 2020

Version 2

Lipoxxygenase activity determination V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.bhqmj5u6>



Neillier Junior¹

¹University of Manitoba



Neillier Junior

University of Manitoba, Universidade Federal de Viçosa

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bhqmj5u6>

Protocol Citation: Neillier Junior 2020. Lipoxxygenase activity determination. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.bhqmj5u6>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 20, 2020

Last Modified: July 17, 2020

Protocol Integer ID: 38381

Keywords: lipoxygenase activity determination lipoxygenase activity on linoleic acid, lipoxygenase activity determination lipoxygenase activity, linoleic acid, formed hydroperoxide

Abstract

Lipoxygenase activity on linoleic acid was determined according to the method described by Axelrod et al. (1981). This method determines the increase in absorbance at 234 nm, resulting from the formation of a conjugated double bond system in the formed hydroperoxide.

Materials

MATERIALS

☒ Sodium phosphate monobasic monohydrate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #S9638**

☒ Sodium phosphate dibasic **Merck MilliporeSigma (Sigma-Aldrich) Catalog #S3264**

☒ Tween 20 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P1379-500ml**

☒ Sodium hydroxide **Merck MilliporeSigma (Sigma-Aldrich) Catalog #S8045**

☒ Linoleic acid **Catalog #L1012**

Safety warnings

! Wear personal protective equipment: gloves, lab coat and mask.

Before start

Organize your workspace

Make sure all solutions and equipment are available.

Reagent Preparation

1 10 mM sodium linoleate stock solution



In a 150 mL Erlenmeyer add:

10 mL distilled water (previously boiled)

78 μ L linoleic acid

90 μ L of tween 20

Keep the solution protected from light by wrapping the Erlenmeyer in aluminum foil.

Mix the solution with the aid of a pipette, taking care not to form bubbles.

Add 0.5 M NaOH until the solution is clarified (approximately 100 μ L).

Transfer the solution to a 25 mL volumetric flask protected from light. Make up the volume to 25 mL.

Divide the stock solution of sodium linoleate into 1.5 mL amber microtubes and store at -20°C .

2 **50.0 mM phosphate buffer, pH 6.0**

Mix sodium phosphate monobasic and dibasic solutions in the proportions indicated below, and dilute to 200 mL with deionized water.

- 6.15 mL of 0.2 M sodium phosphate, dibasic dihydrate ($\text{Na}_2\text{HPO}_4 \cdot 2\text{H}_2\text{O}$ FW = 178.05)
- 43.85 mL of 0.2M sodium phosphate, monobasic, monohydrate ($\text{NaH}_2\text{PO}_4 \cdot \text{H}_2\text{O}$ FW = 138.01)

Note: The dibasic stock sodium phosphate may be somewhat harder to dissolve; adding a little heat may help.

Complete the final volume to 200 mL with deionized water.



Adjust the final pH to 6.0.

Procedure

3 Pipette (in microliters) the following reagents into 1.5 mL microtubes

		Blank	Test
	Phosphate Buffer	1002 μL	1000 μL
	Sodium Linoleate Stock Solution	10.0 μL	10.0 μL
	Enzymatic Extract (sample)	-	2.0 μL

Mix by inversion

4 Zero the spectrophotometer with Blank content at A234 nm.

5 Immediately after enzyme addition (Test), mark the time and pour the contents into a suitable cuvette.

After 30 s of reaction onset, monitor readings at A234 nm for 120 s.

Calculations

6 Calculate velocity from absorbance values obtained

$$V_0 = \Delta A_{234 \text{ nm}} (\epsilon l \Delta t)^{-1}$$



- V_0 : enzymatic activity
- $\Delta A_{234\text{ nm}}$: absorbance variation at 234 nm
- ϵ : molar extinction coefficient of linoleic acid hydroperoxides at 234 nm
- l : optical path
- Δt : time (120 s)