



Feb 01, 2024

Live-cell imaging; cell death assay

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzy1d5lx1/v1>

Minee-Liane Choi¹

¹UCL Institute of Neurology



Minee L Choi

Korea Advanced Institute of Science & Technology, The Franci...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzy1d5lx1/v1>

Protocol Citation: Minee-Liane Choi 2024. Live-cell imaging; cell death assay. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.rm7vzy1d5lx1/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 26, 2022

Last Modified: May 31, 2024

Protocol Integer ID: 63250

Keywords: ASAPCRN, examining cell death, cell death, cell imaging, cell, assay

Abstract

This protocol contains instructions for examining cell death in live-cell imaging.

- 1 Cell death is detected using Propidium iodide (PI, Thermo Fisher Scientific) or SYTOX™ Green (SYTOX, Thermo Fisher Scientific) which is excluded from viable cells but exhibits red (PI) and green (SYTOX) fluorescence following a loss of membrane integrity and Hoechst 33342 (Hoechst, Thermo Fisher Scientific).
 - 2 20 µM PI or 500 nM SYTOX and 10 uM Hoechst are directly added to the dishes.
 - 3 Cells are incubated for 15 min.
 - 4 Live-cell imaging is performed using a confocal microscope.
- Note**
- Live-cell imaging is performed using a confocal microscope (Zeiss LSM 710 or 880 with an integrated META detection system). For confocal microscopes, illumination intensity is limited to 0.1-0.2% of laser output to prevent phototoxicity, and the pinhole is set to allow optical slice at approximately 1-2 µm.
- 5 Hoechst and PI are excited by 405 nm and 543 nm laser lines with the emission between 405 nm to 470 nm and 570 nm to 640nm respectively.
 - 6 SYTOX is excited by a 488 nm laser with emissions between 488 nm and 516 nm.