



Feb 01, 2024

## Live-cell imaging; reactive oxygen species (Superoxide)

DOI

[dx.doi.org/10.17504/protocols.io.4r3l2oyj4v1y/v1](https://dx.doi.org/10.17504/protocols.io.4r3l2oyj4v1y/v1)

Minee L Choi<sup>1</sup>

<sup>1</sup>UCL Institute of Neurology



Minee L Choi

Korea Advanced Institute of Science & Technology, The Franci...

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l2oyj4v1y/v1>

**Protocol Citation:** Minee L Choi 2024. Live-cell imaging; reactive oxygen species (Superoxide). **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.4r3l2oyj4v1y/v1>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

We use this protocol and it's working

**Created:** May 25, 2022



**Last Modified:** May 31, 2024

**Protocol Integer ID:** 63243

**Keywords:** ASAPCRN, measuring superoxide, rate of cytosolic superoxide generation, cytosolic superoxide generation, superoxide generation, rate of superoxide generation, cell imaging, superoxide, reactive oxygen species, oxidized dye, cell, reduced dye, oxidized form, form of the dye, dye

## Abstract

This protocol contains the instruction for measuring superoxide using dihydroethidium (HEt) which allows the rate of superoxide generation to be measured which is present as the ratio of the oxidized form of the dye over the reduced form. HEt allows the rate of cytosolic superoxide generation to be measured, by the slope of the ratio of oxidized dye to reduced dye.

## Troubleshooting

- 1 Cells are washed and loaded [M] 2 micromolar ( $\mu\text{M}$ ) dihydroethidium (HEt; Thermo Fisher Scientific) in the recording buffer.

#### Note

HEt is an indicator of superoxide which exhibits blue fluorescence in the cytosol before oxidation, and the nucleus presents a red fluorescence upon oxidation.

- 2 The recording is performed using an epi-fluorescence inverted microscope equipped with 20x objective after a quick loading (2-3 min) in order to limit the intracellular accumulation of oxidized product, and the dye was present throughout the imaging.

#### Note

Live-cell imaging is performed using an epi-fluorescence inverted microscope equipped with a CCD camera (Retiga; QImaging). For epi-fluorescence inverted microscope, excitation is provided by a xenon arc lamp with the beam passing through a monochromator (Cairn Research) and emission was reflected through a long-pass filter to a cooled CCD camera and digitized to 12-bit resolution (Digital Pixel Ltd, UK).

- 3 Excitation is at 530 nm, and emission recorded above 560 nm is assigned to the oxidized form, while excitation at 380 nm and emission collected from 405 nm to 470 nm is assigned to the reduced form.
- 4 The ratio of the fluorescence intensity, resulting from its oxidized/reduced forms, is quantified.