

Apr 18, 2024

Liver Tissue RNA Isolation - University of Minnesota TMCs

DOI

<https://dx.doi.org/10.17504/protocols.io.n92ld83yxv5b/v1>

Laura J Niedernhofer¹, David A Bernlohr¹

¹University of Minnesota, Minneapolis, MN

Cellular Senescence Net...



Allie Abdellatif

MIBAM, University of Minnesota, Minneapolis, MN

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ld83yxv5b/v1>

Protocol Citation: Laura J Niedernhofer, David A Bernlohr 2024. Liver Tissue RNA Isolation - University of Minnesota TMCs. protocols.io <https://dx.doi.org/10.17504/protocols.io.n92ld83yxv5b/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 26, 2024

Last Modified: April 18, 2024

Protocol Integer ID: 97411

Keywords: liver tissue rna isolation, minnesota tmcs invitrogentm trizoltm plus rna purification kit, rna from liver tissue, rna isolation, total rna, rna, minnesota genomics center, liver tissue, purification

Funders Acknowledgements:

NIH

Grant ID: 5U54AG079754-02

NIH

Grant ID: 5U54AG076041-03

Abstract

Invitrogen™ TRIzol™ Plus RNA Purification Kit provides a simple, reliable, and rapid method for isolating high-quality total RNA from a wide variety of samples, including animal and plant cells and tissue, bacteria, and yeast.

This protocol was used to isolate RNA from liver tissue for BulkRNA sequencing in collaboration with The University of Minnesota Genomics Center.



TFS-Assets-LSG-manuals-Trizol_Pl... 196.1KB



Isolate RNA

- 1
 - a. Lyse and homogenize samples in TRIzol Reagent by adding 1mL TRIzol reagent per 50-100mg tissue to the sample and use a homogenizer
 - b. Incubate for 5 minutes to permit complete dissociate of the nucleoproteins complex
 - c. Add 0.2mL of chloroform per 1mL TRIzol Reagent used for lysis, then securely cap the tube.
 - d. Incubate for 2-3 minutes.
 - e. Centrifuge the sample for 15 minutes at  12000 x g at  4 °C - the mixture separates into a lower red phenol-chloroform, interphase and a colorless upper aqueous phase.
 - f. Transfer ~600 µL of the colorless upper phase containing the RNA to a new tube
 - g. Add an equal volume of 70% ethanol, then mix well by vortexing.
 - h. Invert the tube to disperse any visible precipitate that may form after adding ethanol.

Bind the RNA to the membrane

- 2
 - a. Transfer up to 700 µL of the sample to a spin cartridge with (collection tube).
 - b. Centrifuge  12000 x g for 15 seconds.
 - c. Discard the flow through, then reinsert the spin cartridge into the same collection tube.
 - d. Repeat step 2a -step 2c until the entire sample has been processed.

Wash the RNA on the membrane

- 3
 - a. Add 700 µL of Wash Buffer I to the spin cartridge.
 - b. Centrifuge at  12000 x g for 15 seconds.
 - c. Discard the flow-through, then reinsert the spin cartridge into the same collection tube.
 - d. Add 500 µL of Wash Buffer II to the spin cartridge.
 - e. Centrifuge at  12000 x g for 15 seconds.
 - f. Discard the flow-through, then reinsert the spin cartridge into the same collection tube.
 - g. Repeat step 3d-3f once.

Elute the RNA

- 4
 - a. Centrifuge at  12000 x g for 1 minute to dry the membrane.
 - b. Discard the collection tube, then insert the spin cartridge into a recovery tube.



- c. Add 30 μL -3 $\times$ 100 μL (3 sequential elutions with 100 μL each) of RNase-free water to the center of the spin cartridge.
- Note: If you are performing sequential elutions, collect all eluates in the same tube.
- d. Incubate for 1 minute.
- e. Centrifuge at $>12,000 \times g$ for 2 minutes.
- f. Discard the spin cartridge.

The recovery tube contains the purified total RNA.

- Store the purified RNA on ice if used within a few hours. For long-term storage, store the purified RNA at -80°C .
- If highly pure RNA without genomic DNA contamination is required, perform DNase I treatment after purification (see PureLinkTM RNA Mini Kit User Guide (Pub. No. MAN0000406)).