

Jan 24, 2019

Lysogeny Broth (1L)

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.xf5fjq6>

Brian Smith¹, baltrus@email.arizona.edu micro.tucson¹

¹University of Arizona



Brian A Smith

University of Arizona

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.xf5fjq6>

Protocol Citation: Brian Smith, baltrus@email.arizona.edu micro.tucson 2019. Lysogeny Broth (1L). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.xf5fjq6>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use these protocols commonly in our group and they are working.

Created: January 24, 2019

Last Modified: January 24, 2019

Protocol Integer ID: 19677

Keywords: lysogeny broth, 1l



Fill container aproximately 750mL of H2O

- 1 Add stir bar and set to mix

Add ingredients

- 2
 -  10 g Bacto Tryptone
 -  5 g Bacto Yeast Extract
 -  10 g NaCl

Bring volume to 1L by adding H2O

- 3 Allow to mix again with stir bar.

Autclave

- 4 Add agar to container to be autclaved
 -  15 gThen add media
- 5 Autclave at 121°C for 30min