

Oct 20, 2023

Magnetic resonance imaging (MRI)

 [Nature Communications](#)

 In 3 collections

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l6x3mdlqe/v1>

Núria Peñuelas¹, Ariadna Laguna¹, Miquel Vila¹

¹Vall d'Hebron Research Institute

Vilalab Public

Nuria



Miquel Vila

VHIR-CIBERNED-ASAP

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l6x3mdlqe/v1>

External link: <https://doi.org/10.1038/s41467-024-53168-7>

Protocol Citation: Núria Peñuelas, Ariadna Laguna, Miquel Vila 2023. Magnetic resonance imaging (MRI). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bp2l6x3mdlqe/v1>

**Manuscript citation:**

Laguna, A., Peñuelas, N., Gonzalez-Sepulveda, M. *et al.* Modelling human neuronal catecholaminergic pigmentation in rodents recapitulates age-related neurodegenerative deficits. *Nat Commun* **15**, 8819 (2024). <https://doi.org/10.1038/s41467-024-53168-7>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 19, 2023

Last Modified: November 17, 2023

Protocol Integer ID: 89600

Keywords: magnetic resonance imaging, mouse brain, mri, brain, mouse

Abstract

Magnetic resonance imaging (MRI) of mouse brain

- 1 Prepare the equipment: 7T Bruker BioSpec 70/30USR scanner (Bruker BioSpin GmbH, Ettlingen, Germany) equipped with a mini-imaging gradient set (400mT/m) and using a 72-mm inner diameter linear volume coil as a transmitter and a dedicated mouse brain surface coil as a receiver.
- 2 Prepare the software: Acquire and process all MRI data on a Linux computer using Paravision 5.1 software (Bruker BioSpin GmbH, Karlsruhe, Germany) (RRID:SCR_001964, <http://www.bruker.com/service/support-upgrades/software-downloads/mri.html>).
- 3 Anesthetize the animals (1.5-2% isoflurane in 1 L/min oxygen for maintenance) and place them into an animal bed with bite-bar and ear-bars for optimal head immobilization.
- 4 Control the respiration rate (50-100 bpm) and core body temperature ($37 \pm 1^{\circ}\text{C}$) with an animal monitoring and control system (SA Instruments, Stony Brook, NT).
- 5 Acquire high resolution T1-weighted spin-echo images in the axial plane with and without fat saturation pulses.
- 6 Acquire two slice packages of 7 contiguous slices each containing the SNpc and LC regions and using the following parameter: TE=7 ms; TR= 500 ms; NA=56; FOV=1.92×1.92 cm²; MTX=128×128; ST= 0.25 mm; thus, resulting in a spatial resolution of 150×150×250 μm³; Acqt= 1h per image.