

May 23, 2025

Maintenance of hPSC

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l6xmjrlqe/v1>



Lyn Healy¹, Valeria Fernandez Vallone², Nathalie Lefort³, Katarzyna Ludwik², Tamer Onder⁴, Lisa Pavinato^{5,6}, Fatma Visal OKUR⁷, Harald Stachelscheid²

¹The Francis Crick Institute;

²Core Unit pluripotent Stem Cells and Organoids - Berlin Institute of Health @ Charite, Berlin, Germany;

³Université de Paris, Imagine Institute, iPSC Core Facility, INSERM UMR U1163, F-75015 Paris, France.;

⁴Koc University;

⁵Institute of Oncology Research (IOR), Bellinzona Institutes of Science (BIOS+), Bellinzona, Switzerland;

⁶Faculty of Biomedical Sciences, Università della Svizzera Italiana, Lugano, Switzerland.;

⁷Hacettepe University, Center for Stem Cell Research and Development (PEDISTEM) and Hacettepe University Faculty of Medicine, Department of Pediatrics

CorEuStem



Harald Stachelscheid

Core Unit pluripotent Stem Cells and Organoids - Berlin Inst...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l6xmjrlqe/v1>

Protocol Citation: Lyn Healy, Valeria Fernandez Vallone, Nathalie Lefort, Katarzyna Ludwik, Tamer Onder, Lisa Pavinato, Fatma Visal OKUR, Harald Stachelscheid 2025. Maintenance of hPSC. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.bp2l6xmjrlqe/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 15, 2023

Last Modified: May 23, 2025

Protocol Integer ID: 90972

Keywords: iPSC, hPSC, confluence, maintenance, spontaneous differentiation, maintenance of established hpSC line, maintenance of hpSC, established hpSC line, coated culture vessel, culture vessel, hpSC, maintenance

Funders Acknowledgements:

COST CorEuStem

Grant ID: CA20140

Disclaimer

The reported protocols are based on the authors experience, and may partially differ from the original protocols provided by the companies.

Abstract

This protocol describes maintenance of established hPSC lines in expansion media on matrix coated culture vessels.

Guidelines

Depending on the growth characteristics of specific hPSC line, cells must be passaged every 5-7 days in order to maintain log phase of growth and avoid induction of differentiation.

In general hPSC lines are maintained in a 6 well format, however information on other tissue culture vessel formats are also included in this protocol as a guidance.



Materials

LABORATORY EQUIPMENT AND CONSUMABLES

Use sterile material

- 1/5/10 mL serological pipettes
- 15/50 mL conical tubes
- Cell culture treated plastic vessels of choice e.g. 24, 12 or 6-well plates, T25, T75 flasks, 10cm dishes
- 10/200/1000µL tips and micropipettes (optional)
- Aspirator pump with disposable pipette
- Centrifuge
- Microscope, if available Stereo Microscope
- Disposable scraping tools (pipette tips, glass rods, small size cell scraper, syringe and needle)
- Incubator at 37°C and 5% CO₂ or under hypoxic conditions, 5% O₂/ 5% CO₂
- Class II Biosafety Cabinet

MEDIA AND REAGENTS

Vitronectin

⊗ Vitronectin (VTN-N) Recombinant Human Protein, Truncated **Thermo Fisher Catalog #A14700**

⊗ Vitronectin XF™ **STEMCELL Technologies Inc. Catalog ##07180**

Laminin

⊗ MACSmatrix Laminin 511 **Miltenyi Biotec Catalog #130-136-454**

⊗ rhLaminin-521 **Thermo Fisher Catalog #A29248**

⊗ CellAdhere™ Laminin-521 100 µg **STEMCELL Technologies Inc. Catalog #77003**

Basement Membrane Products

⊗ Geltrex®; LDEV-Free, hESC-Qualified, Reduced Growth Factor Basement Membrane Matrix **Thermo Fisher Catalog #A1413301**

⊗ Matrigel hESC-qualified (Corning Cat# 354277) **Corning Catalog #354277**

Cell Culture Media

⊗ Essential 8™ Medium **Gibco - Thermo Fisher Scientific Catalog #A1517001**

⊗ Essential 8 Flex complete medium (E8) **Gibco - Thermo Fisher Scientific Catalog #A2858501**



- ⊗ StemFlex Medium Thermo Fisher Scientific Catalog #A3349401
- ⊗ mTeSR™1 500 mL Kit STEMCELL Technologies Inc. Catalog #85850
- ⊗ mTeSR plus media kit STEMCELL Technologies Inc. Catalog #100-0276
- ⊗ eTeSR STEMCELL Technologies Inc. Catalog #100-1215
- ⊗ StemMACS™ iPS-Brew XF human: Basal medium and supplement Miltenyi Biotec Catalog #130-104-368

Homemade Cell Culture Media: Essential 8 (E8) Medium (see protocol [E8 media production](#))

Selection of culture conditions

15m

1 The following points should be considered when choosing hPSC culture conditions:

15m

- A number of media will work with an array of different matrices.
- The rate of growth of cell lines will differ between different matrix/media combinations and cell lines may exhibit a transient change in morphology until adapted to the new culture conditions.
- When transitioning cell lines between different media/matrix combinations follow the manufacturers guidance. Usual recommendations: do not switch the media and matrix at the same time. Make a gradual change of media including proportions until full transition.
- hPSC often require a couple of passages to adapt to the new culture conditions.
- Following transition to new culture conditions, perform quality control testing as appropriate for the intended use of the cells.
- For commercial media, prepare, store and use as directed by the manufacturer. For media made in house, use and store under the same conditions established for the validation of the media.

	A	B	C	D	E	F	G
	Media/Matrix	Truncated Vitronectin	Vitronectin XF	Matrigel	Geltrex	Laminin 521	Laminin 511
	Essential E8	XX	X	X	XX	X	
	Essential E8 Flex	XX	X	X	XX	XX	
	mTeSR1	X	XX	XX	X		
	mTeSR Plus		XX	XX			
	StemMACS iPSC-Brew XF			XX	XX	XX	XX
	Stem Flex			X	XX	XX	X

Table 2. Media and matrix possible combinations

XX = recommended by manufacturer and tested by CoreEuStem member

X = tested and recommended by CoreEuStem member

Homemade Cell Culture Media: Essential 8 (E8) Medium (see protocol [E8 media production](#))

**Note**

The matrix media list is not exhaustive. It reflects the media and matrices that are used by the members of CorEuStem for their maintenance protocols.

- 2 Inspect hPSC cultures daily using a microscope to monitor morphology, the presence of spontaneous differentiated cells, and confluence. Based on these observations, determine if the cultures require further action (e.g. removal of differentiated cells, passaging). Use reference images in the protocol: **Reference pictures of hiPSC cultured in defined conditions** to guide your assessment of the cultures. Cell lines may vary in their growth rate and morphology in different media/matrix combinations.

15m

STEP CASE**Culture >70% confluent** 1 step

Culture is >70% confluent and < 10% of the colonies in the culture appear differentiated upon visual inspection.

- 3 If the cells require passaging, refer to the protocol **Non-enzymatic passaging of hPSC** or **Single cell passaging of hPSC**. Otherwise perform a complete media change (refer to Culture <70% confluent tab).

Note

Assessing confluency is inherently subjective. Whenever possible, use imaging systems capable of calculating the surface area of the culture vessel occupied by cells to ensure a consistent and objective evaluation of cell culture confluency.

If such imaging tools are not available, utilize a collection of images captured during the culture process at various stages of confluency to train staff in recognizing and estimating different confluency percentages.

Keep in mind that the growth rate required to achieve confluency is influenced by both the specific cell line and the cell culture system used.

Protocol references

- Ludwig T, Bergendahl V, Levenstein M, et al. Feeder-independent culture of human embryonic stem cells. *Nat Methods*. 2006;3(8):637–646. <https://doi.org/10.1038/nmeth902>
- Chen G, Gulbranson D, Hou Z, et al. Chemically defined conditions for human iPSC derivation and culture. *Nat Methods*. 2011;8(5):424–429. <https://doi.org/10.1038/nmeth.1593>
- Passaniti A, Kleinman HK, Martin GR. Matrigel: history/background, uses, and future applications. *J Cell Commun Signal*. 2022;16(4):621–626. <https://doi.org/10.1007/s12079-021-00643-1>
- Braam SR, Zeinstra L, Litjens S, et al. Recombinant vitronectin is a functionally defined substrate that supports human embryonic stem cell self-renewal via alphavbeta5 integrin. *Stem Cells*. 2008;26(9):2257–2265. <https://doi.org/10.1634/stemcells.2008-0291>
- Rodin S, Antonsson L, Niaudet C, et al. Clonal culturing of human embryonic stem cells on laminin-521/E-cadherin matrix in defined and xeno-free environment. *Nat Commun*. 2014;5:3195. <https://doi.org/10.1038/ncomms4195>
- Nakagawa, M., Taniguchi, Y., Senda, S. *et al*. A novel efficient feeder-free culture system for the derivation of human induced pluripotent stem cells. *Sci Rep* **4**, 3594 (2014). <https://doi.org/10.1038/srep03594>