



Mar 14, 2025

Making Drosophila Artificial Hemolymph

DOI

<https://dx.doi.org/10.17504/protocols.io.q26g7mqg9gwz/v1>

Zachary Freyberg¹

¹University of Pittsburgh



Zachary Freyberg

University of Pittsburgh

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g7mqg9gwz/v1>

Protocol Citation: Zachary Freyberg 2025. Making Drosophila Artificial Hemolymph. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.q26g7mqg9gwz/v1>

Manuscript citation:

Doi: 10.1038/ncomms10652

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 14, 2025

Last Modified: March 14, 2025

Protocol Integer ID: 124407

Keywords: Drosophila, Brain, Dissection, making drosophila artificial hemolymph, drosophila artificial hemolymph, artificial hemolymph, making artificial hemolymph, drosophila brain dissection, imaging, brain,

Abstract

This is a protocol for making artificial hemolymph (AHL) for Drosophila brain dissection and imaging.

Materials

NaCl
KCl
CaCl₂
MgCl₂
NaH₂PO₄
Sucrose
Trehalose
HEPES
ddH₂O
pH meter

Safety warnings

 N/A

Ethics statement

N/A

Before start

Make sure to make the stock solutions ahead of making the final AHL buffer.



Making Drosophila Artificial Hemolymph (AHL)

1 Make stock solutions

1.1 Stock Solutions

3M NaCl	MW: 58.44	87.66g to 500mL
1M KCl	MW: 74.56	37.28g to 500mL
500mM CaCl ₂	MW: 147.00	14.7g to 200mL
500mM MgCl ₂	MW: 203.3	20.33g to 200mL
500mM NaH ₂ PO ₄	MW: 155.99	15.59g to 200mL
*200mM Sucrose	MW: 342.30	34.23g to 500mL
*100mM Trehalose	MW 378.33	3.78g to 100mL
*250mM HEPES	MW: 238.31	11.92 g to 200mL
*Store at 4°		

2 Make 100 mL of AHL from the stock solutions

2.1	108mM NaCl	3.6mL of 3M NaCl
	5mM KCl	.5mL of 1M KCl
	2mM CaCl ₂	.4mL of 500mM CaCl ₂
	8.2mM MgCl ₂	1.64mL of 500mM MgCl ₂
	1mM NaH ₂ PO ₄	.2ml of 500mM NaH ₂ PO ₄
	10mM Sucrose	5mL of 200mM Sucrose
	5mM Trehalose	5mL of 100mM Trehalose
	5mM HEPES	2mL of 250mM HEPES
	4mM NaHCO ₃	.4mL of 1M NaHCO ₃

**add NaHCO₃ last to avoid formation of precipitates

2.2 Add ddH₂O to a final volume of 100 mL

3 Adjust final pH to 7.5

4 Check osmolarity of the solution to ensure that it is at ~265 mOsm



Protocol references

Freyberg, Z, Sonders, MS, Aguilar, JI, Hiranita, T, Karam, CS, Flores, J, Pizzo, AB, Zhang, Y, Farino, ZJ, Chen, A, Martin, CA, Kopajtic, TA, Fei, H, Hu, G, Lin, YY, Mosharov, EV, McCabe, BD, Freyberg, R, Wimalasena, K, Hsin, LW, Sames, D, Krantz, DE, Katz, JL, Sulzer, D & Javitch, JA 2016, 'Mechanisms of amphetamine action illuminated through optical monitoring of dopamine synaptic vesicles in Drosophila brain', *Nature communications*, vol. 7, 10652. <https://doi.org/10.1038/ncomms10652>