

Apr 05, 2023

Matrix Sublimation via In-House Developed Sublimation Apparatus

DOI

<https://dx.doi.org/10.17504/protocols.io.n92ldpyn9l5b/v1>

David Anderson¹, Jeffrey D. Messinger², Jamie Allen¹, Katerina Djambazova¹, Angela Kruse¹, Christine A. Curcio², Kevin L. Schey¹, Richard M. Caprioli¹, Jeff Spraggins¹, Martin Dufrense¹

¹Vanderbilt University; ²University of Alabama at Birmingham

VU Biomolecular Multim...

Human BioMolecular Atl...



Katerina V Djambazova

Vanderbilt University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ldpyn9l5b/v1>

Protocol Citation: David Anderson, Jeffrey D. Messinger, Jamie Allen, Katerina Djambazova, Angela Kruse, Christine A. Curcio, Kevin L. Schey, Richard M. Caprioli, Jeff Spraggins, Martin Dufrense 2023. Matrix Sublimation via In-House Developed Sublimation Apparatus. [protocols.io https://dx.doi.org/10.17504/protocols.io.n92ldpyn9l5b/v1](https://dx.doi.org/10.17504/protocols.io.n92ldpyn9l5b/v1)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: September 28, 2022

Last Modified: October 18, 2023

Protocol Integer ID: 70635

Keywords: MALDI IMS, matrix, Sublimation, high resolution maldi imaging mass spectrometry, matrix sublimation, matrix sublimation via in, house developed sublimation apparatus scope, sublimation device, developed sublimation device, small molecule matrices onto tissue section, small molecule matrice, scope, concentration, matrix type

Abstract

Scope:

This protocol describes the use of an in-house developed sublimation device to deposit small molecule matrices onto tissue sections for high resolution MALDI imaging mass spectrometry (IMS). Concentrations will vary based on matrix type.

Materials

Slides with sample affixed

Matrix of choice

Solvents: acetone, isopropyl alcohol, ethanol

Dry ice



Sample Section Preparation

- 1 Chill three solutions of [M] 150 millimolar (mM) ammonium formate On ice for 30 minutes.
- 2 Wash sectioned tissue slides with three separate solutions of [M] 150 millimolar (mM) ammonium formate for 00:00:45 each. 45s
- 3 Dry tissue using a stream of nitrogen gas.
- 4 Store in a vacuum desiccator for 00:20:00 . 20m


Cold Trap Preparation

- 5 Ensure the cold sleeve is at least 2/3 full of ethanol or isopropyl alcohol when trap is fully seated.
- 6 Carefully add dry ice chips to the ethanol bath in the sleeve and allow chips to sublime off.
Repeat until minimal sublimation of dry ice is observed (ethanol has reached equilibrium temperature).



Sublimation Apparatus Set up

- 7 Ensure that the vacuum line is securely installed between the cold trap and the apparatus lid.
- 8 Ensure that the electrical connections between the heater controller and the passthrough in the apparatus base are secure and remain dry.
- 9 Place sample slides in the sample holder indentations.
Use copper tape or magnetic holder as needed to secure samples as close to the center of the holder as possible.
- 10 Place sample holder in the lid using the embedded magnets.

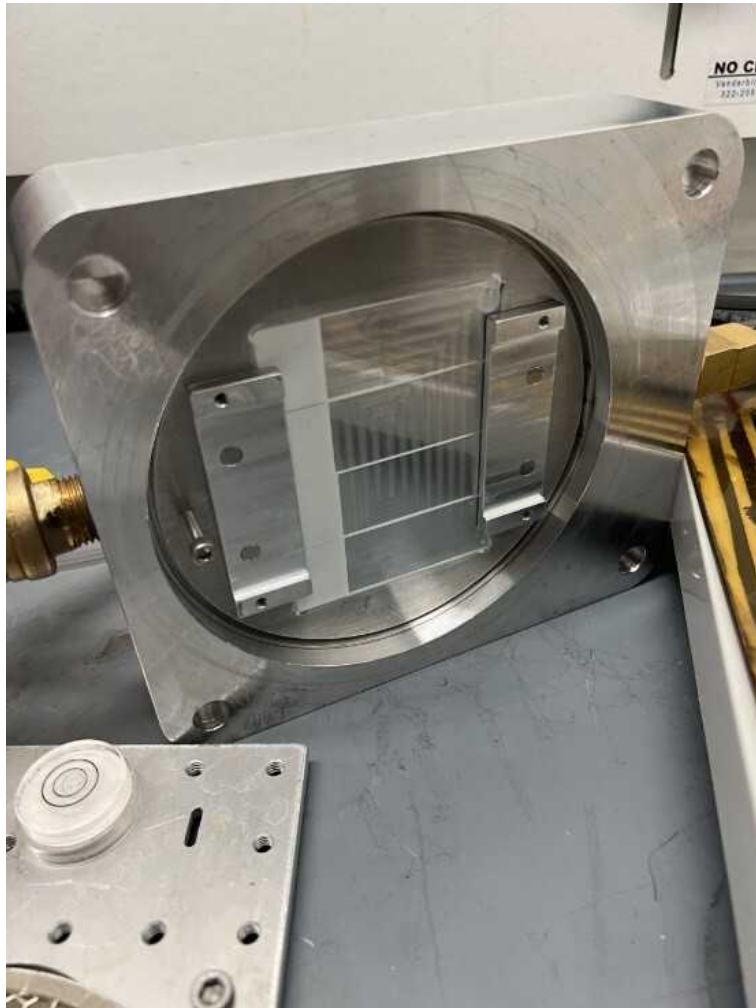

Matrix Preparation

5m

- 11 Dissolve 5mg of matrix in  2 mL of acetone for a total matrix concentration of [M] 2.5 mg/mL .


Note

Concentration will vary depending on matrix used.

- 12 Place matrix pan on a level surface and evenly distribute the matrix solution by dispensing the solution in an equilateral triangle around the center of the matrix pan.

Note

Dispense three solutions of 0.667 mL each.

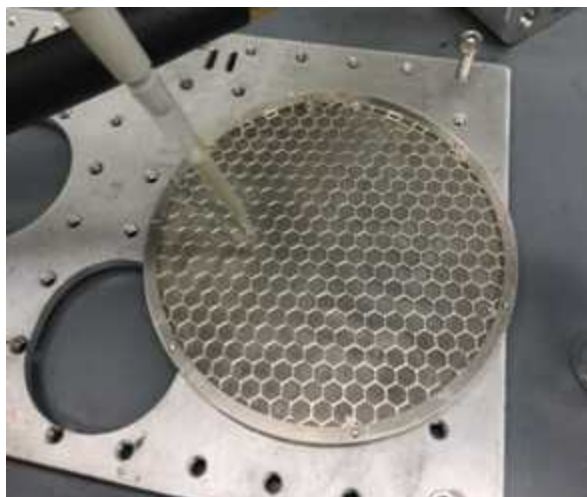

- 13 Wait for solvent to dry (⌚ 00:03:00).

3m

Sublimation Procedure

30m

- 14 Place matrix pan on the heat plate in the apparatus base.
- 15 Check that mating surfaces of the lid, base, and the O-ring are clean, and carefully seat the O-ring in the groove in the base.
- 16 Place apparatus lid on the base, ensure that base is fully seated in the indentions on the induction heater top cover. Start pulling vacuum.
- 17 Once it is clear that vacuum is established, ensure that bath drain stopcock is closed. Add a small amount of acetone (🧴 20 mL) into the well than fill with dry ice and allow to cool for ⌚ 00:03:00 .
- 18 Wait for the pressure to fall to < 50 mTorr measured at the pump (usually ~25 mTorr). Note the pressure at the pump and at the apparatus.

3m



- 19 Turn on internal thermocouple.
Turn on induction heater, set heater to 390F temperature setting, and start a Start
 00:10:00 timer.

10m

Note

Temperature will fluctuate. Make sure ice bath is stocked with dry ice and that there is sufficient melt to cover the bath surface.

- 20 Turn off the heater and remove dry ice chips from the ice bath. Replace ice bath with warm water and open to drain excess water. Repeat warm water exchanges until the apparatus is at room temperature.

Note

A heat gun can be used to help warm up the valve prior to turning it.
The entire process may take up to 15 minutes to complete.

- 21 Turn off the vacuum pump and break vacuum with the leak valve on the vacuum manifold and remove matrix-coated samples.

Clean Up

5m

- 22 Clean matrix pan and sample holder with acetone.

Annealing

1m

- 23 Place samples on a heating plate (130°C) for 30 seconds.