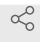


Oct 14, 2019

MC38 and B16-F10 culture media

Elinor Gottschalk¹, Bulent Arman Aksoy¹, Pinar Aksoy¹, Jeff Hammerbacher¹¹Medical University of South Carolina

1 Works for me

 Sharedx.doi.org/10.17504/protocols.io.2y2gfye Elinor Gottschalk

DOI

dx.doi.org/10.17504/protocols.io.2y2gfye

PROTOCOL CITATION

Elinor Gottschalk, Bulent Arman Aksoy, Pinar Aksoy, Jeff Hammerbacher 2019. MC38 and B16-F10 culture media. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.2y2gfye>

LICENSE

————— This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

CREATED

May 17, 2019

LAST MODIFIED

Oct 14, 2019

PROTOCOL INTEGER ID

23290

MATERIALS TEXT

MATERIALS

[EMD Millipore™ Stericup™ Sterile Vacuum Filter Units Fisher](#)

Scientific Catalog #SCGPU05RE

[Fetal Plus® Atlas](#)

Biologicals Catalog # FP-0500-A

[Penicillin-Streptomycin Gibco - Thermo](#)

Fisher Catalog #15140122

[HyClone™ Dulbeccos High Glucose Modified Eagles Medium Fisher](#)

Scientific Catalog #SH30022FS

[Trypsin 0.05% 1X Solution VWR](#)

Scientific Catalog #16777-202

[CytoOne T75 filter cap TC flask USA](#)

Scientific Catalog #CC7682-4875

Mix components and sterile filter

1. Working in the hood, remove 50 ml of DMEM from the 500 ml bottle ([HyClone™ Dulbecco's High Glucose Modified Eagles Medium](#)).
2. To the remaining 450 ml DMEM add: 50 ml FBS ([Fetal Plus](#)) (10%) and 5 ml [penicillin-streptomycin](#) (1%).
3. Sterile filter into a [500 ml 0.22 filter bottle](#).
4. Store at **4 °C** in .