

Jan 11, 2026

Measurement of Lysosomal pH Using LysoSensor Yellow/Blue DND-160

DOI

<https://dx.doi.org/10.17504/protocols.io.q26g7nrx1lwz/v1>

Ali Ghoochani^{1,2,3,4}, Monther Abu-Remaileh^{1,2,3,4}

¹Department of Chemical Engineering, Stanford University, Stanford, CA 94305, USA;

²Department of Genetics, Stanford University, Stanford, CA 94305, USA;

³The Institute for Chemistry, Engineering and Medicine for Human Health (Sarafan ChEM-H), Stanford University, Stanford, CA 94305, USA;

⁴Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 20815, USA

asap



Ali Ghoochani

Stanford University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g7nrx1lwz/v1>

Protocol Citation: Ali Ghoochani, Monther Abu-Remaileh 2026. Measurement of Lysosomal pH Using LysoSensor Yellow/Blue DND-160. protocols.io <https://dx.doi.org/10.17504/protocols.io.q26g7nrx1lwz/v1>



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 08, 2025

Last Modified: January 11, 2026

Protocol Integer ID: 222007

Keywords: measurement of lysosomal ph, quantitative lysosomal ph value, lysosomal ph, ratiometric dye lysosensor yellow, using lysosensor yellow, lysosensor yellow, fluorescence measurement, ph calibration curve, fluorescence, dye incubation

Abstract

This protocol describes the measurement of lysosomal pH using the ratiometric dye LysoSensor Yellow/Blue DND-160. The method includes dye incubation, fluorescence measurement, and the use of a pH calibration curve to derive quantitative lysosomal pH values.

Materials

- LysoSensor Yellow/Blue DND-160 (Thermo Fisher Scientific)
- 1x DPBS
- MES calibration buffers (pH 3.5–7.0)
- eBioscience Monensin Solution (1000X) (Thermo Fisher Scientific, Cat# 00-4505-51)
- Nigericin sodium salt (Sigma, Cat# SML1779)
- Microplate reader capable of dual excitation/emission readings



Materials

1

LysoSensor Yellow/Blue DND-160 (Thermo Fisher Scientific)
1X DPBS
MES calibration buffers (pH 3.5–7.0)
eBioscience Monensin Solution (1000X) (Thermo Fisher Scientific, Cat# 00-4505-51)
Nigericin sodium salt (Sigma, Cat# SML1779)
Microplate reader capable of dual excitation/emission readings

MES Calibration Buffer Composition:

2

- 20 mM MES
- 20 mM NaCl
- 110 mM KCl
- 1x monensin
- 20 μ M nigericin

Procedure:

20m

- 3 Culture SH-SY5Y cells in complete media. Once cells become 90% confluence, wash cells with 1XDPBS and add serum-depleted DMEM for 72 hours.

Note

This is a general protocol for measuring lysosomal pH. It can be applied to a variety of cell types, with or without serum starvation depending on the specific experimental conditions.

- 4 Wash cells once or twice with 1x DPBS.

- 5 Incubate cells (experimental cells and cells for generating pH calibration curve) with 5 μ M LysoSensor Yellow/Blue DND-160 in DMEM for  00:10:00 at  37 °C

10m

- 6 Optional: Wash cells once with 1x DPBS.



- 7 Generate a calibration curve by equilibrating LysoSensor-loaded cells in MES calibration buffers with pH values ranging from 3.5 to 7.0.
- 8 Incubate cells in each buffer for  00:10:00 at  37 °C before measuring fluorescence. 10m
- 9 Measure fluorescence using a microplate reader at the following settings:
 - Excitation at 340 nm → Emission at 440 nm
 - Excitation at 380 nm → Emission at 540 nm
- 10 Calculate the emission ratio (440/540 nm), which correlates with lysosomal pH.
- 11 Use the calibration curve to interpolate lysosomal pH from measured emission ratios.