



Jan 11, 2026

Measurement of Oxygen Consumption Rate (OCR) in iPSC-Derived Neurons and iPSCs Using Seahorse XFe96 Analyzer

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l214o3g1y/v1>

Ali Ghoochani^{1,2,3,4}, Monther Abu-Remaileh^{1,2,3,4}

¹Department of Chemical Engineering, Stanford University, Stanford, CA 94305, USA;

²Department of Genetics, Stanford University, Stanford, CA 94305, USA;

³The Institute for Chemistry, Engineering and Medicine for Human Health (Sarafan ChEM-H), Stanford University, Stanford, CA 94305, USA;

⁴Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 20815, USA

asap



Ali Ghoochani

Stanford University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l214o3g1y/v1>



Protocol Citation: Ali Ghoochani, Monther Abu-Remaileh 2026. Measurement of Oxygen Consumption Rate (OCR) in iPSC-Derived Neurons and iPSCs Using Seahorse XFe96 Analyzer. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.4r3l214o3g1y/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 22, 2025

Last Modified: January 11, 2026

Protocol Integer ID: 223019

Keywords: mitochondrial oxygen consumption rate, measurement of oxygen consumption rate, using seahorse xfe96 analyzer, oxygen consumption rate, seahorse xfe96 analyzer this protocol, seahorse xfe96 analyzer, ipsc, ocr, undifferentiated ipsc, derived neuron

Abstract

This protocol describes the procedure to measure mitochondrial oxygen consumption rate (OCR) in iPSC-derived neurons and undifferentiated iPSCs using the Seahorse XFe96 Analyzer.

Materials

- Seahorse XFe96 microplates (Agilent Technologies)
- Poly-D-Lysine (for neurons)
- rhLaminin (for iPSCs)
- Seahorse XF Base Medium (Agilent Technologies)
- DMEM/F12 (Thermo Fisher Scientific)
- Neurobasal-A (Thermo Fisher Scientific)
- MEM Non-Essential Amino Acids (Thermo Fisher Scientific)
- GlutaMAX (Thermo Fisher Scientific)
- N2 Supplement (Thermo Fisher Scientific)
- B27 Supplement with/without antioxidant (Thermo Fisher Scientific)
- NT-3 (10 ng/ml)
- BDNF (10 ng/ml)
- Mouse Laminin (1 µg/ml)
- Doxycycline hydrochloride (2 µg/ml)
- Pyruvate (1 mM)
- Glucose (20 mM)
- Oligomycin (1 µM final)
- 2,4-Dinitrophenol (DNP, 50 µM final)
- Rotenone (0.5 µM final)
- StemFlex™ Medium (Thermo Fisher Scientific)
- Seahorse XFe Report Generator
- Protein quantification reagents



Maintain iPSCs

- 1 Culture iPSCs in StemFlex™ Medium on rhLaminin-coated plates (2.5 µg/mL in 1x DPBS).
- 2 Replace medium every other day.

Initiate Pre-Differentiation (Day -3 to 0)

- 3 Detach iPSCs using ReLeSR™ and centrifuge to pellet.
- 4 Resuspend in N2 Pre-Differentiation Medium:
 - Knockout DMEM/F12
 - 1x MEM Non-Essential Amino Acids
 - 1x N2 Supplement
 - 10 ng/mL NT-3
 - 10 ng/mL BDNF
 - 1 µg/mL Mouse Laminin
 - 10 µM Thiazovivin
 - 2 µg/mL doxycycline
- 5 Seed 1.5×10^6 cells/well in rhLaminin-coated 6-well plates with 2 mL/well of medium.

Pre-Differentiation (Day -3 to 0)

- 6 Maintain cells without medium change.
- 7 On Day 0, dissociate with Accutase and centrifuge.

Neuronal Differentiation and Maturation (Days 0–21)

- 8 Plate 5×10^4 pre-differentiated cells per well into **Poly-D-Lysine-coated Seahorse XFe96 microplates.**



- 9 Culture in Classic Neuronal Medium (**1:1 DMEM/F12 and Neurobasal-A**) supplemented with:
 - 1x MEM NEAA
 - 0.5x GlutaMAX
 - 0.5x N2
 - 0.5x B27 with/without antioxidant (Based on the experiment)
 - 10 ng/mL NT-3
 - 10 ng/mL BDNF
 - 1 µg/mL Mouse Laminin
 - 2 µg/mL doxycycline
- 10 Day 7: Half-medium change, no doxycycline.
- 11 Day 14: Repeat half-medium change with same composition (no doxycycline).
- 12 Day 21: Cells ready for Seahorse assay.

Preparation for Seahorse Assay

- 13 Preparation of Sensor Cartridge (a day before experiment):

- 13.1 Hydrate the sensor cartridge:

The day prior to measuring OCR, hydrate the sensor cartridge with Seahorse XF Calibrant  200 µL /well and  400 µL PBS/ edge well and keep the sensor cartridge at  37 °C in a non-CO2 incubator overnight



Preparing the Assay Medium

- 14 On the day of the assay, replace the growth medium with pre-warmed assay medium ( 180 µL) containing Seahorse XF base medium, 1 mM pyruvate, and 20 mM glucose.
- 15 Incubate the microplate in a non-CO2 incubator at  37 °C for  01:00:00 to equilibrate the temperature and pH.



Preparing Modulating Agents

- 16 In the meantime, prepare following modulating agents (10X) in assay medium for injection:

A	B	C	D
Port Loading Design	10X concentration	Add to Port volume	Final well concentration (μM)
Port A: Oligomycin	10 μM	20 μL	1 μM
Port B : 2,4-Dinitrophenol (DNP)	500 μM	25 μL	50 μM
Port C : Rotenone	5 μM	27 μL	0.5 μM

- 16.1 Load 10X concentration of agents into port of the cartridge (From step 13)

Calibrating the Analyzer

- 17 Place the loaded sensor cartridge into analyzer and run the program. The calibration usually takes  00:15:00 .

Measuring the OCR

- 18 After calibration is done, transfer the Seahorse XFe96 cell microplate (from step 3.1) to the analyzer and press "continue"

Post-Assay Procedures

- 19 Protein Normalization:

- 19.1 After the assay, aspirate the assay medium and lyse the cells with a suitable lysis buffer and measure protein concentration using a BCA assay.



19.2 Use this data for normalizing the OCR results.

OCR Measurement in iPSCs (Undifferentiated)

20 Plate 5×10^4 iPSCs per well into rhLaminin-coated Seahorse XFe96 microplates in StemFlex™ Medium.

21 Incubate 48 h at 37°C, 5% CO₂.

22 **Seahorse Assay :**

22.1 follow same steps as in **Steps 13–19** above.